

MINI-PAM-II

Photosynthesis Yield Analyzer

Manual

2.163/10.2024

6. Edition, October 2024

MINI_PAM_II_07.docx

© Heinz Walz GmbH, 2022

Heinz Walz GmbH · Eichenring 6 · 91090 Effeltrich · Germany

☎ +49 9133 7765-0 · 📠 +49 9133 5395

✉ info@walz.com · ⭐ www.walz.com

Contents

1	Safety Instructions	7
1.1	General Safety Instructions	7
1.2	Special Safety Instructions	7
1.3	Handling of Fiberoptics	8
2	Introduction.....	9
2.1	Overview	10
3	Components and Setup	13
3.1	Extent of Delivery (Basic System)	13
3.2	Optoelectronic Unit	14
3.2.1	Batteries.....	19
3.3	2010-A Distance Clip 60°	19
3.4	Accessories	21
3.4.1	MINI-SPEC/MP Miniature Spectrometer	21
3.4.2	2035-B Leaf-Clip Holder	26
3.4.3	2054-L External LED Source	29
3.4.4	2035-B/RLC Conversion Kit for 2030-B/2035-B Leaf Clips	31
3.4.5	2060-B Arabidopsis Leaf Clip	32
3.4.6	Fiberoptics Adapter 90° 2030-B90.....	33
3.4.7	Leaf Positioning Setup DUAL-BA	33
3.4.8	2060-A Fiberoptics Holder for Surfaces.....	34
3.4.9	2065-M Mini Quantum/Temp.-Sensor	35
3.4.10	DLC-8 Dark Leaf Clip.....	35
3.4.11	KS-2500 Suspension Cuvette.....	37
3.4.12	Oxygen Package	37
3.4.13	BCS-9590 Barcode Scanner	38
3.4.14	MINI-PAM/F1 Miniature Fiberoptics.....	39
3.4.14.1	Calibration of the internal light sensor using the MINI-SPEC/MP	40
3.4.15	MQS-B/A Adapter Set for Thin Fiberoptics	41
4	Touchscreen Operation	43

Contents

4.1	Calibration.....	43
4.2	Top-level Windows	44
4.2.1	Basic Data	46
4.2.2	Primary Data.....	50
4.2.3	Porometer.....	50
4.2.4	Quenching Analysis	50
4.2.5	Ft-Chart.....	51
4.2.6	Spectrometer	52
4.2.7	Actinic + Yield	54
4.2.8	Induction Curve.....	56
4.2.9	Light Curve	57
4.2.10	Recovery.....	58
4.2.11	Actinic Light List.....	61
4.3	Main Menu	66
4.3.1	PAM Settings.....	69
4.3.1.1	Meas. Light	69
4.3.1.2	Meas. Light Sett.....	70
4.3.1.3	Gain	71
4.3.1.4	Damping	72
4.3.1.5	ETR-Factor	72
4.3.1.6	Fo' Mode.....	72
4.3.1.7	Adjust F-Offset.....	73
4.3.2	Light Sources.....	74
4.3.2.1	Actinic Light	74
4.3.2.2	Actinic Intensity.....	74
4.3.2.3	Actinic Factor	74
4.3.2.4	Far-red	74
4.3.2.5	Far-red Sett.....	76
4.3.2.6	Light Panel Sett.	76
4.3.2.7	SAT Settings.....	77
4.3.3	Program/Clock.....	78
4.3.3.1	Clock.....	78
4.3.3.2	Clock Interval	79
4.3.3.3	Clock Item.....	79

Contents

4.3.3.4	Actinic + Yield	80
4.3.3.5	Induction Curve.....	81
4.3.3.6	Light Curve	82
4.3.3.7	Yield + # Poro	82
4.3.3.8	Select Y(II), Y(NPQ), Y(NO) or Y(II), NPQ, qL	83
4.3.4	Sensors.....	83
4.3.4.1	External PAR Sensor.....	85
4.3.4.2	Porometer	85
4.3.4.3	O2/pH Sensor Settings	85
4.3.4.4	Spectrometer	88
4.3.4.5	Leaf Clip.....	91
4.3.4.6	Internal PAR	95
4.3.5	MINI-PAM-II Settings	97
4.3.6	Memory	99
4.3.7	Info	100
4.3.7.1	MINI-PAM-II Info	101
4.3.7.2	Sensor Info	101
4.3.7.3	Firmware Info	101
4.3.8	Reset	103
4.3.8.1	Reset Settings	103
4.3.8.2	Reset Sys. Settings	105
4.3.8.3	Reset Porometer.....	106
5	WinControl-3 Installation	107
5.1	Installation process	107
5.2	WinControl-3 Program Group	108
5.2.1	PAM Firmware Update	109
5.2.2	Uninstall WinControl-3.....	111
5.2.3	WinControl-3.....	111
5.2.4	WinControl-3 Network Mode.....	112
5.2.5	WinControl-3 Offline Mode	112
6	WinControl-3 Operation	113
6.1	Chart Window	114
6.1.1	Menus and Tabs	115
6.1.2	Buttons.....	118

Contents

6.1.3	Chart.....	119
6.1.4	Val./Rec./Chan.....	121
6.1.5	Results.....	124
6.1.6	Status Bar.....	124
6.2	Induct. Curve/Light Curve Windows.....	128
6.3	SAT-Chart Window.....	131
6.4	Spectrum Window.....	132
6.5	Report Window.....	136
6.6	Memory Window.....	139
6.6.1	Clear Memory.....	141
6.6.2	Download Memory.....	141
6.7	Batch Window.....	142
6.7.1	Add Command.....	142
6.7.2	Record Macro.....	144
6.7.3	Options.....	144
6.8	Control Window.....	146
6.9	Sensors Window.....	148
6.10	Settings Window.....	150
6.10.1	Title Bar.....	151
6.10.2	Light Intensity.....	151
6.10.3	Actinic Light List.....	152
6.10.4	F-Offset.....	152
6.10.5	System Parameter.....	153
6.10.6	Programs.....	153
6.10.7	Clock.....	154
6.10.8	Indicators.....	154
6.11	System Settings Window.....	155
6.11.1	Title Bar.....	157
6.11.2	Custom PAR Gain.....	157
6.11.3	Int. PAR sensor.....	157
6.11.4	Ext. PAR/T sensors.....	159
6.11.5	PAM Trim Val.....	159
6.11.6	Internal Temperature Sensor.....	160
7	Hints & Troubleshooting.....	161

Contents

7.1	Instrument Settings.....	161
7.2	Default settings	161
7.3	F ₀ Fluorescence.....	162
7.5	Signal Noise.....	164
7.6	System Hangs	164
7.7	External PAR Sensor is not Responding	164
7.8	Intensity of external lamp cannot be adjusted	164
8	Saturation Pulse Analysis.....	165
8.1	Pulse-amplitude Modulated (PAM) Fluorescence	165
8.2	Saturation Pulse Analysis	168
8.3	Relative Electron Transfer Rate (ETR)	172
8.3.1	Reviews on Saturation Pulse Analysis of Photosystem II.....	173
9	Specifications MINI-PAM-II	177
9.1	Basic System	177
9.1.1	Optoelectronic Unit	177
9.1.2	Fiberoptics MINI-PAM/F	179
9.1.3	Power Supply MINI-PAM-II/N	179
9.1.4	Battery Charger 000190101101	180
9.1.5	Distance Clip 60° 2010-A	180
9.1.6	Complementary Items.....	180
9.1.7	Software WinControl-3.....	180
9.1.8	Transport Case MINI-PAM/T	181
9.2	Accessories	181
9.2.1	2035-B Leaf-Clip Holder	181
9.2.2	Fiberoptics Adapter 90° 2030-B90	182
9.2.3	External LED Source 2054-L	182
9.2.4	Dark Leaf Clip DLC-8.....	183
9.2.5	Arabidopsis Leaf Clip 2060-B	183
9.2.6	Conversion Kit for 2030-B/2035-B Leaf Clips 2035-B/RLC	184
9.2.7	Fiberoptics Holder for Surfaces 2060-A.....	184
9.2.8	Mini Quantum/Temp.-Sensor 2065-M	185
9.2.9	Miniature Spectrometer MINI-SPEC/MP	185
9.2.10	Flat Entrance Optics SPEC/P	186
9.2.10.1	Fluorescence and Reflectance Optics SPEC/R.....	186

Contents

9.2.11	PAR Calibration Block 000160101439	186
9.2.12	Suspension Cuvette KS-2500.....	187
9.2.13	Magnetic Stirrer with Fiberoptics Holder MKS-2500.....	187
9.2.14	Compact Tripod ST-2101A.....	187
9.2.15	Barcode Scanner BCS-9590	188
9.2.16	MINI-PAM/F1 Miniature Fiberoptics.....	188
10	Guarantee	189
10.1	Manufacturer's Guarantee	189
10.2	Conditions.....	189
10.3	Instructions	190
10.4	Applicable law	191
11	Index	193
12	List of Figures.....	201
13	List of Tables.....	205

1 Safety Instructions

1.1 General Safety Instructions

- Read safety instructions and the operating instructions prior to operation of the device and its accessories.
- Pay attention to all safety warnings.
- Keep device and its accessories away from water or high moisture areas.
- Keep the device and its accessories away from dust, sand, and dirt.
- Do not put the device and its accessories near sources of heat.
- Ensure that neither liquids nor foreign bodies get inside the device or its accessories.
- Ensure sufficient ventilation.
- Connect the device only to the power source indicated in the operating instructions or on the device. If the device is not in use, remove the mains plug from the socket.
- The device and its accessories should only be repaired by qualified personnel.

1.2 Special Safety Instructions


- The MINI-PAM-II is a highly sensitive instrument which should be only used for research purposes, as specified in this manual. Follow the instructions of this manual to

avoid potential harm to the user and damage to the instrument.

- The MINI-PAM-II can emit very strong light! To avoid harm to your eyes, never look directly into the light port of the MINI-PAM-II or its fiberoptics.
- Switch off MINI-PAM-II before connecting or disconnection 2054-L External LED Source.

1.3 Handling of Fiberoptics

The MINI-PAM/F Fiberoptics of the MINI-PAM-II is made of delicate glass fibers. Each glass fiber extents over the entire length of the fiberoptics. Do not sharply bend, stretch, or crush the fiberoptics as this can break the glass fibers. Each broken fiber reduces light transmission by the fiberoptics. If many fibers are broken, the fiberoptics must be replaced. Place protection caps on fiber ends when the fiberoptics is not in use.

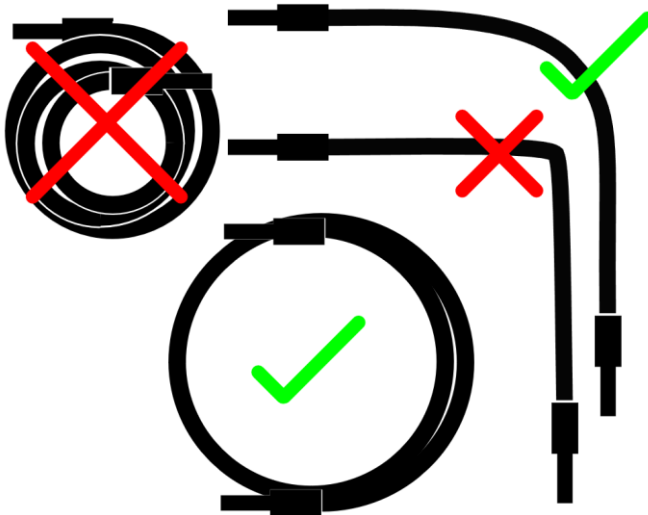

2 Introduction

- The “Photosynthesis Yield Analyzer MINI-PAM-II” has been designed for highly sensitive saturation pulse analysis of photosystem II (PS II) in the field as well as in the laboratory. The automatically calculated parameters are F_V/F_M (maximum photochemical yield), $Y(II)$ (effective photochemical yield) and its complementary yields $Y(NPQ)$ and $Y(NO)$, as well as parameters of photochemical (q_L , q_P) and non-photochemical quenching (q_N , NPQ) (see Table 29, page 174).
- The instrument continues the tradition of the preceding MINI-PAM chlorophyll fluorometer. The major technical advancements of the MINI-PAM-II are the consistent use of energy-efficient LEDs, an internal PAR sensor, and stand-alone operation by a touchscreen which is well readable under natural light conditions. Also, a far-red LED has been added for selective excitation of photosystem I.
- A further technical progress is the newly designed leaf clip sensor (2035-B) which measures photosynthetically active radiation (PAR) at leaf level with high accuracy and, thus, provides reliable light intensity data for calculations of electron transport rates (ETR).
- A variety of add-ons make the MINI-PAM-II a highly versatile measuring system which can be configured to meet the needs of your research goal. The range of accessories includes a multi-colored lighting unit (Section 3.4.3, page 29), an optical oxygen sensor (Section 3.4.12, page 37) and a miniature spectrometer (Section 3.4.1, page 21).
- For long-term field campaigns, the memory capacity has been upgraded to keep data of more than 27,000 saturation pulse

analyses. The fluorometer is powered by of-the-shelf AA (Mignon) batteries which are easily replaceable even under field conditions. One set of batteries lasts for up to 1000 saturation pulse analyses.

- The MINI-PAM-II can be operated in the stand-alone mode or by the well-proven WinControl-3 software. WinControl-3 has been introduced with the JUNIOR-PAM fluorometer and now operates many other fluorometers like the DIVING-PAM, MONITORING-PAM, and WATER-PAM. In addition to the features available in the stand-alone mode, the software allows evaluations of light-response curves by a non-linear fitting routine and automatic execution of custom-designed experiments using the built-in batch file feature.

2.1 Overview

The MINI-PAM-II fluorometer provides a vast range of settings and protocols for measuring fluorescence. To make full use of these opportunities, become acquainted with terminology and principles of saturation pulse analysis. The present manual provides a chapter dealing with the basics of saturation pulse analysis (Chapter 8, page 165).

Chapter 8 also provides a short list of review papers on PAM chlorophyll fluorescence and saturation pulse analysis (page 173). Help with technical problems is provided by Chapter 7 (page 161).

In the field, the MINI-PAM-II is mostly operated in stand-alone mode via its touch screen. Chapter 4 (page 43) provides detailed instructions on how to use the touchscreen interface including advice on fluorescence induction and light curve programs.

In the laboratory, the WinControl-3 software is normally used to operate the MINI-PAM-II. Chapters 5 and 6 (page 103 and 113, respectively) introduce the features and functions of the WinControl-3 software.

This manual also includes sections on safe handling of the MINI-PAM-II (Chapter 1, page 7), and on the extent of delivery of the basic fluorescence system and its accessories (Chapter 3, page 13). Technical information is provided by Chapter 9 (page 177), and guarantee conditions are given in Chapter 10 (page 189).

The manual includes a keyword index (Chapter 11, page 193), as well as a list of the illustrations and tables used (Chapter 12, page 201, and Chapter 13, page 205, respectively).

3 Components and Setup

3.1 Extent of Delivery (Basic System)

Item	Order Code
Optoelectronic Unit	MINI-PAM-II/B or -/R
Fiberoptics	MINI-PAM/F
Power Supply	MINI-PAM-II/N
Battery charger	190101101
Battery (12 x)	160101990
USB Cable Type A to Mini B	130606252
Distance Clip 60°	2010-A
Sloped Plexiglas Rack	240313614
Stylus	160201311
Carrying Strap	150401922
Software WinControl-3	
MINI-PAM-II Manual	

3.2 Optoelectronic Unit

Sockets, fiberoptics port and on/off switch of the rear side of the MINI-PAM-II are outlined and numbered in Fig. 1, their properties and functions are explained in Table 1, page 15. Fastening of the carrying belt is illustrated in Fig. 2, page 16.

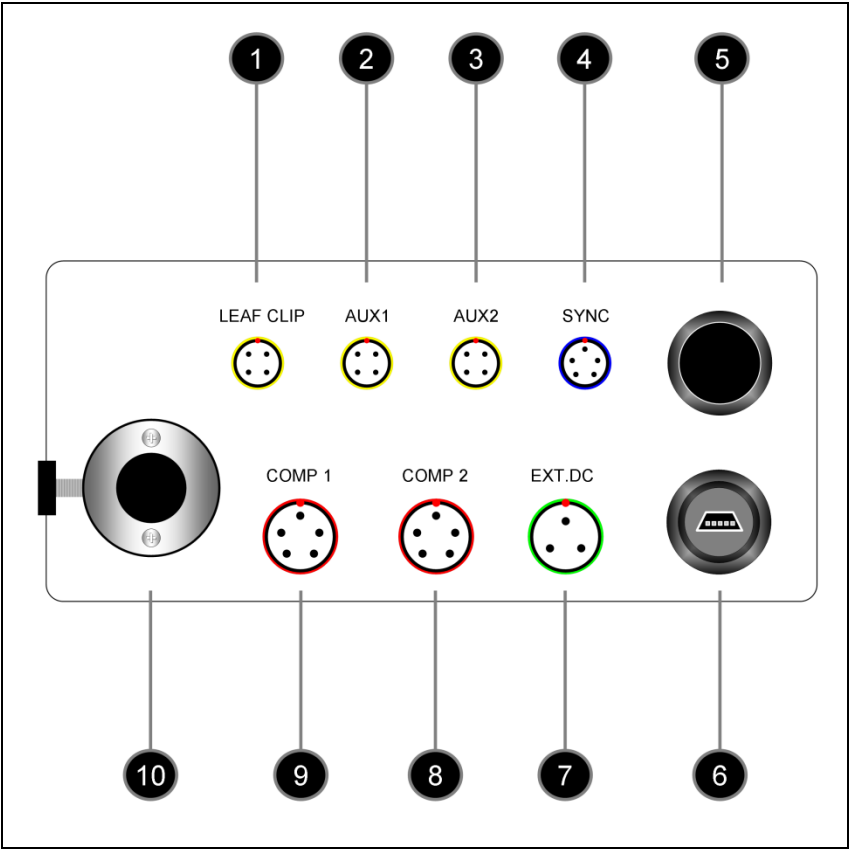


Fig. 1: Rear Panel of MINI-PAM-II

Table 1: Rear Panel of MINI-PAM-II.

Numbering refers to Fig. 1, page 14	Function
1	LEAF CLIP Socket for 2035-B Leaf-Clip Holder
2	AUX 1 Electronically configured as LEAF CLIP socket
3	AUX 2 Electronically configured as LEAF CLIP socket
4	SYNC Socket for external light source which emits synchronized with MINI-PAM-II measuring light
5	ON/OFF MINI-PAM-II switch
6	USB SOCKET Receptacle for MINI-B USB plug.
7	EXT. DC Socket for Power Supply MINI-PAM-II/N
8	COMP 2 Prepared for peripherals mastering RS232 communication (e.g. bar code scanner)
9	COMP 1 Electronically configured as COMP 2
10	LIGHT PORT Port for Fiberoptics MINI-PAM/F

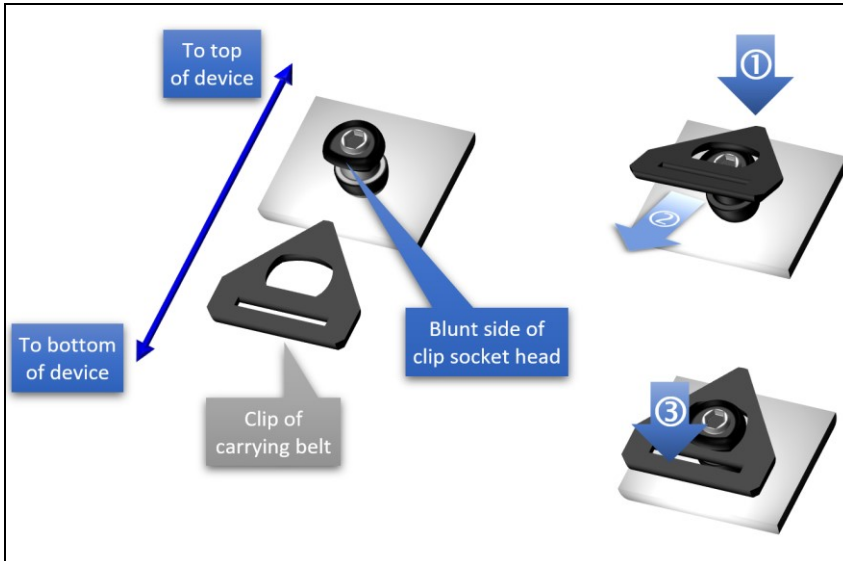


Fig. 2: Carrying Belt Fastening

Note Noses for the clips of the carrying belt are mounted on each side of the MINI-PAM-II device.
From top to bottom, the nose consists of a bluntly shaped head, a circular groove, and a cylindrical base.


Instruction

- ① Position apex of belt clip above the groove on the opposite of the blunt site.
- ② While keeping the position of the apex, pull base of the belt clip to further move the clip apex into the groove.
- ③ While maintaining the pull, press down the base of the clip.

Note Great caution should be exercised to prevent dirt or foreign matter from entering the ports or sockets of the MINI-PAM-II. Do not force a plug into the wrong socket. Orientate each plug so that the red dot on the plug coincides with the red dot of the socket. Do not try to disconnect a plug by pulling at the cable. Disconnect plug by pulling at the rippled bushing of the plug.

Fig. 3: Self-locking Connector

To lock connection, align red dots of plug and socket and push in. To disconnect, hold the knurled sleeve and pull out. Applies for connection numbers 1 - 4 and 7 - 9 Fig. 1.

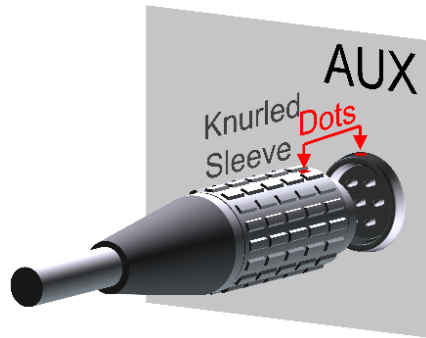

Marked End of Fiberoptics



Fig. 4: Fiberoptics

Insert the end marked "TO= PAM" into fiberoptics port.

Two versions of the MINI-PAM-II fluorometer are available. The MINI-PAM-II/B fluorometer is equipped with a blue LED which in the MINI-PAM-II/R fluorometer is replaced by a red LED. This LED is electronically driven to act as measuring and as actinic light source. In addition, both versions of the MINI-PAM-II offer a far-red LED. Normalized spectra of blue, red, and far-red LEDs are shown in Fig. 5.

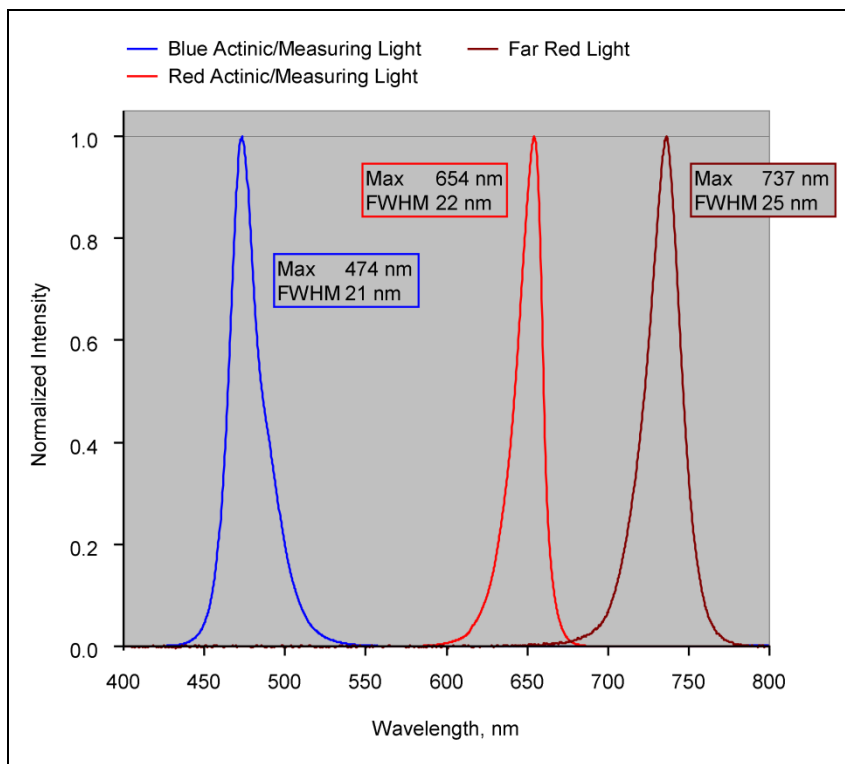


Fig. 5: Emission Spectra of MINI-PAM II LEDs

Normalized emission spectra of blue LED (MINI-PAM-II/B), red LED (MINI-PAM-II/R), and far-red LED (MINI-PAM-II/B and R).

3.2.1 Batteries

When operated independently, the system is powered by six AA (Mignon) rechargeable batteries (1.2 V/2 Ah). The system can also be powered by non-rechargeable batteries. The battery compartment of the opto-electronic unit does not have a charging function. Therefore, the device can be connected to line power even in the presence of non-rechargeable batteries.

The battery compartment is closed by an aluminum plate. Its locking mechanism functions properly if the label “INNER FACE” on the aluminum plate faces the batteries.

3.3 2010-A Distance Clip 60°

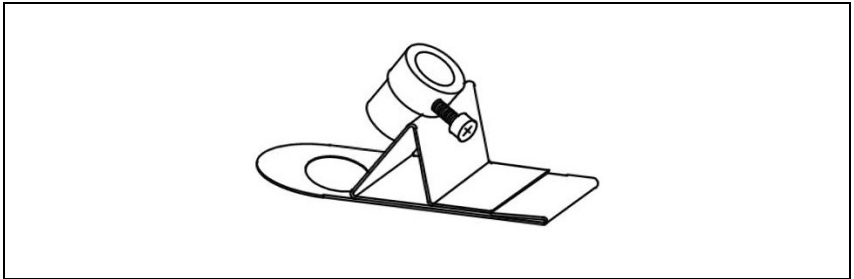


Fig. 6: Distance Clip 60° 2010-A

The 2010-A clip positions the fiberoptics end-piece relative to the sample. The axis of the end-piece is positioned at a 60° angle relative to the sample plane. Two different spacer rings may be used to increase the distance between fiberoptics and sample.

In case of relative thick leaves, or when lichens and mosses are examined, the sample may be placed below the hole of the 2010-A clip. Normal leaves are usually examined above this hole. In the

latter case, the leaf can be held between the folded parts of the clip.

The distance between fiberoptics exit plane and sample has considerable influence on signal amplitude and effective light intensities (Fig. 7, page 20). With a 60° angle between sample plane and fiberoptics, the distance between leaf surface and fiberoptics varies. Hence, the leaf surface is exposed to slightly heterogeneous light intensities when actinic light is applied *via* the fiberoptics. A much more pronounced intensity gradient exists inside the leaf due to shading by the top chloroplast layers. In essence, the measured signal will be dominated by that part of the leaf which receives maximal intensity, as this also is most strongly excited by the measuring light and emits most of the fluorescence which is received by the fiberoptics.

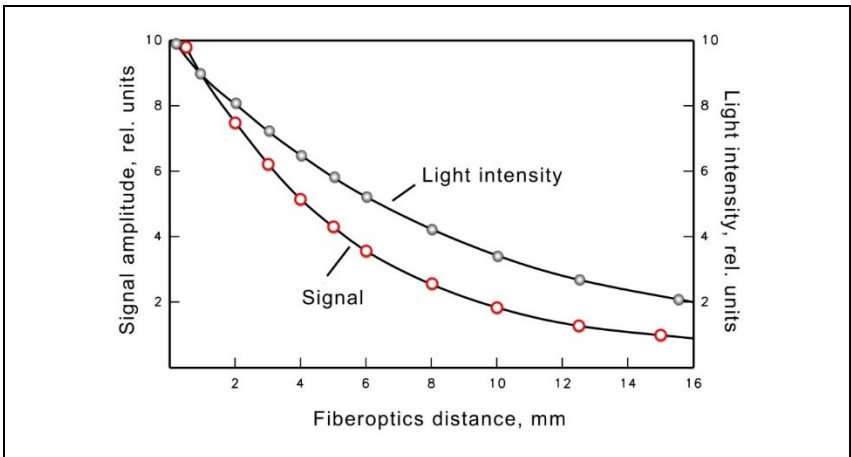


Fig. 7: Measuring Light and Signal

Signal amplitude and light intensity as a function of the distance between fiberoptics and sample level

3.4 Accessories

3.4.1 MINI-SPEC/MP Miniature Spectrometer

Originally, the spectrometer MINI-SPEC/MP (Fig. 11, page 25) has been introduced as accessory for the DIVING-PAM-II. For this reason, the MINI-SPEC/MP possesses an underwater-type connector. Proper connection to the MINI-PAM-II requires that the cable plug is completely inserted before the screw is tightened (see Fig. 8).

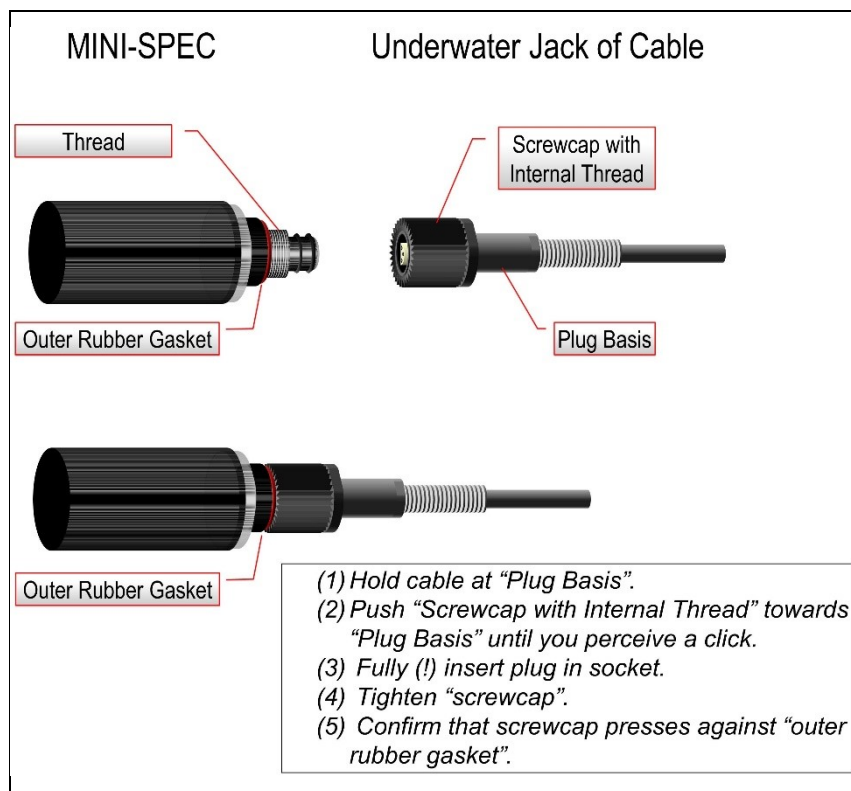


Fig. 8: Connection of MINI-SPEC/MP

The spectrometer is calibrated to measure spectra of quantum fluxes. Integration of these spectra over the visible range (400 – 700 nm) yields PAR data equivalent to those recorded by Walz quantum sensors.

Like the PAR sensor of the 2035-B or 2065-M devices (Section 3.4.1, page 21 and Section 3.4.11, page 37), the spectrometer can be employed to calibrate the internal PAR sensor of the MINI-PAM-II (see Section 4.3.4.6, page 95). To this aim, the MINI-PAM-II light guide and the entrance optics of the spectrometer are inserted in the PAR calibration block (Fig. 9). The light guide can be inserted either in the 60° or the 90° port according to the two possible orientations of the light guide in the Leaf Clip Holder 2035-B. With both pieces fully inserted, the distance between fiberoptics end and diffusing disk of the spectrometer matches the corresponding standard distances between fiberoptics end and sample level in the Leaf Clip Holder 2035-B.


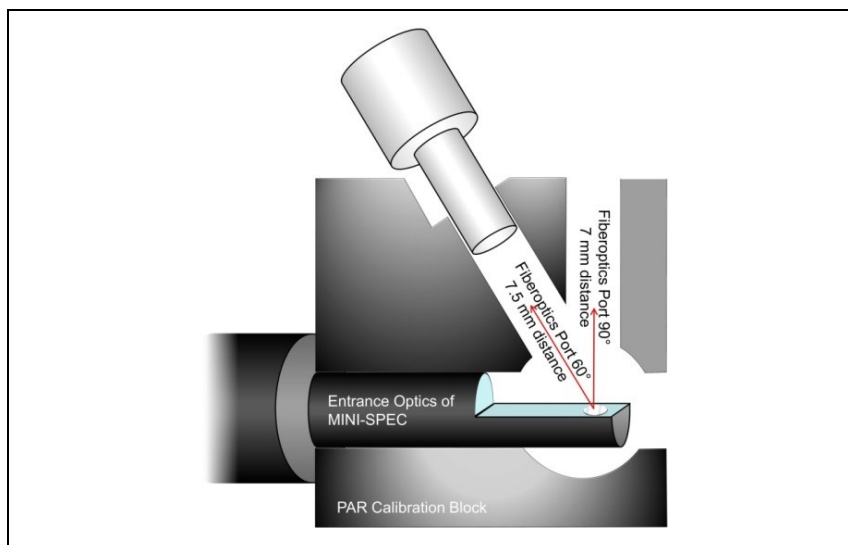


Fig. 9: PAR Calibration Block

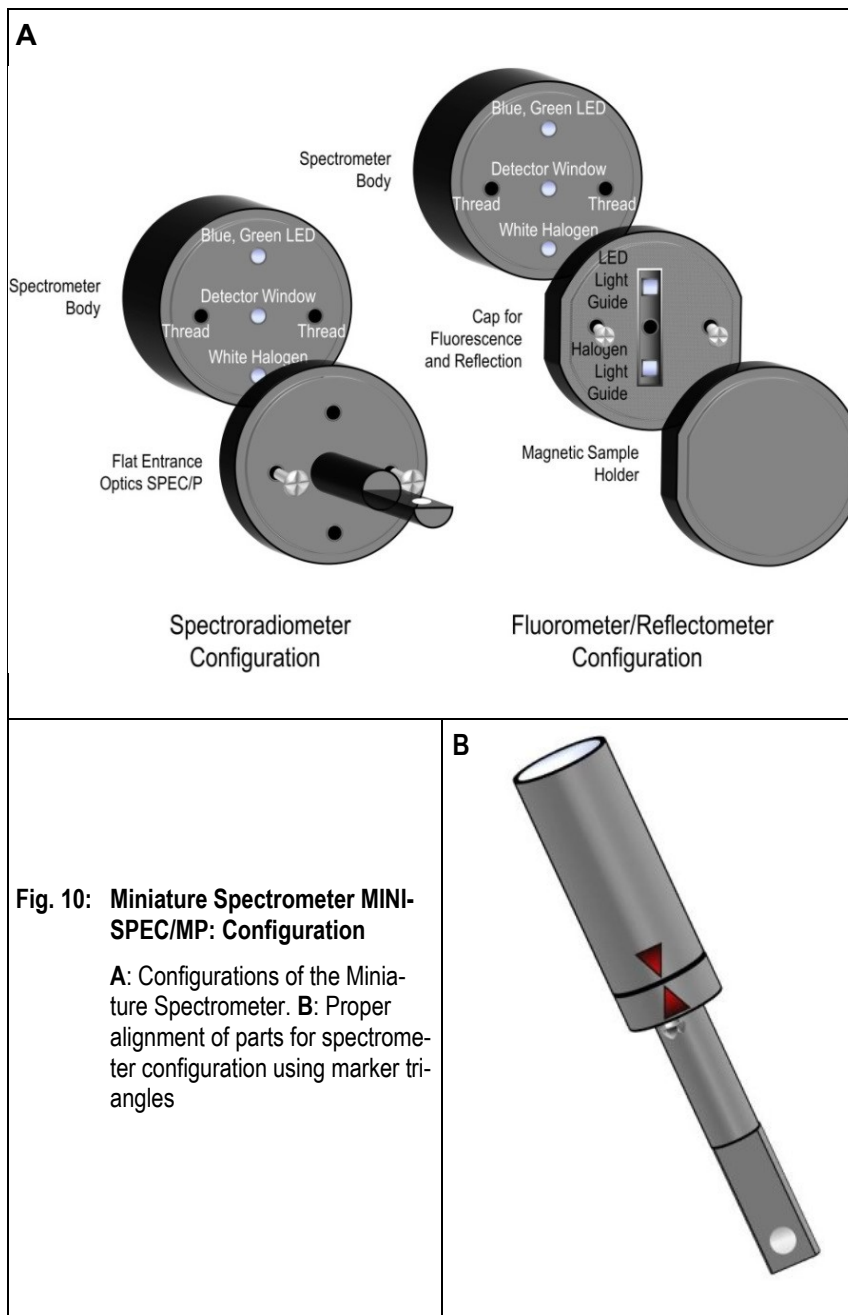

Replacing the light entrance optics by the cap for fluorescence and reflectance (Fig. 10, page 23) extends the range of spectral information attainable by the miniature spectrometer.

In the fluorescence mode, a light guide in the cap leads light from a blue or a green LED located inside the spectrometer body to the sample. The excited fluorescence reaches the spectrometer body through a central hole in the cap. Green light penetrates deeper into photosynthetic tissue than blue or red light (Terashima et al (2009) *Plant Cell Physiol* 50:684-697). Therefore, the two excitation colors permit probing the effect of varying depths of penetration of excitation light on the shape of fluorescence spectra. The spectrometer automatically compensates for the spectral sensitivity of the detector.

For reflectance measurements, a halogen lamp is employed whose emission is transferred through another light guide to the sample (Fig. 10, page 23).

From reflectance measurements, reflectance (R) is derived by dividing a sample spectrum (r_{sample}) by the spectrum of a white light-diffusing reflector consisting of a fluoropolymer with very high diffuse reflectance in the entire visible range ($r_{\text{reference}}$):

$$R(\lambda) = \frac{r_{\text{sample}}(\lambda)}{r_{\text{reference}}(\lambda)}$$

The reference material is part of delivery. To maintain its optical properties, do not touch the white surface, keep away dirt, dust, and humidity, and store the reflectance standard in a closed container when not used.


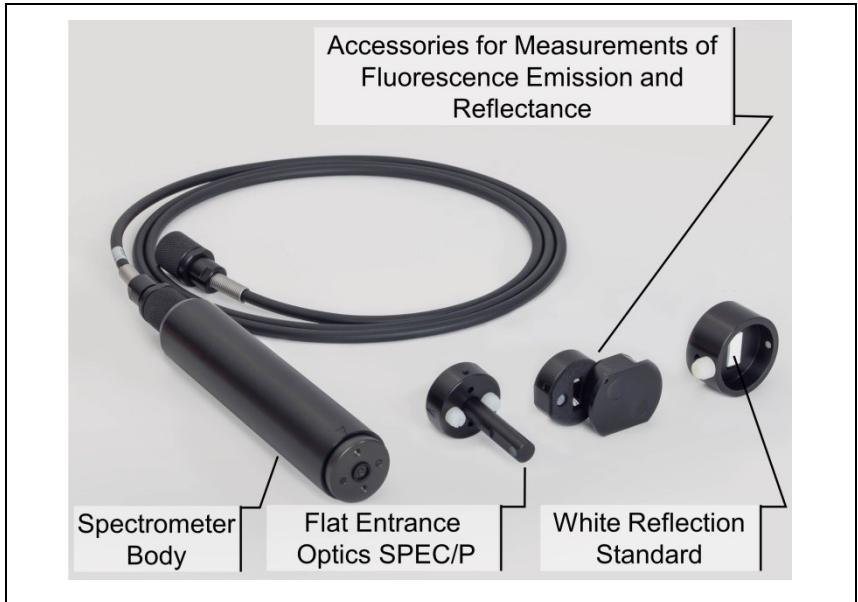


Fig. 11: MINI-SPEC/MP

In addition to the items shown, the MINI-SPEC/MP includes a mounting ring (order number 000244905514) into which the spectrometer can be inserted.

Fig. 12: Emission Spectra of Miniature Spectrometer Light Sources

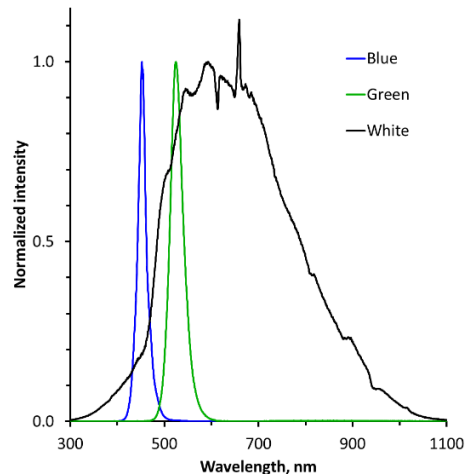

3.4.2 2035-B Leaf-Clip Holder

The Leaf-Clip Holder 2035-B must be connected to the LEAF CLIP socket (Fig. 1, page 14) to record PAR, leaf temperature and ambient humidity in parallel with chlorophyll fluorescence. In the stand-alone mode, readings of environmental data are taken with every saturation pulse analysis, but these data can be continuously recorded when the MINI-PAM-II is operated by the WinControl-3 software.

The Leaf-Clip Holder 2035-B is almost indispensable for field investigations, when ambient conditions may vary considerably. It substitutes for the standard "Distance Clip" (2010-A) as a device for defined positioning of the fiberoptics relative to the leaf plane.

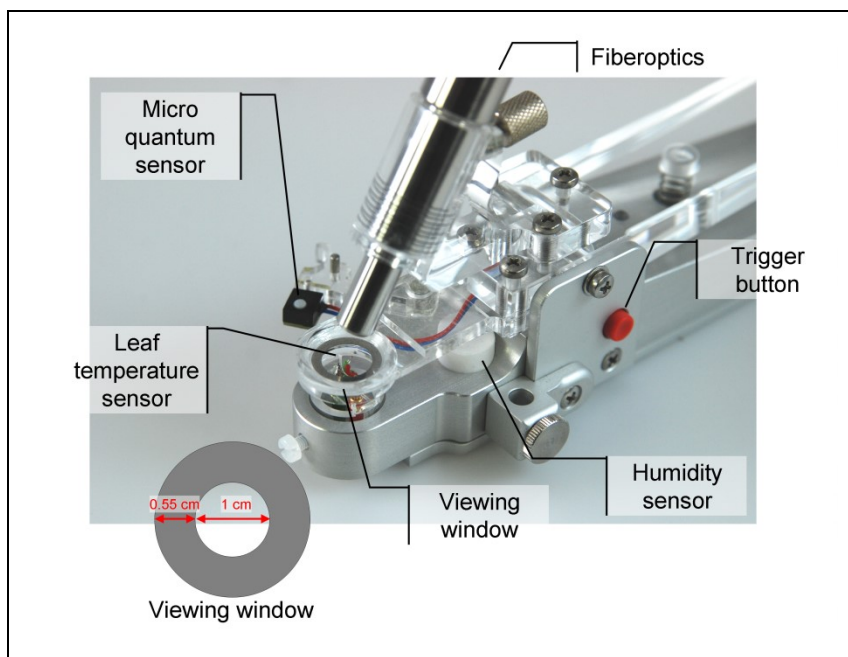


Fig. 13: 2035-B Leaf-Clip Holder

Further, using the PAR sensor of the leaf clip, the internal PAR sensor of the MINI-PAM-II can be readily calibrated.

In the 2035-B holder, the leaf is resting on a Perspex tube with widened crest. The tube can be vertically adjusted to account for different leaf thicknesses. The fiberoptics axis forms a 60° angle with the leaf plane. Optionally, a 90° fiberoptics adapter (2030-B90) is available for applications requiring homogenous illumination by actinic light applied *via* the fiberoptics (see Fig. 17, paged 32). The distance between fiberoptics and leaf can be varied. For most applications the minimal distance is recommended (maximal signal). Larger distances can be defined by spacer rings. The illuminated leaf area is limited by a steel ring with 10 mm Ø opening.

At the bottom of the Leaf-Clip Holder 2035-B, a tripod mounting thread is provided. Mounting the device on a tripod (e. g. Compact Tripod ST-2101A) facilitates long term measurements with the same plant.

The handle of the Leaf-Clip Holder 2035-B features a red push-button for remote control of the MINI-PAM-II. Pressing the button triggers a saturation pulse and associated measurements of fluorescence levels for “fluorescence quenching analysis”.

Micro-Quantum-Sensor

A micro quantum sensor is integrated into the Leaf-Clip Holder 2035-B to monitor the photosynthetic active radiation (PAR, between 400 and 700 nm) to which the sample is exposed. The micro-quantum-sensor measures light intensity in $\mu\text{mol quanta m}^{-2} \text{s}^{-1}$. The $\mu\text{mol quanta m}^{-2} \text{s}^{-1}$ is the unit of photon flux density. Hence, the micro-quantum-sensor measures photosynthetic photon flux density (PPFD).

The optoelectronic elements of this micro-quantum-sensor are:

- A 3 mm \varnothing diffusing disk.
- High stability silicon photovoltaic detector with filter set for PAR correction, magnetically attached to 2035-B Leaf Clip Holder.
- Cosine response characteristics (Angular dependence: Error < 3 % for angle between -30 ° and +30 ° from normal axis).

The sensor is factory calibrated and calibration factors are stored in the internal memory of the 2035-B leaf clip. The stability of calibration depends on keeping the diffuser clean. It is advisable to check calibration regularly by comparison with a standard quantum sensor. Any deviation can be corrected by entering a new calibration factor in WinControl-3 (Section 6.11.4, page 159) or on the touch screen (Section 4.3.4.1, page 85). A substantial increase of the calibration factor from its original value indicates dirt-deposition on the diffuser, which may be reversed by gentle cleaning using a cotton tip applicator, moistened with some diluted ethanol.

Thermocouple

A NiCr-Ni thermocouple is mounted in the Perspex tube on which the leaf area is resting. The thermocouple is forming a loop that gently presses against the lower surface of the leaf. This arrangement results in effective temperature equilibration between leaf and thermocouple and protects the thermocouple from direct sun radiation.

The reference couple is located on the circuit board, near the thermovoltage amplifier, enclosed in the bottom part of the holder. The relationship between thermovoltage and temperature is almost linear. With decreasing temperatures there is a small decline of $\Delta V / ^\circ\text{C}$. Calibration was performed at 25 °C. At 0 °C or -15 °C the deviation amounts to 0.5 or 0.8 °C, respectively.

Humidity Sensor

A calibrated, capacitive-type humidity sensor measures humidity conditions close to the sample surface.

Data Display

All sensor data are displayed on the touchscreen window “Primary Data” (Fig. 31, page 49).

Signal LED

See Table 2.

Table 2: Signal Code of LED on 2035-B Leaf Clip Holder.

LED action	Status
Flashing green	Normal operation.
Continuous green	Communication from MINI-PAM-II to 2035-B clip interrupted. This happens temporarily during firmware update of MINI-PAM-2.
Flashing red	(a) Broken thermocouple: inspect and ask for repair kit. (b) Internal error on 2035-B EPROM: contact Walz.
Continuous red	As “flashing red” plus communication interrupted.

3.4.3 2054-L External LED Source

For experiments requiring different actinic light colors, we offer an external light source which can be attached to the 2035-B leaf clip (Fig. 14, page 30). The light source consists of four four-chip LED RGBW sources each capable of emitting red, green, blue, and white light. Total intensity and color composition can be regulated by the software of the MINI-PAM-II or by WinControl-3. The maximum PAR of each light quality is $1500 \mu\text{mol m}^{-2} \text{s}^{-1}$.

Note Switch off MINI-PAM-II before connecting or disconnecting 2054-L External LED Source.

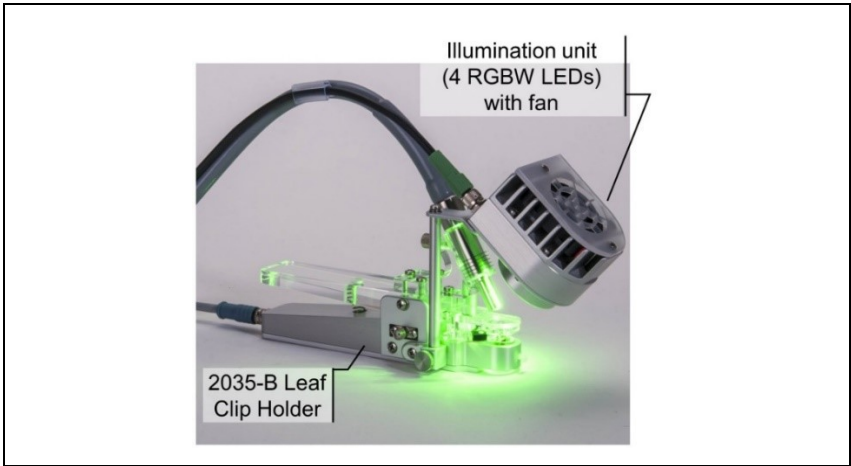


Fig. 14: 2054-L External LED Source

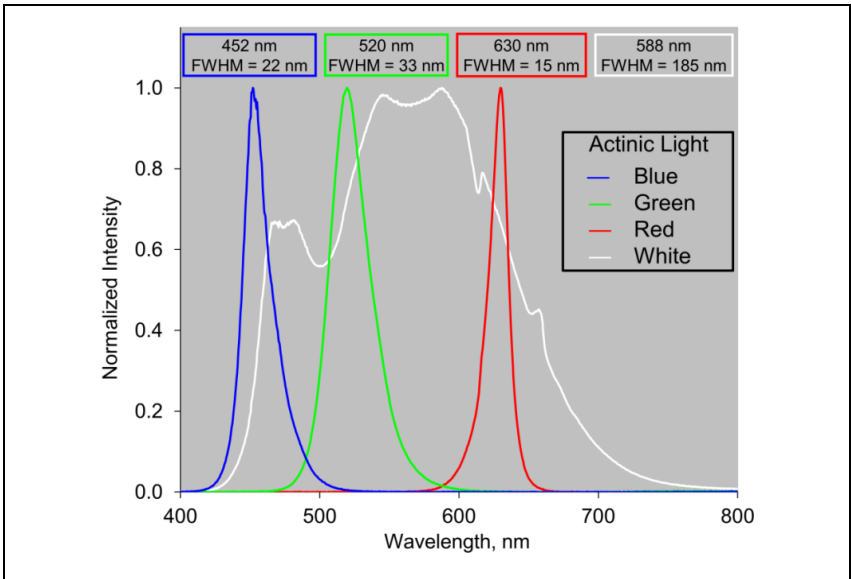


Fig. 15: Emission spectra of the 2054-L External LED Source

3.4.4 2035-B/RLC Conversion Kit for 2030-B/2035-B Leaf Clips

The 2035-B/RLC kit shields the sample area of 2030-B/2035-B Leaf Clips from external light (Fig. 16). The accessory has been designed for field experiments with the fluorometer's internal light source. The 2035-B/RLC kit is especially suitable for rapid light curve experiments (RLC) in outdoor studies.

See Fig. 17 (page 32) for instruction to mount the fiberoptics duct "Fiberoptics Adapter 90° 2030-B90".



Fig. 16: 2035-B/RLC Conversion Kit

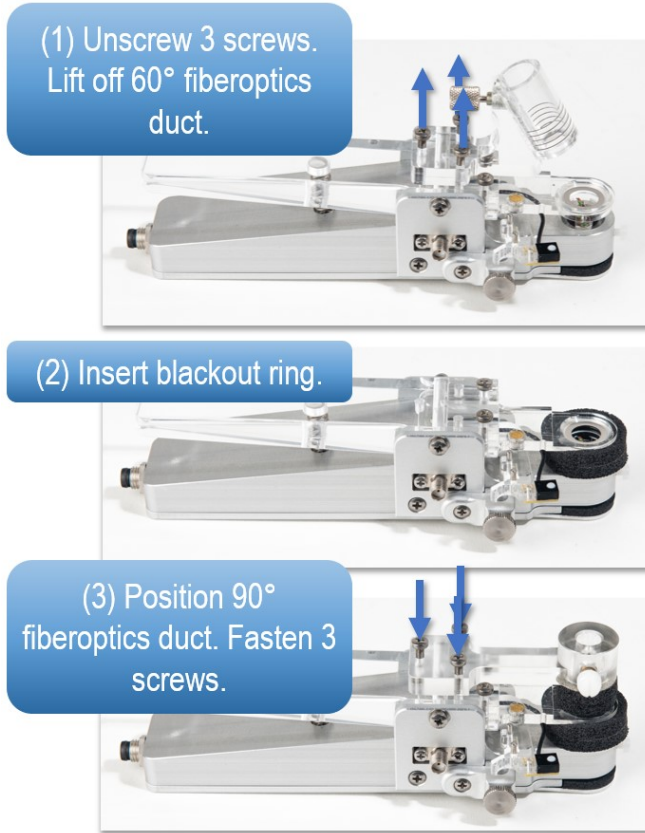


Fig. 17: Fiberoptics Adapter 90° 2030-B90: Mounting

3.4.5 2060-B Arabidopsis Leaf Clip

This leaf clip is made to position small samples in the beam of the fiberoptics of the MINI-PAM-II. Usually, the 2060-B clip is combined with the 2065-M Mini Quantum/Temp.-Sensor to measure PAR at sample level and lower leaf temperature (see Fig. 18, page 33).

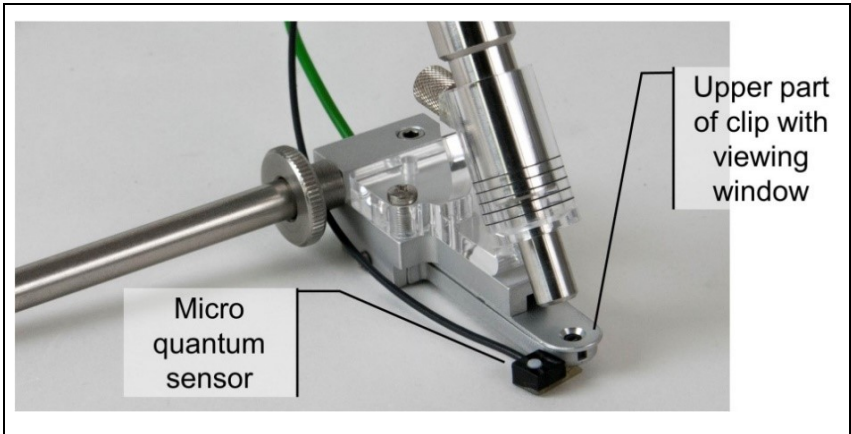


Fig. 18: 2060-B & 2065-M Arabidopsis Clip with Quantum/Temp.-Sensor

3.4.6 Fiberoptics Adapter 90° 2030-B90

The fiberoptics adapter 90° can be attached to the leaf clips 2035-B and 2060-B to position the fiberoptics of the MINI-PAM-II at 90° angle relative to leaf plane. See Section 3.4.4 for instructions on mounting the Fiberoptics Adapter 90° 2030-B90.

3.4.7 Leaf Positioning Setup DUAL-BA

The DUAL-BA has been designed to position the fiberoptics of a PAM fluorometer easily and quickly on attached leaves. The accessory is compatible with the DUAL-PAM/F, MINI-PAM-II and PAM-2500 fluorometers. It includes a duct with laterally mounted neodymium magnet and a screw to lock the fiberoptics tip. An angle bracket made of a spring steel strip positions the leaf in front of the fiberoptics, where the steel strip is held by the magnet of the duct. The DUAL-BA includes a fiberoptics guide with metal rod to mount the guide on a stand.



Fig. 19: Leaf Positioning Setup DUAL-BA

3.4.8 2060-A Fiberoptics Holder for Surfaces

The holder positions the fiberoptics of the MINI-PAM-II on bulky samples. When combined with the 2065-M Mini Quantum/Temp.-Sensor, temperature and PAR of the surface can be measured (Fig. 20).

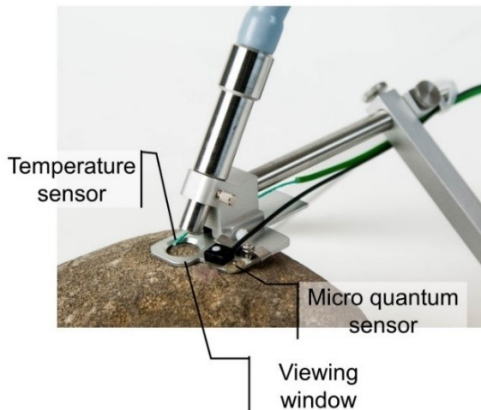


Fig. 20: 2060-A & 2065-M, Surface Holder with Quantum/Temp.-Sensor

3.4.9 2065-M Mini Quantum/Temp.-Sensor

The light and temperature sensors of the 2065-M device can be mounted on both the 2060-B Arabidopsis Leaf Clip and the 2060-A Fiberoptics Holder for Surfaces. Sensors and electronics of the 2065-M device are identical to the 2035-B clip (Section 3.4.1, page 21).

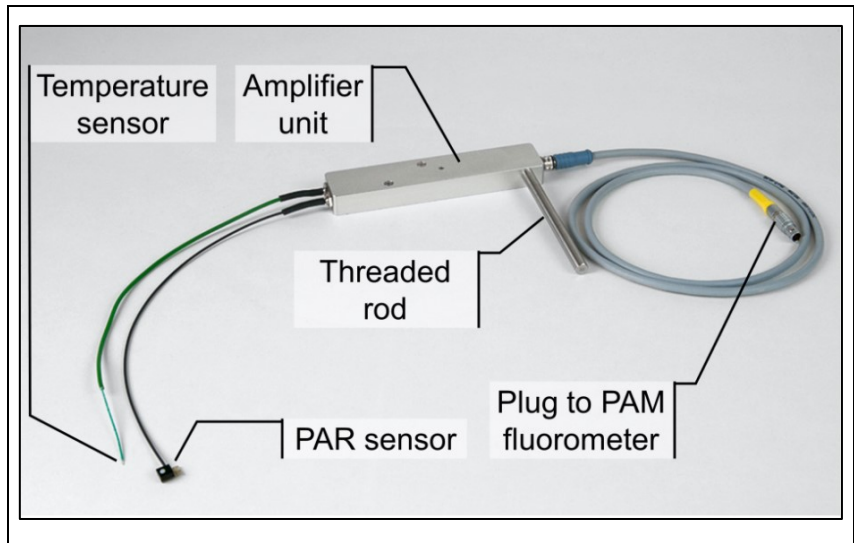


Fig. 21: Mini Quantum/Temp.-Sensor 2065-M

3.4.10 DLC-8 Dark Leaf Clip

The DLC-8 leaf clip permits dark-acclimation of small leaf areas in the field which is essential for proper determination of the maximal quantum yield F_v/F_m and for recording of dark-light induction kinetics. The Dark Leaf Clip DLC-8 weighs approx. 4 g and, hence, can be attached to most types of leaves without any detrimental effects.



Fig. 22: DLC-8 Dark Leaf Clip

The tip of the fiberoptics of the MINI-PAM-II fits snugly into the DLC-8 port. With the fiber tip inserted, the sliding shutter of the DLC-8 can be opened so that F_0 and F_M level fluorescence can be measured without interference of ambient light.

Using the Dark Leaf Clip DLC-8, the fiberoptics is positioned at right angle with respect to the leaf surface at the relatively short distance of 7 mm. Consequently, signal amplitude is distinctly higher (factor of 2.4) compared to the Leaf-Clip Holder 2035-B with 60° fiberoptics angle. To avoid signal saturation, the settings of measuring light intensity and gain have to be lowered with respect to the standard settings (Fig. 45, page 69).

When the shutter is still closed and the measuring light is on, an artifactual F_t signal is observed. This signal is due to a small fraction of measuring light which is reflected from the closed shutter to the photodetector. However, this background signal is of no concern as the reflection is much smaller when the shutter is opened, and the measuring light is strongly absorbing by the leaf sample instead of being reflected by the metal surface of the shutter.

3.4.11 KS-2500 Suspension Cuvette

The suspension cuvette includes a 400 μ l sample compartment made of stainless steel with POM exterior. The cuvette is equipped with a 7 mm fiberoptics window adapter, an injection port for microliter syringes, and nozzles for connecting an external flow-through water-bath for temperature control.

NOTE: Avoid corrosion by thoroughly cleaning and drying the inside of the cuvette.



Fig. 23: KS-2500 Suspension Cuvette and MKS-2500 Magnetic Stirrer

3.4.12 Oxygen Package

The oxygen package consists of an optode sensor, an oximeter (FSO2-1) and an interface (FSO2-AK). The sensor is inserted through the needle port of the KS-2500 cuvette (Fig. 23, Fig. 24). The oximeter controls the sensor and transmits the oxygen signal via the interface to the MINI-PAM-II. A separate manual is provided for this package.



Fig. 24: O₂ measurement. System Overview

3.4.13 BCS-9590 Barcode Scanner

The barcode scanner is the ideal add-on when many differed samples are probed. Simply convert your sample IDs into barcodes, print labels, and mark your samples. Then, the BCS-9590 Barcode Scanner writes your sample ID into the memory of the MINI-PAM-II or the report data in WinControl-3.

In Report table, barcodes are printed in the Aux1 column in the data line of the subsequently executed saturation pulse analysis. Writing barcode and measuring results in the same line facilitates data sorting when data of many samples are collected.



Fig. 25: **BCS-9590 Barcode Scanner**

3.4.14 MINI-PAM/F1 Miniature Fiberoptics

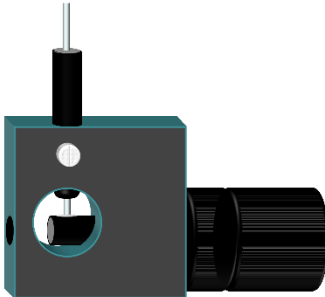
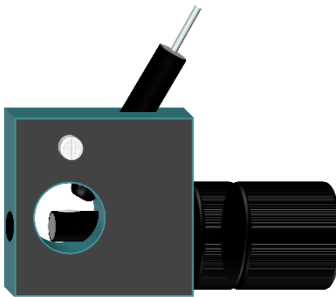
The MINI-PAM/F1 fiberoptics is useful for investigation of small areas. It consists of a single coated plastic fiber with an active diameter of 2 mm.



Fig. 26: **Miniature Fiberoptics MINI-PAM/F1**

3.4.14.1 Calibration of the internal light sensor using the MINI-SPEC/MP

Table 3: Miniature Fiberoptics MINI-PAM/F1

What is measured?	PAR at fiberoptics tip	PAR at sample level (Standard geometry, distance to sample level 7.5 mm, angle 60°. See Fig. 9.)
Setup	 <p>Fiber touches center of the diffusing disk of the MINI-SPEC.</p>	 <p>Fiber flush with adapter which is fully inserted in 60° port of the calibration block.</p>
Calibration of internal light sensor	<p>Spec. Cal. Factor = 3.01.</p> <p>See 4.3.4, page 83. The Spec. Cal Factor can only be entered on the MINI-PAM-II screen.</p>	<p>Spec. Cal. Factor = 1.22</p>
Calibration of actinic light list	<p>Go to "Actinic Light List" on MINI-PAM-II display (Section 4.2.11, page 61) or Settings Window of WinControl-3 (Section 4.3.2.1, page 74), execute "Calibrate".</p>	

3.4.15 MQS-B/A Adapter Set for Thin Fiberoptics

The MQS-B/A Adapter Set is designed to measure the light intensity of the MINI-PAM/F1 fiberoptics using an MQS-B light sensor together with a data logger having a high-impedance BNC input, for example, the Walz ULM-500 Universal Light Meter & Data Logger.

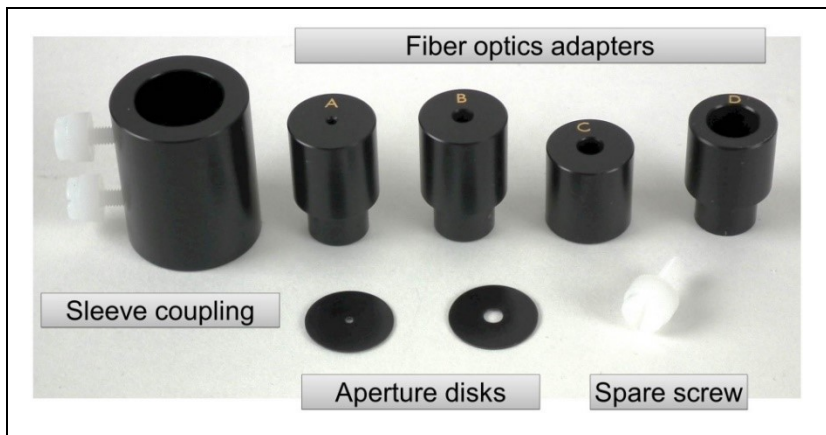


Fig. 27: MQS-B/A Adapter Set

4 Touchscreen Operation

Touchscreen operation of the MINI-PAM-II allows saturation pulse analysis of photosynthesis independent from a computer. Using the touchscreen, single F_V/F_M or $Y(II)$ measurements as well as more complex experiments like induction and light curves are feasible. Continuous recording of steady state fluorescence, however, requires operation of the MINI-PAM-II by WinControl-3 running on an external computer.

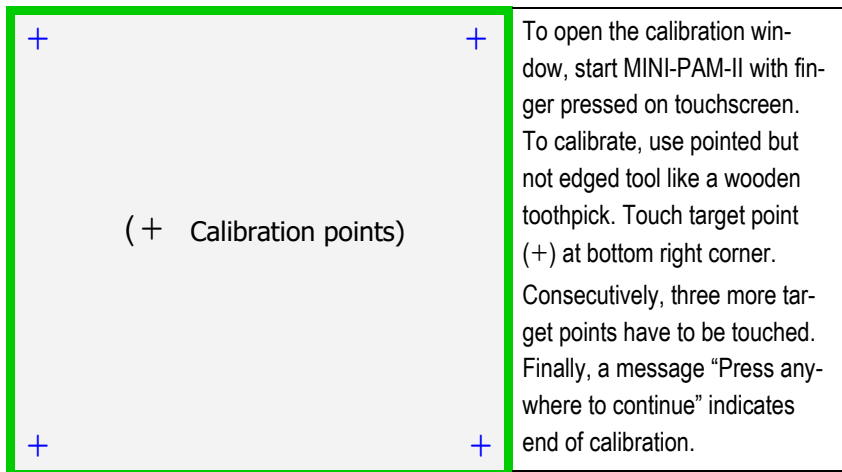


Fig. 28: Touch Screen Calibration

4.1 Calibration

The MINI-PAM-II can be operated using its resistive touchscreen. The touchscreen is factory calibrated. If touch operation is imprecise, calibration can be performed as described in Fig. 28.

4.2 Top-level Windows

The MINI-PAM-II provides various windows for control and data display. Frequently used commands, fluorescence data and the intensity list of actinic light (PAR list) are presented in 11 “top-level” windows (Table 4, below).

Table 4: Overview of Top-level Windows

	Window	Content
1	Basic Data	Minimum data set and basic action keys for filed operation.
2	Primary Data	Data of last saturation pulse analysis and current levels of fluorescence and additional measured parameters.
3	Porometer	Measurements, porometer status and commands (only when a MINI-PAM-II/POROMETER is connected).
4	Quenching Analysis	Fluorescence levels of last saturation pulse analysis and corresponding fluorescence ratios ($Y(II)$, F_v/F_m , ...).
5	Ft-Chart	Fluorescence signal over 25 or 125 s (see Fig. 71, page 98).
6	Spectrometer	Control and display of light, fluorescence and reflectance spectra (only when a MINI-SPEC/MP is connected).
7	Actinic + Yield	Short illumination program with saturation pulse analysis.
8	Induction Curve	Graphic of saturation pulse analyses of the last induction curve experiment.
9	Light Curve	Graphic of saturation pulse analyses of the last light response curve.
10	Recovery Curve	Graphic of saturation pulse analyses of the last dark phase following either an induction or a light response.
11	Actinic Light List	Set values of photosynthetically active radiation PAR in $\mu\text{mol m}^{-2} \text{s}^{-1}$.

Accessible from all top-level windows is the Main Menu and its submenus which allow adjusting settings of the MINI-PAM-II and its peripherals (Fig. 42, page 66; Fig. 43, page 67).

Main Panel			Side Panel	Side Panel
Window Title	100	A	SAT	The panel provides a button for saturation pulses (SAT) which is in some windows replaced by a start/stop button for automated experiments or by a CAL button to start PAR calibration. Navigation keys, a "MENU" key and action keys are located on the side panel.
Data, Graph			START	
			CAL	
			Navigation	
Action Keys			MENU	
			Act.L.	
Information Line			START	
			MEM	

Main Panel

From top to bottom: Window title and currently selected character to mark measurements, numerical and/or graphical data, action keys and information line. Action keys are absent or differently arranged in some windows. The information line displays date + time, or PAM activity, or informs on critical fluorometer states. Alphanumeric data are displayed in the "Cursor Mode" graphics windows

Fig. 29: Principal Screen Layout of Top-level Windows

Generally, all top-level windows consist of the "main panel" and the "side panel" (Fig. 29, page 45). The top line of the main panel displays window title, the battery charge percentage, and a capital letter which is added as a mark to each saturation pulse analysis. When the MINI-PAM-II/POROMETER is connected, the GPS status is indicated (Table 5).

Table 5: GPS-Status



GPS status. Left, ON. Right, ON and location determined.
(Only when the MINI-PAM-II/POROMETER is connected.)

The bottom of the main panel provides various action keys. Depending on the window, these action keys trigger saturation pulse analysis, control light conditions or affect graphic display. The “Information Line” at the bottom edge of the main panel shows current date and time. The Information Line can also include alphanumeric data when graphics windows are in the “Cursor Mode” which is started by touching the graph area.

The side panel provides arrow keys to change windows, control keys for fluorometer functions, and a **MENU** key to access the main menu. Control keys are different between windows.

An example is the top button on the sidebar. It triggers saturation pulses on the first three top-level windows (cf. Table 4, page 44), but the same button starts an experimental routine in the windows Actinic + Yield, Induction Curve, Light Curve, and Recovery Curve. Sections 4.2.1 to 4.2.11 will introduce all top-level windows in detail, except the Porometer window which is introduced in a separate manual:

https://www.walz.com/products/chl_p700/mini-pam-ii/downloads.html

4.2.1 Basic Data

The Basic Data window (Fig. 30, page 47) displays a reduced set of four data for fast sampling under field conditions. From these data, the Ft and PAR represent current measurements but the Y(II), or F_v/F_m , and the ETR are derived from the last saturation pulse analysis.

The bottom of the window provides keys for frequently used commands: **Rec** starts a new data set (Record), **Fo,Fm** determines maximum PS II photochemical yield, **Clock** starts repetitive triggering of saturation pulses, or an automated experimental routine,

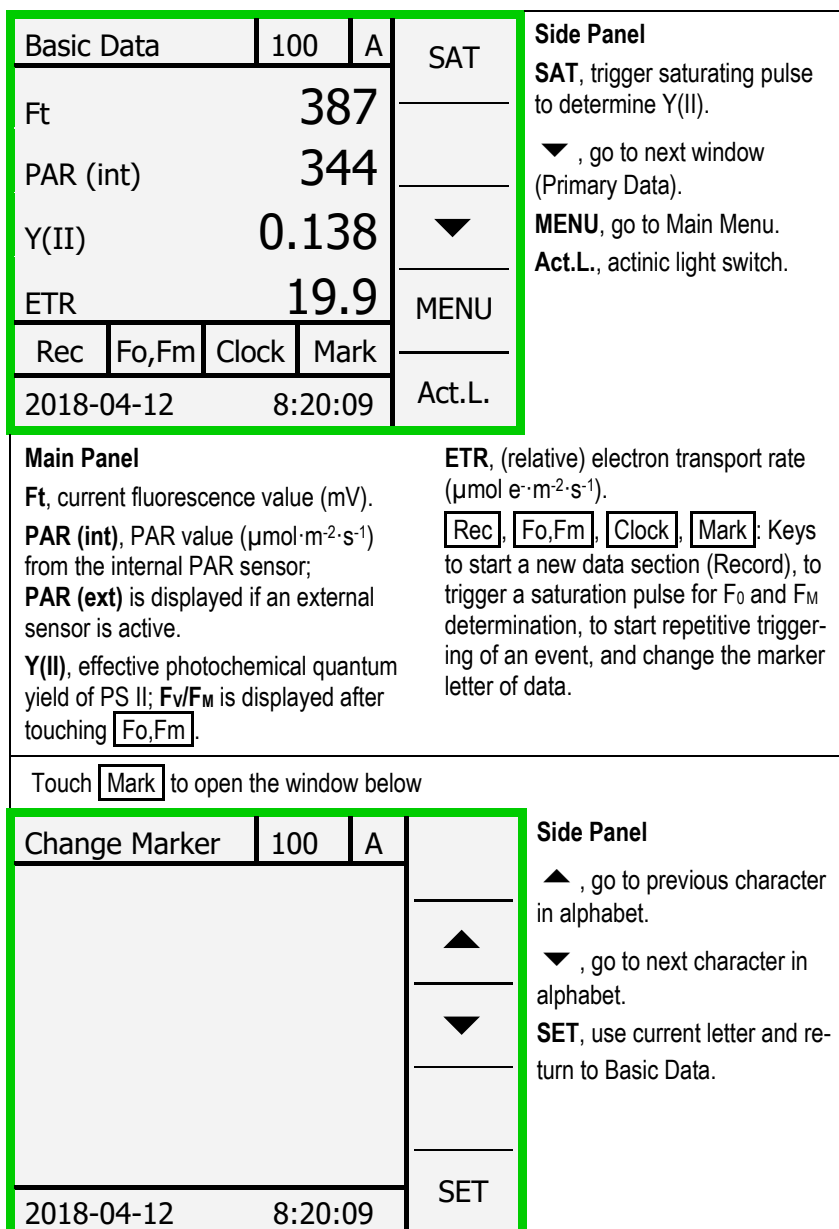


Fig. 30: Basic Data, Change Marker

and **Mark** opens the window “New Marker”. A “Mark” is a letter saved with each saturation pulse analysis.

To select a marker letter, use up and down keys in the window “Change Marker”. The current marker is shown on the top edge of the window. Touching the **SET** key confirms the current selection and returns to the Basic Data window.

The side panel of the Basic Data window offers four commands: the **SAT** key determines effective PS II photochemical yield, and the **Act.L.** represents a switch for actinic light. The **MENU** key leads to the Main Menu window as in all other top-level windows. The downward arrow key switches to the next window. Two arrow keys (up and down) are present in all other top-level windows except the last (Actinic Light List; Table 4, page 44) which requires only the upward key.

Comment on **Rec (New Record)**

Starting a new Record disconnects current saturation pulse analyses from that of the previous Record. Hence, without F_0 and F_M determinations after **Rec**, a **SAT** command will only result in data of $Y(II)$, q_P , and q_L , because these parameters can be calculated without F_0 and F_M (see Table 29, page 174).

However, the calculation of q_P and q_L requires a F_0' determination. F_0' can be measured by the routine “ F_0' mode”. If the “ F_0' mode” is inactive, F_0' can be calculated according to Oxborough and Baker (see Chapter 8, page 165). These calculations, however, require F_0 and F_M data. Note that all calculated F_0' values are preceded by a swung dash (tilde).

If a Record holds more than one F_0 , F_M determination, the latest F_0 , F_M will be used to calculate fluorescence coefficients. All calculations using both **Fo,Fm** and **SAT** data are only valid if the

two measurements were made with the same sample and under the same measuring conditions. Violating these conditions sometimes results in fluorescence ratio parameters exceeding their valid range (compare Table 29, page 174). Invalid data will be displayed in grey on the touch screen.

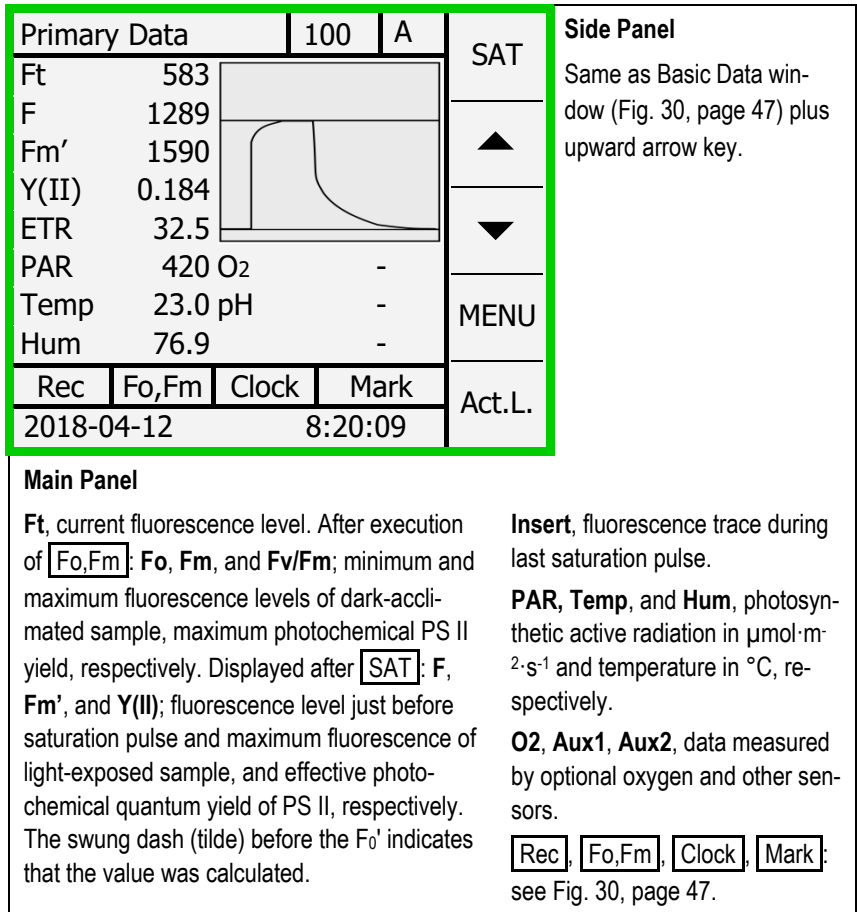


Fig. 31: Primary Data

4.2.2 Primary Data

The insert in the Primary Data window displays the fluorescence trace of the last saturation pulse. In addition, the window repeats the data of the previous one (F_t , F_v/F_m or $Y(II)$, ETR, PAR). New information of the Primary Data window is the F_0 and F_m signal levels (after **Fo,Fm** was pressed) or the F and F_m' signal levels (after **SAT** was pressed). Additionally, temperature, °C (Temp) and relative humidity, % (Hum), as measured by the MINI-PAM-II leaf clip (2035-B), are displayed.

4.2.3 Porometer

See separate manual.

4.2.4 Quenching Analysis

The window “Quenching Analysis” provides a complete overview on fluorescence levels and the fluorescence ratio quotients calculated by the MINI-PAM-II. Data line 2 to data line 4, compare data of the light exposed sample (left) with data of the dark-acclimated sample (right).

Quenching Analysis				100	A	SAT	Side Panel Same as Basic Data window (Fig. 30, page 47) plus upward arrow key.
F	919					▲	
Fm'	1250	Fm	2965				
Fo'	~454	Fo	575				
Y(II)	0.265	Fv/Fm	0.806			▼	
Y(NPQ)	0.425	qP	0.416				
Y(NO)	0.310	qL	0.206				
NPQ	1.372	qN	0.667		MENU		
Rec	Fo,Fm	Clock	Mark			Act.L.	
2018-04-12				8:20:09			

Main Panel

Numerical data: see Fig. 97, page 167 and Table 29, page 174.

Rec, Fo,Fm, Clock, Mark: see Fig. 30, page 47.

Fig. 32: Quenching Analysis

4.2.5 Ft-Chart

The Ft-Chart displays a 25 or 125 s interval of Ft where the right-most level of the graph corresponds to the current Ft value. The X axis interval can be adjusted in the menu “MINI-PAM-II Settings (Fig. 71, page 98).

Continuous Ft values are not saved when the MINI-PAM-II is operated in the stand-alone mode. Recording of Ft requires operation of the MINI-PAM-II by WinControl-3.

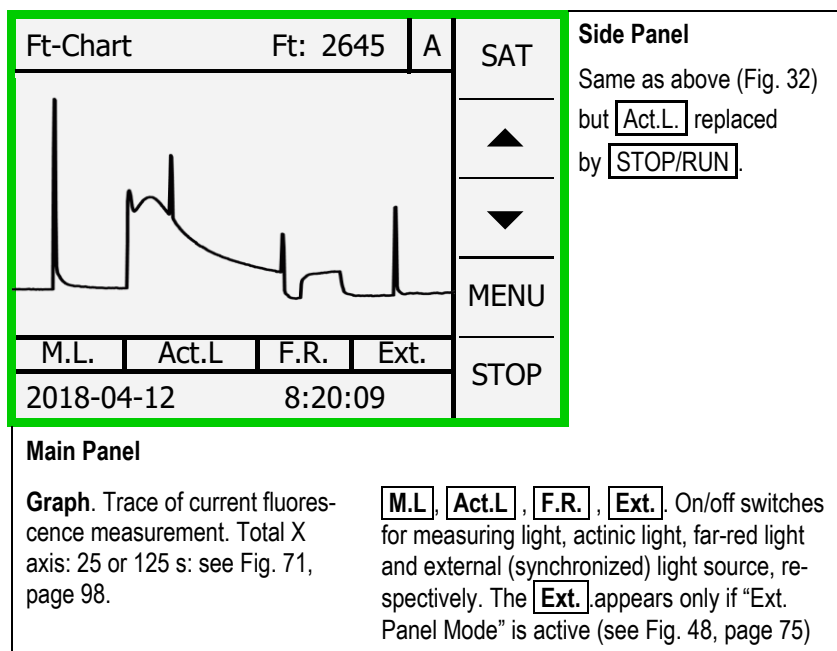



Fig. 33: Ft Chart

4.2.6 Spectrometer

The **SPEC** command of the Spectrometer window triggers a spectrometric measurement. In the standard mode, the spectrometer is equipped with a tube containing the optical entrance for external radiation (Fig. 10, page 23).

Spectra of photon flux densities are measured in units of $\text{nmol photons m}^{-2} \text{s}^{-1} \text{nm}^{-1}$. PAR corresponds to the integral of the spectrum from 400 to 700 nm. PAR is derived from the spectrometer when "Ext. PAR Sensor" is set to "On" in menu "Sensor Settings" (available by selecting "Sensors" in the Main Menu), and the spectrometer is selected as external PAR sensor (Main Menu →

Sensors → External PAR sensor → Spectrometer; Fig. 56, page 85).

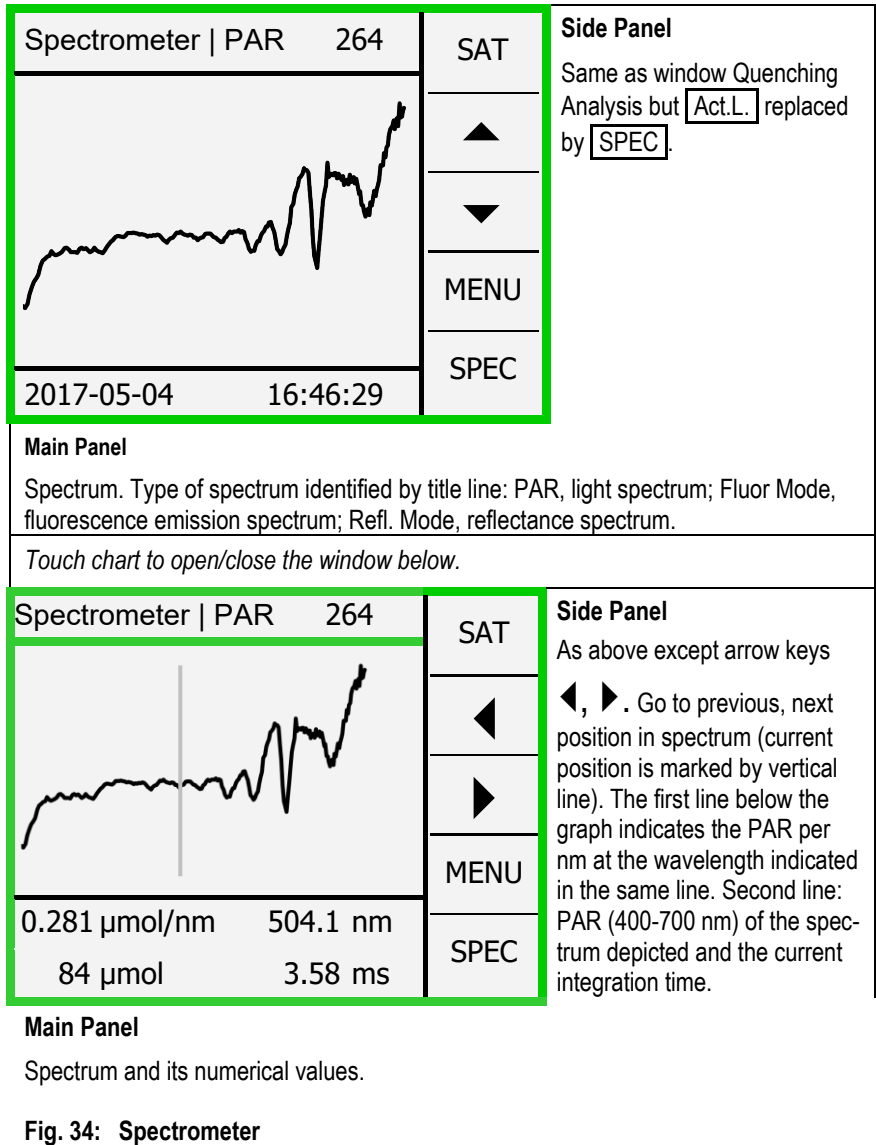

Note that the spectrometer measures the PAR of the MINI-PAM-II internal light correctly only in the CAL mode (Fig. 39, page 61) because only in the CAL mode the LEDs are operated continuously. In the measuring mode, actinic light is pulse-width modulated and the PAR value is fluctuating displayed.

To record fluorescence or reflectance spectra, change spectrometer configuration as described in Fig. 10 (page 23). Then select operation mode: Main Menu → Sensor → Spectrometer → Operation Mode.

In menu Operation Mode, the items “Fluores. (blue)”, “Fluores. (green)” and “Reflectance” set the spectrometer mode to fluorescence spectra with blue excitation, fluorescence spectra with green excitation and reflectance spectra, respectively.

For reflectance measurements, first measure dark current in complete darkness, then the 100% reflectance signal with the white reference material (Fig. 11, page 25), and finally the sample.

In the Spectrometer window, touching the graph display area show a vertical cursor line as well as the x-y data of the intersection between cursor line and spectrum. Use the arrow keys to navigate through the spectra. Touching the graph display again returns to the original function of arrow keys.

4.2.7 Actinic + Yield

The “Actinic + Yield” window is the first of four windows for automated measuring routines. The parameters of the Actinic + Yield routine can be set in the menu “Actinic + Yield Settings” (Fig. 51, page 80) which is accessible via “Program/Clock” of the Main Menu.

The routine illuminates a sample with actinic light of a defined period. Depending on settings, saturation pulse analysis is carried out prior and after actinic light exposure, or only after actinic light exposure. Because the Actinic + Yield experiment is rather short, it is frequently employed when response to light exposure needs to be evaluated for many samples.


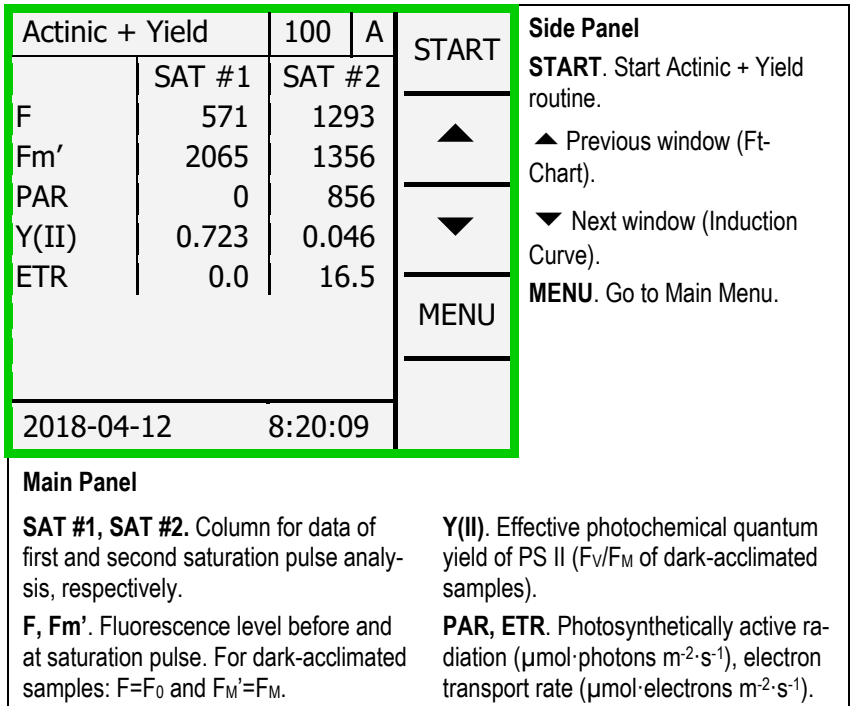


Fig. 35: Actinic + Yield

4.2.8 Induction Curve

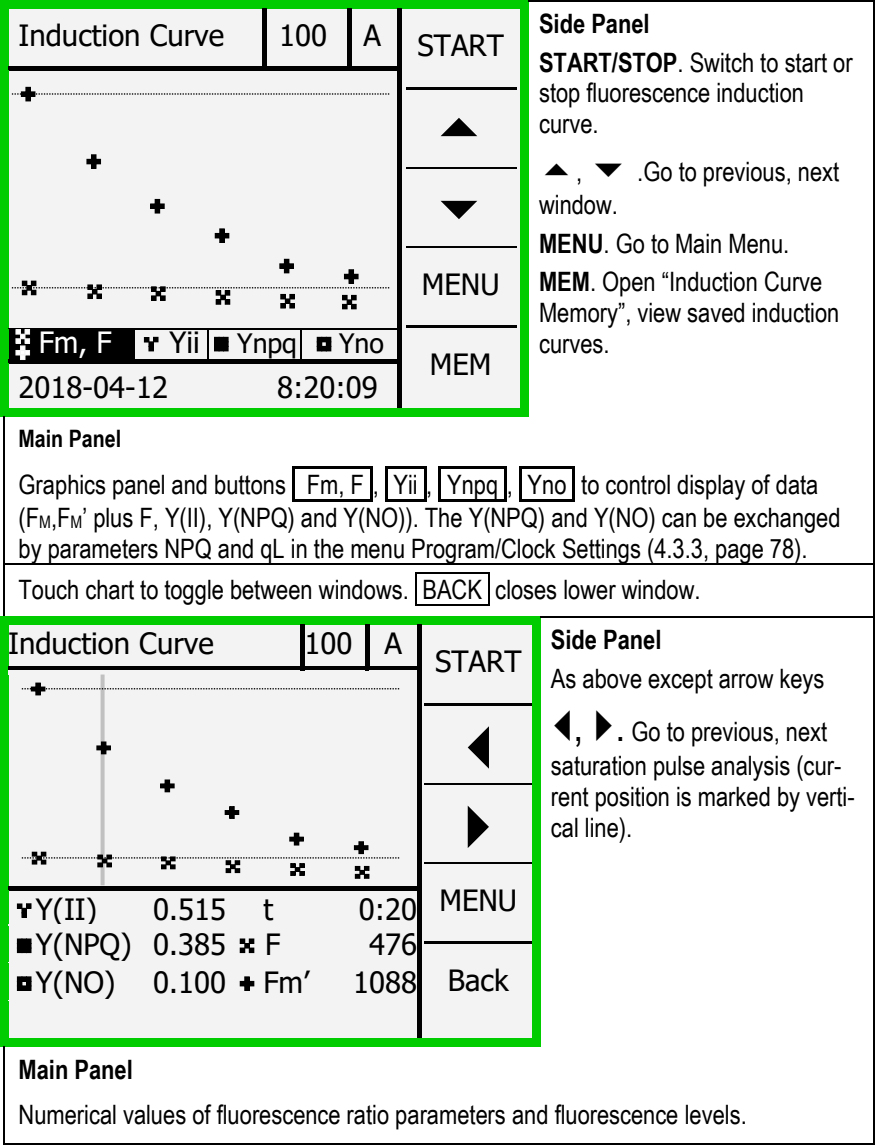

The “Induction Curve” window controls fluorescence induction curve experiments (Fig. 36, page 56). The graphics panel provides a qualitative picture of induction curve data.

The window also provides numerical data of fluorescence ratio parameters and fluorescence levels. These numerical data appear by touching the graphics panel. The navigation keys move the cursor (grey vertical line) from one saturation pulse analysis to another. The cursor indicates the position within the induction curve of the currently displayed set of numerical data.

The parameters of an induction curve experiment, number of saturation pulse analyses and interval between them, can be adjusted in the menu “Induction Curve Settings” (Fig. 52, page 81) which is available over the Program/Clock line of the Main Menu (Fig. 42, page 66). In the “Induction Curve Settings” menu, you can choose to continue fluorescence monitoring after the induction curve in the dark (recovery curve). The **MEM** of the Induction Curve window opens the “Induction Curve Memory” window. In this window, the navigation keys scroll through stored induction curves.

4.2.9 Light Curve

In analogy to the previous window (Induction Curve), the “Light Curve” window provides buttons to start and stop light curves and to survey them (Fig. 37, page 59). Also, the Light Curve window provides numerical data of saturation pulse analysis which can be accessed by touching the graphics panel. Navigation between different saturation pulses analyses and selection of displayed data works as described for the previous window (Induction Curve).

Light curve parameters (number and duration of light steps, initial PAR, and recovery curve) can be adjusted in the menu “Light Curve Settings” (Fig. 53, page 82). Touching the **MEM** button and scrolling using the arrow keys on the side bar permits viewing stored light curves.

4.2.10 Recovery

Often, fluorescence recovery experiments are automatically appended to an induction or a light curve (cf. 0 and 4.2.9). In these cases, the recovery curves are started automatically. Otherwise, recovery curves can be initiated (and cancelled) manually using the **START/STOP** button on the side panel of the Recovery window (see Fig. 38, page 60).

As in previous Induction and Light Curve windows, recovery data are represented graphically and numerically. The time course of recovery curves is fixed: each curve last 39 min during which 7 saturation analysis are carried out. In case of a preceding induction or light curve, the last saturation pulse analysis of the induction or light curve corresponds to the first one of recovery. The interval between neighboring saturation pulse analysis roughly doubles with time (Table 6, page 15).

In the Recovery window, the **MEM** button opens the window “Recovery Memory” in which all recovery kinetics can be viewed independent if they are connected to an induction or light curve or if they represent separate experiments.

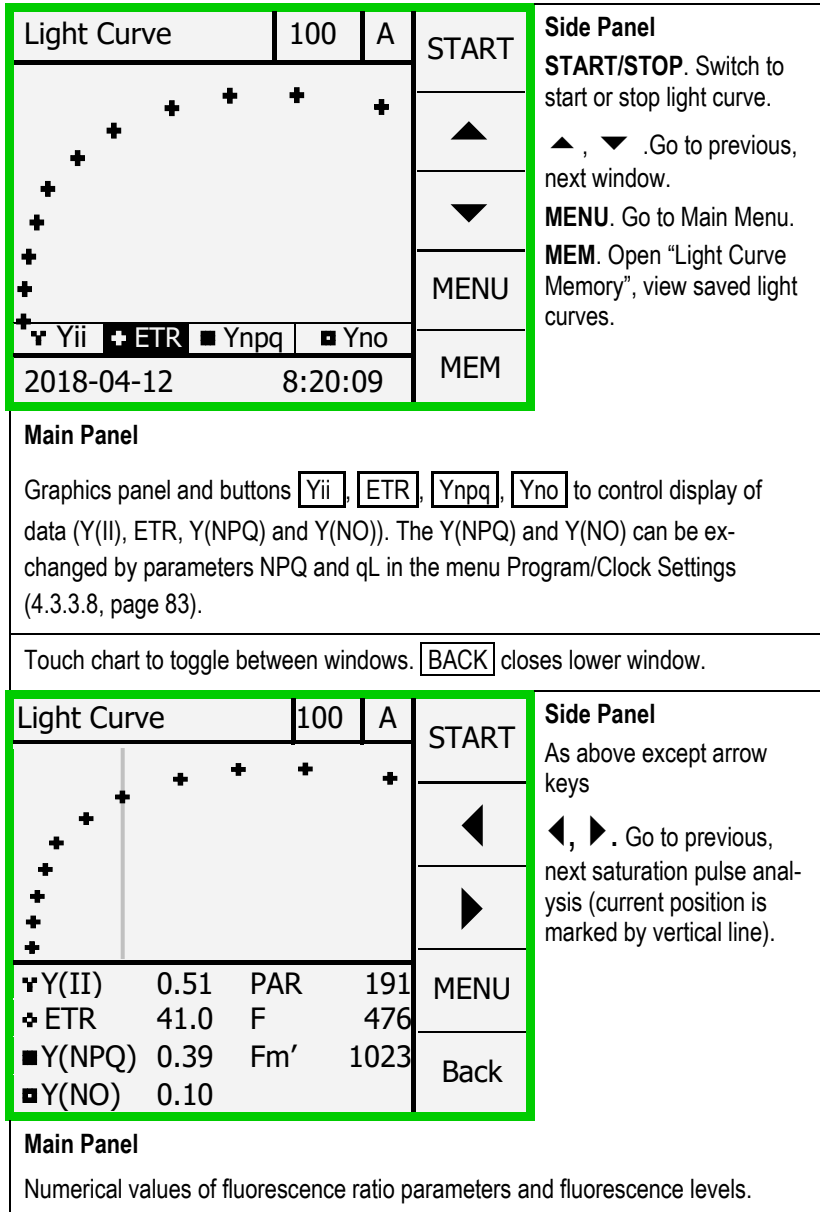


Fig. 37: Light Curve

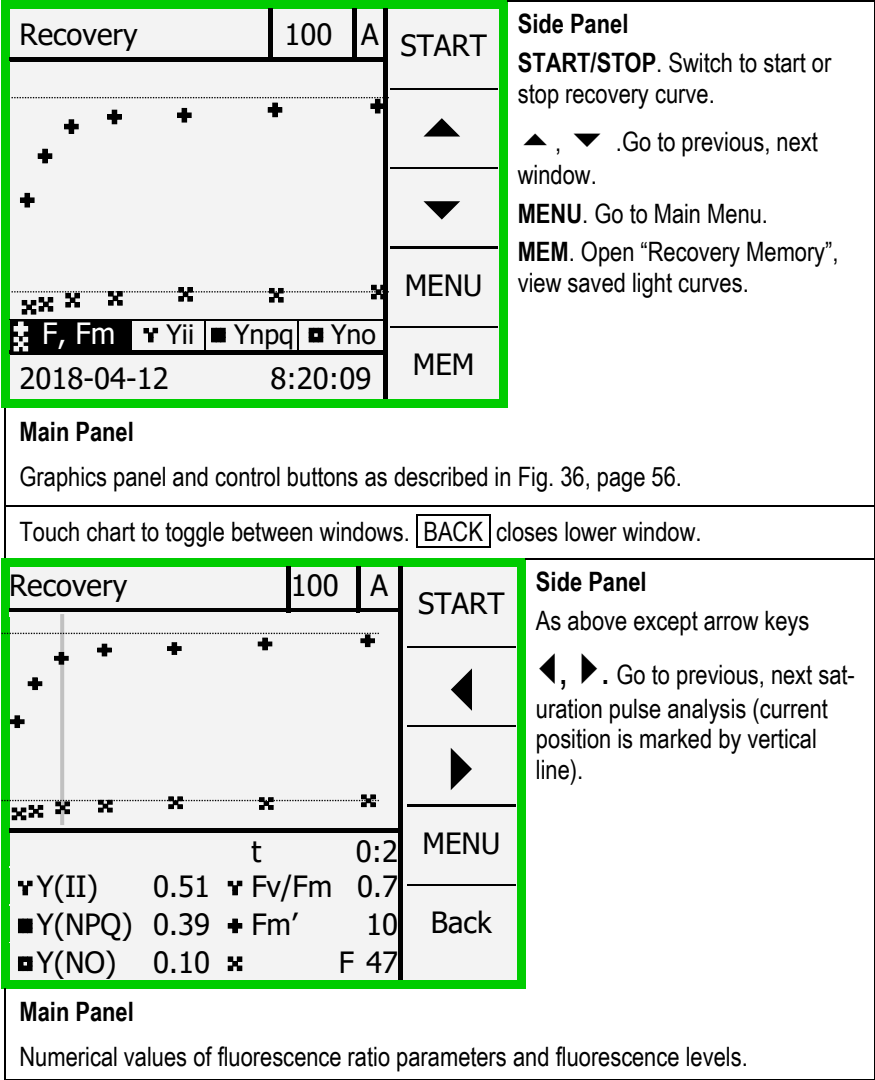


Table 6: Sequence of Saturation Pulse Analyses in a Recovery Curve

SAT number	Time in darkness, min
1	0:00
2	0:30
3	1:30
4	4:00
5	9:00
6	19:00
7	39:00

4.2.11 Actinic Light List

Actinic Light List				100	A	CAL	Side Panel CAL. Calibrate Actinic Light List.
1:	25	7:	285				
2:	45	8:	420			▲	▲ Go to previous window.
3:	65	9:	625				MENU , go to Main Menu.
4:	90	10:	820				
5:	125	11:	1150				
> 6:	190	12:	1500			MENU	
External/Internal Actinic Light							
2018-04-12 8:20:09							

Main Panel

PAR target values ($\mu\text{mol m}^{-2} \text{s}^{-1}$) of the 12 settings for actinic light intensity. The PAR target values are valid for an actinic light factor of 1.00 (see menu “Light Sources” Fig. 47, page 75). The “>” indicates the currently selected PAR level.

Fig. 39: Actinic Light List

The window “Actinic Light List” displays PAR values (in $\mu\text{mol m}^{-2} \text{s}^{-1}$) which increase roughly exponentially with intensity settings from 1 to 12. The entire PAR list can be multiplied by the internal actinic factor (“Int. Act. Factor, Fig. 47, page 75).

In the factory, the internal actinic light is adjusted so that the PAR values of the Actinic Light List apply to the sample level of the 2035-B leaf clip with the MINI-PAM-II fiberoptics fully inserted (i.e. distance between fiberoptics tip and sample level of 7 mm; angle between end piece of fiberoptics and sample level of: 60°). Because of the same spatial situation, these factory settings also apply to the 2060-B Arabidopsis Leaf Clip when distance rings 2 mm and 4 mm are used, and the 2010-A Distance Clip 60°.

External versus Internal Actinic Light

The main panel of the window Actinic Light List (Fig. 39, page 61) indicates the light source used for actinic illumination.

External Actinic Light is displayed when:

- An external lamp (2054-L External LED Source) is connected to the SYNC port of the MINI-PAM and “PAM Mode (Act)” is selected as operation mode (see 4.3.2.6, page 76).

Internal Actinic Light is displayed when:

- An external lamp (2054-L External LED Source) is connected to the SYNC port of the MINI-PAM and “Ext. Panel Mode” is selected as operation mode (see 4.3.2.6, page 76)..
- An external lamp is not connected.

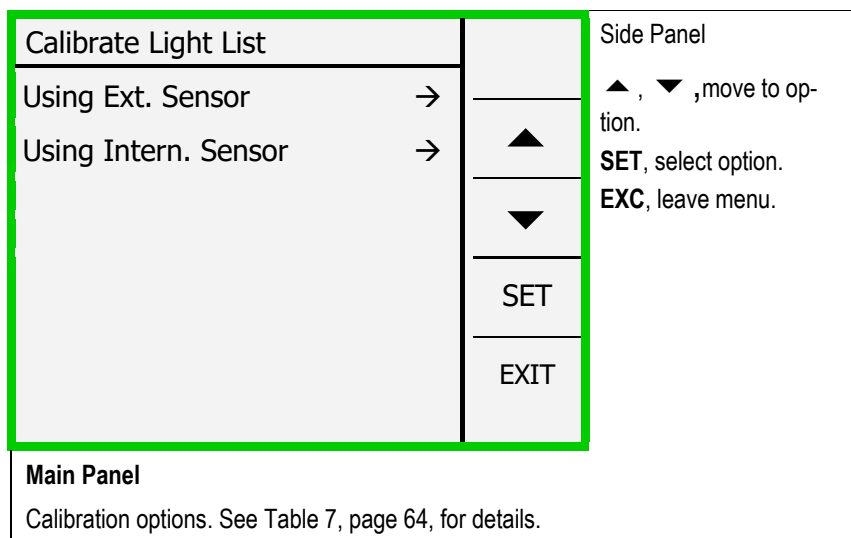


Fig. 40: Calibrate Light List

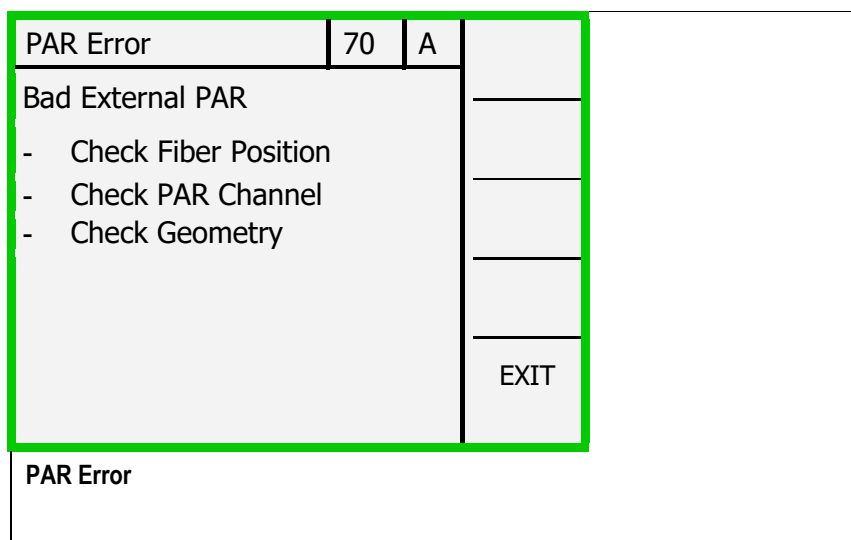


Fig. 41: PAR Error

The **CAL** command adjust internal or external actinic light so that the PAR at sample level corresponds to the “Actinic Light List”.

Clicking **CAL** opens the window “Calibrate Light List” in which an external or internal light sensor can be selected (Fig. 40). The option “Use External Light Sensor” requires that an external light sensor is connected. If “Use External Light Sensor” is attempted in the absence of an external sensor, an error message pops up (Fig. 41, page 63). The same error message appears when the external sensor is not exposed to the actinic light, or when the external sensor to be used was not set as active sensor (see Fig. 56, page 85). The option “Use Internal Light Sensor” cannot be used to calibrate an external light source. Both the external and internal light sensor options are only available for the internal actinic light source.

Table 7: The CAL Command

Light source being calibrated.	Ext. PAR Sensor (e.g., 2035-B Leaf-Clip Holder).	External Light Source (e.g., LED Light Source 2054-L).	Action.
Internal actinic light.	Connected	Absent OR Present and “Ext. Panel Mode” active.	If “Using Ext. Sensor” selected: Calibration of internal light sensor. Then, calibration of the Actinic Light List with the internal light sensor.
Internal actinic light.	Absent or connected	As above.	If “Using Intern. Sensor” selected: Calibration of the Actinic Light List with the internal light sensor.
External actinic light.	Connected	Present and “Ext. Panel Mode” active.	Calibration of the Actinic Light List with the external light sensor.*

Using the MINI-SPEC/MP Miniature Spectrometer as external PAR sensor: See Section 3.4.1, page 21.

Calibration of the internal light sensor: See Section 4.3.4.6, page 95.

Calibration of an external light source. If “External Actinic Illumination” is displayed by the window “Actinic Light List”, the **CAL** command adjusts the PAR emitted by the 2054-L External LED Source to the PAR values of the Actinic Light List. Typically, the external sensor is the PAR sensor of 2035-B leaf clip.

To separately adjust each of the four light types of the 2054-L lamp, **CAL** command switches the PAR sensor of the 2035-B leaf clip automatically to the “Mixed Mode” (see Section 4.3.4.5, page 91). After calibration, the measuring mode of the PAR sensor is returned to its previous mode. When checking the light intensities of the 2054-L lamp, make sure that the Mixed Mode is selected.

Internal Light Sensor. The internal sensor is exposed to a small and constant fraction of internal actinic light. The properly calibrated internal PAR sensor measures continuously internal light intensity. In the factory, the internal sensor has been adjusted to indicate the PAR values at sample level of the 2035-B leaf clip when the fiberoptics is fully inserted.

When the sample level differs from the standard spatial arrangement (2035-B leaf clip), the internal PAR sensor must be recalibrated. For this purpose, a calibrated PAR sensor is positioned at the selected sample level. Note that the PAR sensor of the 2035-B leaf clip is magnetically attached and can easily be disconnected to adjust the internal PAR sensor for various measuring configurations.

If a calibrated PAR sensor feeds its data in the MINI-PAM-II (as is the case with the PAR sensor of 2035-B leaf clip), the internal PAR sensor can be calibrated automatically. To do so, select “Calibrate” in window “Int. PAR Sensor Settings” and start automatic adjustment of the calibration factor (see Section 4.3.4.6, page 95).

If the calibrated PAR sensor is not compatible with the MINI-PAM-II, switch light on, and vary in the window “Int. PAR Sensor Settings” (Fig. 69, page 96) the “Calibration Factor” until internal PAR matches external PAR.

4.3 Main Menu

The “Main Menu” (Fig. 42, page 66) is the central hub to access settings, calibration data, hardware information and the memory of the MINI-PAM-II. The Main Menu consists of eight items. An item is selected by the arrow keys of the side panel followed by the **SET** command, or by directly touching a menu line. Most of these items lead to submenus that themselves link to lower-level menus. The complex architecture of the Main Menu and its submenus is outlined in Fig. 43, page 67.

Main Menu	A		Side Panel
PAM Settings	→		▲, ▼. Up/down keys move cursor.
Light Sources	→		SET . Open submenu of selected line.
Program/Clock	→	▲	EXIT . Return to previous (top-level) window.
Sensors	→		
MINI-PAM-II Settings	→	▼	
Memory	→		
Info	→	SET	
Reset	→		
		EXIT	

Fig. 42: Main Menu

PAM Settings	
Meas. Light	On/Off
Meas. Light Sett.	→
Gain	#
Damping	#
ETR-Factor	#
Fo' Mode	On/Off
Adjust F-Offset	→

Meas. Light Settings	
Meas. Light Int.	#
Meas. Light Freq.	#
ML-F High	On/Off

Light Sources	
Actinic Light	On/Off
Actinic Intensity	#
Actinic Factor	#
Far Red	On/Off
Far Red Sett.	→
Light Panel Sett.	→
SAT Settings	→

Far Red Settings	
Far Red Int.	#
SAT Red Width (s)	#

Light Panel Settings	
Intensity	#
Red	#
Green	#
Blue	#
White	#
Light Panel	#
Operation Mode	→

SAT Pulse Settings	
SAT Intensity	#
SAT Width (s)	#

Light Panel Oper. Mode	
PAM Mode (Act)	
Ext. Panel Mode	X

Program/Clock Sett.	
Clock	On/Off
Interval (mm:ss)	#
Clock Item	→
Actinic + Yield	→
Induction Curve	→
Light Curve	→
Yield + # Poro Only	→
Y(II), Y(NPQ), Y(NO)	X
Y(II), NPQ, qL	

Actinic+Yield Settings	
Width (mm:ss)	#
Initial Pulse	Yes/No

Ind. Curve Settings	
Delay (mm:ss)	#
Width (mm:ss)	#
Length	#
With FoFm Pulse	
Add Recovery	Yes/No

Light Curve Settings	
Width (mm:ss)	#
Initial Intensity	#
Length	#
With FoFm Pulse	On/Off
Add Recovery	Yes/No

Clock Item	
SAT Pulse	X
Actinic + Yield	
Light Curve	
Light Curve + R	
Induction Curve	
Ind. Curve + R	
Poro Only	
Yield + Poro Only	

Fig. 44: First Three Submenus in Detail

4.3.1 PAM Settings

The menu item “PAM Settings” includes adjustments of the way how the MINI-PAM-II acquires PAM fluorescence. The PAM Settings menu contains seven items (Fig. 45, page 69) but also displays the current Ft value and the currently active offset, which is automatically subtracted from the raw signal to obtain Ft.

PAM Settings		A	OFF	Side Panel ON/OFF. Switch for measuring light. ▲ , ▼ . Up/down keys move cursor. SET. Open submenu (→) or select parameter, change parameter value by up/down keys. EXIT. Return to Main Menu.
Meas. Light	On/Off		▲	
Meas. Light Sett.	→			
Gain	1		▼	
Damping	2			
ETR-Factor	0.84			
Fo' Mode	On/Off		SET	
Adjust F-Offset	→			
Ft	421		EXIT	
Current F-Offset	70			

Fig. 45: PAM Settings

4.3.1.1 Meas. Light

On/Off switch for measuring light (weak excitation light consisting of μs pulses). Measuring light can be switched by selecting “Meas. Light” and touching **SET**. The **OFF/ON** key turns measuring light off or on, independent of the menu item selected.

4.3.1.2 Meas. Light Sett.

The item “Measuring Light Settings” opens a menu for configuring the measuring light. The variables are measuring light intensity (Meas. Light Int.), measuring light frequency (Meas. Light. Freq.), and the option to set measuring light frequency to high (MF-F High). To adjust, select variable by the Up/Down arrows and press **SET**. If several options are available, select desired setting by Up/Down arrows followed by **SET**.

Measuring Light Settings	A		Side Panel
Meas. Light Int.	6	▲	ON/OFF. Switch for measuring light.
Meas. Light Freq.	3		▲ , ▼ . Up/down
ML-F High	On/Off	▼	keys move cursor.
		SET	SET. Open submenu (→) or select parameter, change parameter value by up/down keys.
		EXIT	EXIT. Return to Main Menu.

Fig. 46: Measuring Light Settings

Measuring light intensity: At constant frequency, measuring light intensity can be considered as proportional to settings 1 to 12. The fluorescence signal increases with measuring light intensity. High measuring light intensities can interfere with the correct F_0 fluorescence measurements.

The average PAR of measuring light at highest frequency and highest intensity setting was measured to be $1.5 \mu\text{mol m}^{-2} \text{s}^{-1}$ by

the PAR sensor of the 2035-B leaf clip and MINI-PAM-II fiber in the fully inserted position. For the same geometrical arrangement, average measuring light intensities can be estimated using the subsequent equation:

$$I_{ML} \left(\frac{\mu\text{mol}}{\text{m}^2 \cdot \text{s}} \right) = 1.5 \left(\frac{\mu\text{mol}}{\text{m}^2 \cdot \text{s}} \right) \cdot \frac{f(\text{Hz})}{100(\text{Hz})} \cdot \frac{\text{Int. Sett.}}{12}$$

where I_{ML} , f and *Int. Sett.* is the current measuring light intensity in $\mu\text{mol m}^{-2} \text{s}^{-1}$, the current measuring light frequency, and the current intensity setting for measuring light, respectively.

Measuring light frequency: Five frequency levels are available (see Table 8, page 72). Increasing the measuring light frequency improves the signal quality but can also be actinic and therefore interfere with F_0 fluorescence measurements. Measuring light frequency does not affect the frequency of acquisition of F_t by the WinControl-3 software.

High measuring light frequency: The highest measuring light frequency is 100 Hz. The “ML-F High” command overrides settings made under “Meas. Light. Freq.” High measuring light frequency improves signal quality but bears the risk that its higher intensity drives photosynthesis, that is, the measuring light becomes actinic. In this case, the F_0 may be overestimated. Measuring light frequency is automatically switched to “high” for saturation pulse analysis.

4.3.1.3 Gain

Selecting “Gain” by the SET gives access to four electronic amplification factors (1 to 4) which can be adjusted by the arrow keys.

Table 8: Measuring Light Frequencies

Setting	Frequency, Hz
1	5
2	10
3	15
4	20
5	25
high	100

4.3.1.4 Damping

Damping is a software-based filter that specifically suppresses high frequency noise and, thus, can improve signal quality. Changing damping settings uses the same principle as described for “Gain”. Default setting for damping is 2 (two). Changing damping to higher values can make the MINI-PAM-II response slow.

4.3.1.5 ETR-Factor

This factor is used for ETR calculations and corresponds to the fraction of incident PAR absorbed by a leaf; its default value is 0.84 (cf. Section 8.3, page 172).

4.3.1.6 Fo’ Mode

After saturation pulses, the “Fo’ Mode” replaces actinic light by far-red light to quickly open PS II reaction centers (cf. Chapter 8, page 165). The measured Fo’ is the minimum Ft during this far-red period. Interval and intensity of far-red illumination can be adjusted in menu “Light Sources” (Fig. 47, page 75).

4.3.1.7 Adjust F-Offset

The “Adjust F-Offset” command determines the background signal for subtraction from the total signal. Background signals must possess the modulation characteristics of measuring light to be recognized by a PAM fluorometer. These signals can arise from:

- Fluorescence from suspension media or detector filter excited by measuring light.
- Traces of modulated excitation light transmitted by the detector filter.
- Non-optical modulated "electronic noise".

Usually, the background signal increases with measuring light intensity and signal amplification (gain). Therefore, the Adjust F-Offset command determines the background signal for all measuring light intensities and all gain settings. The currently active offset is displayed in the bottom line of the PAM Settings window (Fig. 45, page 69).

Procedure

- Choose dim environment.
- Switch off any flickering light sources like fluorescent lamps or computer screens.
- Point fiber tip away from any objects, keep fiber tip clear.
- Run “Adjust F-Offset”

4.3.2 Light Sources

4.3.2.1 Actinic Light

On/off switch for actinic illumination. Instead via the menu item, actinic light can be switched on and off by the **OFF/ON** key on the side panel of the window (Fig. 47, page 75).

4.3.2.2 Actinic Intensity

Intensity regulation for actinic light. Select menu item by **SET** and choose setting using arrow keys. Settings 1 to 12 are available. PAR information of settings is shown in window “Actinic Light List” (Fig. 39, line 61).

4.3.2.3 Actinic Factor

Factor multiplying target PAR values in window “Actinic Light List”. Factor range is 0.5 to 2.0. Maximum intensity might be constraint by LED limits.

4.3.2.4 Far-red

On/off switch for far-red light.

Light Sources			Side Panel ON/OFF , actinic light switch. ▲ , ▼ . Up/down keys move cursor. SET . Open submenu (→) or select setting. To change parameter value, use up/down keys. EXIT . Return to Main Menu.
Actinic Light	on	OFF	
Actinic Intensity	6		
Actinic Factor	1.00	▲	
Far Red	off		
Far Red Sett.	→	▼	
Light Panel Sett.	→		
SAT Settings	→	SET	
		EXIT	

Fig. 47: Light Sources

Light Panel Settings			Side Panel ▲ , ▼ . Up/Down keys to move to a menu line. SET . Select setting, up/down keys change setting value. EXIT . Return to Light Sources menu.
Intensity	-		
Red	0		
Green	1	▲	
Blue	0		
White	0	▼	
Light Panel Operation Mode	→	SET	
		EXIT	

Fig. 48: Light Panel Settings

4.3.2.5 Far-red Sett.

Opens the menu to adjust intensity and duration of far-red illumination. These settings are active for F_0' determinations (Section 8.2, page 168).

4.3.2.6 Light Panel Sett.

Use **SET** to enter submenu “Light Panel Settings” (Fig. 48, page 75). The function of this window depends on the “operation mode”. The operation mode is selected in the last line in this window.

Two operation modes are available: “PAM Mode (Act.)” and “Ext. Panel Mode”. In the first mode (PAM Mode (Act.)), actinic light is provided exclusively by an external source (e.g., the 2054-L External LED Source, Fig. 14, page 30). In the second mode (Ext. Panel Mode), the MINI-PAM-II internal light is used as actinic light to which an external light source may be added (see Table 23, page 147).

Note: Operation of the external light source (2054-L External LED Source) requires line power.

Note: In absence of an external light source, “PAM Mode Act.” is not available.

For **PAM Mode (Act.)**, proper calibration of the external light source is mandatory. In this mode, the item “Intensity” and the Light Panel On/Off switch in Fig. 48 are not available.

In the “PAM Mode (Act.)” mode, the total intensity of all four LED colors is adjusted as described above for internal actinic light, that is, 12 intensity settings are available with the intensities defined in

the “Actinic Light List” (Fig. 39, page 61), and the intensities can be modified by the Actinic Factor (Section 4.3.2.3, page 74).

In the Light Panel Settings window (Fig. 48, page 75), the numbers entered for the four LED colors (red to white) determine the fraction with which each LED color contributes to the actinic light. For example, if red=1, green=1, blue=3 and white=0, the actinic light contains 20% each of red and green, 60% of blue and no white.

In the **“Ext. Panel Mode”**, in the window Light Panel Settings (Fig. 48), the item “Intensity” and the Light Panel On/Off switch are available. The intensity scales range from 0% to 100% (maximum intensity). The typical maximum value (100%) for the individual LEDs is about $1500 \mu\text{mol m}^{-2} \text{s}^{-1}$. When all four LED intensities are set to 100%, the maximum total intensity is $> 5500 \mu\text{mol m}^{-2} \text{s}^{-1}$. Exact PAR values of the external panel are not available in the Ext. Panel Mode.

4.3.2.7 SAT Settings

Moving to “SAT Settings”, and then **SET** opens the submenu “SAT Pulse Settings” (see below) in which relative intensity (1 to 12) and duration (width, 0.2 to 2.0 s) of saturation pulses can be set.

At sample level, at intensity setting 12 the SAT intensity corresponds to $6000 \mu\text{mol m}^{-2} \text{s}^{-1}$ under the geometrical conditions of a 2035-B leaf clip (distance fiberoptics tip to sample level: 7 mm; angle between end piece of fiberoptics and sample level: 60°). The saturation pulse intensity can be adjusted at increments of $500 \mu\text{mol m}^{-2} \text{s}^{-1}$.

4.3.3 Program/Clock

The menu “Program/Clock Settings” provides all options to configure automated experimental routines (Actinic + Yield, Induction Curve, and Light Curve) well as repetitive triggering of single measurements and experimental routines.

Program/Clock Settings				Side Panel
Clock	On/Off			▲ , ▼ . Up/Down keys to move to a menu line. SET . Open submenu (→) or adjust setting using the Up/Down keys. EXIT . Return to Main Menu.
Interval (mm:ss)	→		▲	
Clock Item	→			
Actinic + Yield	→		▼	
Induction Curve	→			
Light Curve	→		SET	
Yield + # Poro	→			
Y(II), Y(NPQ), Y(NO)	X		EXIT	
Y(II), NPQ, qL				

Fig. 49: Program/Clock Settings

4.3.3.1 Clock

On/off switch of clock. The “Clock” triggers repetitively an event at a defined interval. The interval is specified in “Clock Interval” and the event in “Clock Item”.

4.3.3.2 Clock Interval

Adjust clock interval between 10 s and 60 min by selecting “Interval” (up/down keys and **SET**, respectively) and adjusting time interval (up/down keys).

4.3.3.3 Clock Item

Saturation pulse analyses, and the programs Actinic + Yield, Induction Curve and Light Curve can be repetitively performed under clock control. Recovery experiments can be appended to induction and light curves (item Light Curve + R and Induction Curve + R, respectively). Porometer measurements can be triggered without (Poro Only) or in combination with saturation pulse analysis (Yield + Poro Only). To select one of the eight items in menu “Clock Item” (Fig. 50), move cursor to the item of interest and touch **SET**. The selected item is then marked by an X.

Clock Item		Side Panel
SAT Pulse	X	▲ , ▼ . Up/Down keys to move to a menu line.
Actinic + Yield	▲	SET . Open submenu (→) or adjust setting using the Up/Down keys.
Light Curve	▼	EXIT . Return to Main Menu.
Light Curve + R	SET	
Induction Curve		
Induction Curve + R		
Poro Only		
Yield + Poro Only	EXIT	

Fig. 50: Clock Item

4.3.3.4 Actinic + Yield

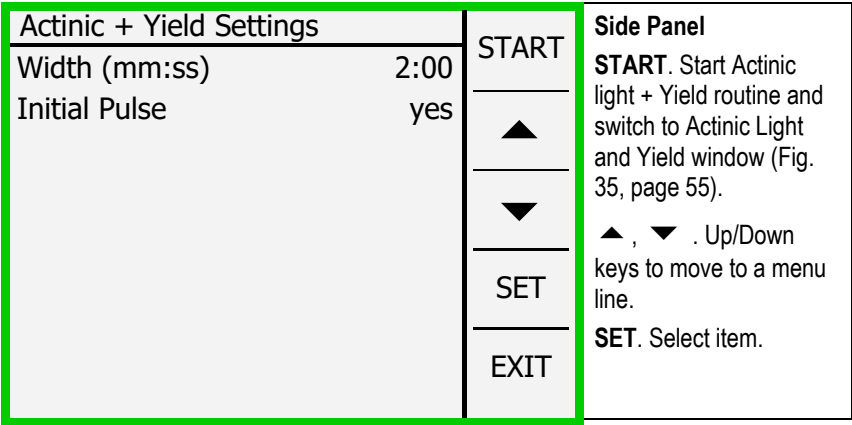


Fig. 51: Actinic + Yield Settings

The behavior of the Actinic + Yield program is defined by two factors (Fig. 51, page 80): the duration (width) of actinic illumination (possible settings from 5 s to 5 min) and the option to start actinic illumination without preceding saturation pulse analysis (Initial pulse). The width of actinic illumination is adjusted as described above for clock interval and initial pulse is selected by the **SET** command. Actinic light intensity is adjusted in the window “Light Sources” (Fig. 47, page 75).

An Actinic + Yield routine can be started using the **START** button on the side panel of Fig. 51. In this case, the screen display will automatically switch to the Actinic + Yield experimental window (Fig. 35, page 55).

4.3.3.5 Induction Curve



Induct. Curve Settings		START   SET EXIT	Side Panel START. Start induction curve experiment and switch to Induction Curve window (Fig. 36, page 56). ▲ , ▼ . Up/down keys move cursor. SET. Selects item which then can be changed by arrow keys.
Delay (mm:ss)	1:00		
Width (mm:ss)	0:30		
Length	10		
With FoFm pulse	on		
Add Recovery	no		

Fig. 52: Induction Curve Settings

Induction curve experiments are configured in the window “Induction Curve Settings” (Fig. 52, page 81).

Delay (range 5 s to 10 min) defines the dark interval between saturation pulse analysis with the dark-acclimated sample (F_0 , F_M determinations, Chapter 8, page 165) and beginning of actinic illumination.

Width (range 5 s to 10 min) is the time interval between two successive saturation pulse analyses during illumination.

Length is the number of saturation pulse analyses carried out during actinic illumination. Thus, the duration of actinic illumination is “Length – 1” times Width.

With FoFm Pulse performs a saturation pulse analysis before actinic light is switched on.

Add recovery appends a recovery curve to an induction curve (see Section 4.2.10, page 58 for information on recovery times).

4.3.3.6 Light Curve

Light Curves are defined in “Light Curve Settings” (Fig. 53, page 82). Properties of Width and Add Recovery are identical for light and induction curves.

Intensity specifies the actinic intensity setting for the first light step (range 1 to 5, for PAR values see Fig. 39, page 61).

Length is the number of light steps which can range from 2 to 12. If length = 5 and intensity = 2, 5 light steps with intensity settings 2, 3, 4, 5, and 6 will be performed. The time required for a light curve is Length times Width.

Light Curve Settings		Side Panel	
Width (mm:ss)	1:00		START
Initial Intensity	3		▲
Length	8		▼
With FoFm pulse	on		▲ , ▼ . Up/down keys move cursor.
Add Recovery	yes		SET. Selects item which then can be changed by arrow keys.
			SET
		EXIT	

Fig. 53: Light Curve Settings

4.3.3.7 Yield + # Poro

This item allows porometer measurements without yield determination between combined yield and porometer measurements. The number entered corresponds to the number of measurements

made by the Porometer alone. For instance, if the “Yield + # Poro” is set to 4, each combined yield plus porometer measurement is followed by 4 porometer measurements without yield determination.

4.3.3.8 Select Y(II), Y(NPQ), Y(NO) or Y(II), NPQ, qL

The last two lines of the menu “Program/Clock Settings” affect graphical and numerical presentation of data in windows Induction Curve, Light Curve, and Recovery Curve. Selecting “Y(II), Y(NPQ) Y(NO)” displays three yield parameters that are used in analyzing energy partitioning. Choosing Y(II), NPQ, and qL displays the classical NPQ parameter and a parameter for indicating the reduction state of PS II (qL). Selection between lines works as described for “Clock Item” (Fig. 50, page 79). See Table 29, page 174 for definitions of fluorescence parameters.

4.3.4 Sensors

Sensor Settings		<div>Side Panel</div> <div>▲ , ▼ . Up/down keys move cursor.</div> <div>SET. Selects item which either open a submenu (lines with →) or select the PAR sensor (selection marked by X).</div> <div>EXIT. Return to Main Menu.</div>
Ext. PAR Sensor	On/Off	
Add int. PAR	On/Off	
Sel. Ext. PAR Sensor	→	
SETTINGS		
Porometer	→	
O ₂ /pH Sensor	→	
Spectrometer	→	
Leaf Clip	→	SET
Internal PAR	→	EXIT

Fig. 54: Sensor Settings

The Main Menu item “Sensors” leads to the window “Sensor Settings” (Fig. 54). The external PAR sensor (if connected) can be set as the light sensor via the first window item (Ext. PAR Sensor). The second window item allows adding the intensity of the internal light source to the external measurement. In addition, “Sensor Settings” includes eight links leading to submenus (distinguished by →). An overview on Sensor Settings and its complex structure of submenus gives Fig. 55.

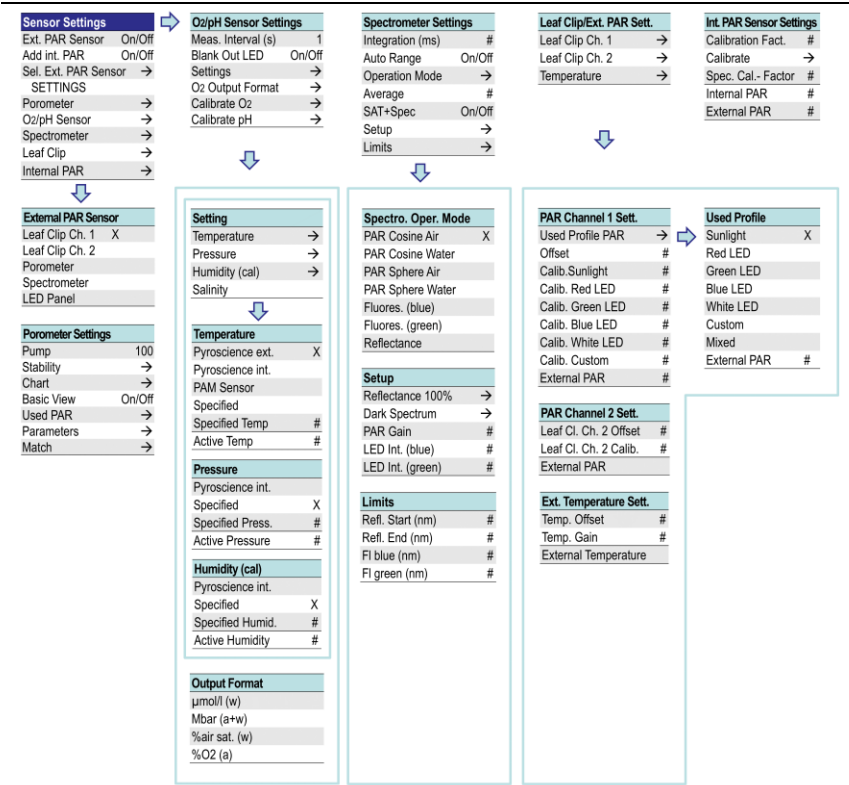


Fig. 55: Sensor Settings Menu and its Submenus

4.3.4.1 External PAR Sensor

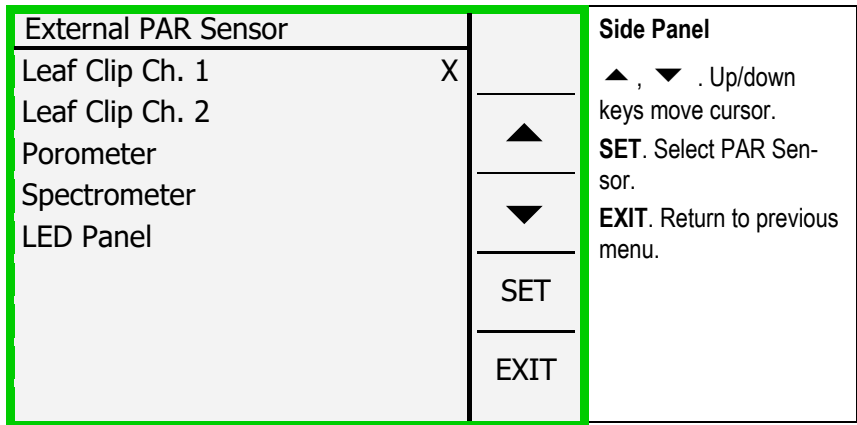


Fig. 56: External PAR Sensor

Ext PAR Sensor opens a list of PAR sensors (Fig. 56, page 85): PAR sensors 1 and (optional) 2 of 2035-B leaf clip, the PAR sensor of the porometer, the miniature spectrometer, and a PAR sensor connected to the LED-Panel RGBW-L084. Select PAR sensor via Up/down keys and **SET**. In the default configuration, Leaf Clip Ch 1 is selected.

4.3.4.2 Porometer

See separate Manual.

4.3.4.3 O2/pH Sensor Settings

The window “O2/pH Sensor Settings” is dedicated to optode sensors measuring either oxygen concentration or pH. (At the time of

printing this manual, the market launch of the pH sensor for the MINI-PAM-II device is in preparation.)

Here, the subject “O2/pH Sensor Settings” is shortly introduced. Details on O₂ optode measurements are provided by a separate manual:

https://www.walz.com/products/chl_p700/mini-pam-ii/downloads.html

and by the manual for the FireStingO2 oxymeter:

(<https://www.pyroscience.com/en/>).

In the window “O2/pH Sensor Settings”, “Meas. Interval” defines the frequency of oxygen readings and “Blank Out LED” switches of MINI-PAM-II internal or external light in the case that it interferes with oxygen measurements.

“Settings” allows entering temperature, pressure, and humidity data for correct sensor calibration. For each of these three parameters, an extra submenu is provided. “Output Format” opens a menu for selection of the parameter used for oxygen representation in the chart.

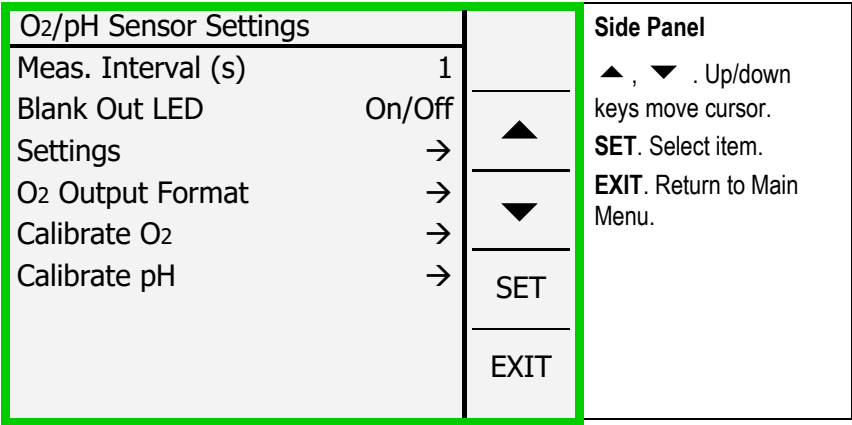


Fig. 57: Oxygen Sensor Settings

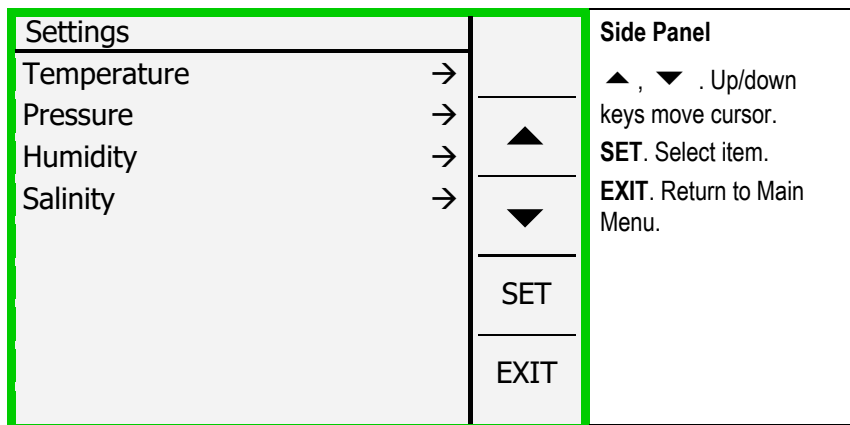


Fig. 58: Settings

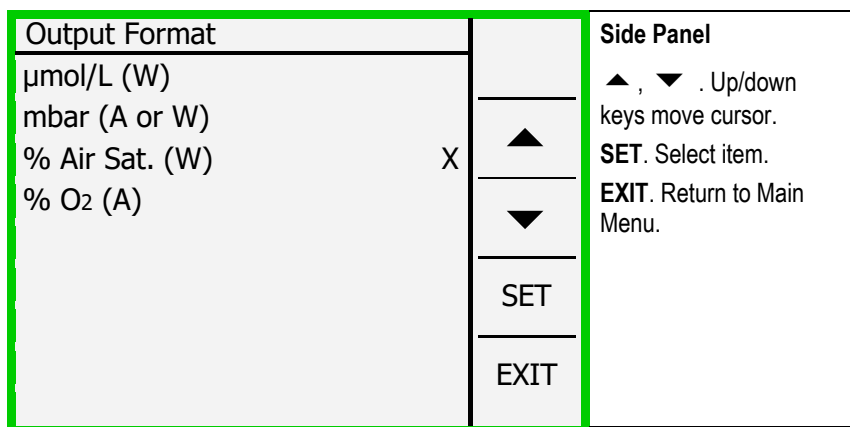


Fig. 59: Output Format

The last two menu items (Calibrate O₂ and Calibrate pH) give access to the windows for calibration of the oxygen and the pH sensor, respectively.

4.3.4.4 Spectrometer

The item “Spectrometer” in menu “Sensor Settings” leads to the window “Spectrometer Settings (Fig. 60, page 88) which is dedicated to configuration on the miniature spectrometer.

Spectrometer Settings		Side Panel	
Integration (ms)	5.00	▲ , ▼	Up/down keys move cursor.
Auto Range	on		SET. Select item.
Operation Mode	→	▼	EXIT. Return to previous menu.
Average	1		
SAT + Spec	off		
Setup	→	SET	
Limits	→		
External PAR		EXIT	

Fig. 60: Spectrometer Settings

Integration time (ms) determines the integrated measuring time in ms for a single spectrum.

Auto Range optimizes the integration time for a spectrum depending on incoming light. Auto Range ON is the default setting. Auto Range OFF is for special applications. When spectra look unrealistic, switch on Auto Range.

Operation Mode opens another menu (Fig. 61, page 89) in which the type of spectrum (light, fluorescence or reflectance) and, for light spectra, the entrance optics (flat=cosine versus spherical) and the environment in which measurements are performed (air or water) can be specified.

The cosine configuration employs a diffusing disk as light entrance (Fig. 10). This configuration shows an approximate cosine response toward incoming radiation. The sphere configuration measures light from all directions with similar weight. At the time of writing of this manuscript, the spherical sensor is under development.

Average determines the number of measurements used to calculate the final spectrum.

SAT+Spec active: each saturation pulse analysis is followed by recording of a spectrum.

Setup opens a menu for recording of the dark spectrum of the miniature spectrometer (Dark Spectrum →). The dark current of the spectrometer at room temperature is measured in the factory and stored on the flash memory of the device. To newly establish the dark current, fully darken the entrance optics (cv. Fig. 10, page 23).

The reference spectrum for reflectance measurements with the white standard is recorded by selecting “Refl. 100%”. The “ampli-

Spectro. Oper. Mode		Side Panel
PAR Cosine Air	X	
PAR Cosine Water		
PAR Sphere Air	▲	
PAR Sphere Water	▼	
Fluores. (blue)	SET	
Fluores. (green)	EXIT	
Reflectance		

Fig. 61: Spectro. Oper. Mode

fication factor “Gain” is automatically adjusted depending on operation mode of the spectrometer. For example, switching the operation mode from air to water increases the Gain from 1.00 to 1.03 to consider the lower light flow to the detector under water. The last two items in the Setup menu allows adjusting the intensity of the blue and green LED used as excitation source for fluorescence emission spectra.

Limits: In the menu “Limits” short and long wavelength borders of reflectance spectra can be defined, and also the short wavelength limits of fluorescence emission spectra.

Setup		Side Panel	
Refl. 100%	→	▲ , ▼	Up/down keys move cursor.
Dark Spectrum	→	SET	SET. Select item.
Gain	1.00	▼	EXIT. Return to previous menu.
LED Int. (blue)	200		
LED Int. (green)	600		
		SET	
		EXIT	

Fig. 62: Setup

Limits			Side Panel
Refl. Start (nm)	400.00		▲ , ▼ . Up/down
Refl. End (nm)	850.00	▲	keys move cursor.
Fl. Blue (nm)	600.00		Change value.
Fl. Green (nm)	640.00	▼	SET . Select item.
		SET	EXIT . Return to previous
		EXIT	menu.

Fig. 63: Limits

4.3.4.5 Leaf Clip

In the Sensors menu, the item “Leaf Clip” opens the Window “Leaf Clip / Ext. PAR Sett.” which deals with the calibration of the Leaf Clip Holder 2035-B (Fig. 64). The factory calibration data are stored on the flash memory of the Leaf Clip Holder 2035-B.

Leaf Clip / Ext. PAR Sett.			Side Panel
Leaf Clip Ch. 1 →			▲ , ▼ . Up/down
Leaf Clip Ch. 2 →		▲	keys move cursor.
Temperatur →			Change value.
		▼	SET . Select item.
		SET	EXIT . Return to previous
		EXIT	menu.

Fig. 64: Leaf Clip / Ext. PAR Sett.

Leaf Clip Channel 1 Settings		Side Panel	
Used Profile	→	▲ , ▼	Up/down
Offset	0		keys move cursor.
Calib. Sunlight	158	▲	Change value.
Calib. Red LED	149	▼	SET. Select item.
Calib. Green LED	181		EXIT. Return to previous
Calib. Blue LED	217	SET	menu.
Calib. White LED	161		
Calib. Custom	150		
External PAR	1007	EXIT	
Used Calibration	149		

Fig. 65: Leaf Clip Channel 1 Settings

Leaf Clip Ch. 1 opens “Leaf Clip Channel 1 Settings” which contains various calibration factors for PAR channel 1, that is, the mini quantum sensor of Leaf Clip Holder 2035-B. The different calibration factors (“Calib. Sunlight to Calib. White”, see Fig. 65) were factory-established and are characteristic for each PAR sensor.

By considering the spectral variations in sensitivity of the PAR sensor, these calibration values are optimized to measure various light qualities. Specifically, Calib. Sunlight is optimized to measure sunlight under clear skies.

Calib. Red, Calib. Green, Blue and Calib. White LED applies to measurements of light from Red, Green, Blue, and White LEDs. Calib. Custom can be chosen by the user. The “Offset” in the window “Leaf Clip Channel 1 Settings” applies to all PAR measurements.

The item Used Profile in the “Leaf Clip Channel 1 Settings window” opens another menu (Used Cal. Profile, Fig. 66, page 93) in which a calibration factor or a calibration profile can be selected. In “Used

Cal. Profile”, the items “Sunlight” to “Custom” correspond to comparable items in the previous window., For example, selecting profile “Sunlight” activates the factor “Calib Sunlight” for PAR measurements.

The ‘Mixed’ profile is designed for the 2054-L external LED light source. It adjust the calibration factor dynamically depending on the contributions of the red, green blue and white LEDs.

Leaf Clip Ch. 2 in window “Leaf Clip/Ext. PAR Sett” (Fig. 67, page 94) opens the calibration menu for a second PAR sensor, which can be connected to the SMA socket on the side of the 2035-B leaf clip. This menu consists of slope (Calib.) and offset of the calibration line. Also, the PAR readout of the active PAR sensor is displayed.

Used Profile		Side Panel
Sunlight	X	
Red LED		
Green LED		
Blue LED		
White LED		
Custom		
Mixed		
External PAR	1007	<p>▲ , ▼ . Up/down keys move cursor. SET. Open submenu. EXIT. Return to previous menu.</p>
Used Calibration		

Fig. 66: Used Cal. Profile

Leaf Clip Channel 2 Settings		<div><div>▲</div><div>▼</div><div>SET</div><div>EXIT</div></div>	Side Panel
Chan. 2 Offset	0		▲ , ▼ . Up/down keys move cursor.
Chan. 2 Calib.	142		SET . Open submenu.
External PAR	1007		EXIT . Return to previous menu.

Fig. 67: Leaf Clip Channel 2 Settings

Temperature leads to the calibration data for the leaf temperature sensor of the 2035-B leaf clip (Fig. 68, page 94). This window lists slope (Gain) and offset of the calibration line.

External Temp. Sensor Sett.		<div><div>▲</div><div>▼</div><div>SET</div><div>EXIT</div></div>	Side Panel
Offset	0.0		▲ , ▼ . Up/down keys move cursor.
Gain	1.00		SET. Select calibration constant; to change calibration constant use up/down keys.
External Temp.	21.6		EXIT. Return to Leaf Clip/Ext. PAR Settings.

Fig. 68: External Temperature Sensor Settings

4.3.4.6 Internal PAR

“Internal PAR” of the “Sensors” menu opens the “Internal PAR Sensor Settings” window (Fig. 69, page 96). This window is dedicated to the calibration of the MINI-PAM-II internal PAR sensor. The window displays the current “Calibration Factor” of the internal PAR sensor. If the miniature Spectrometer is used for calibration of the internal PAR sensor, the spectrometer calibration factor (Spec. Cal. Factor) is important. This factor corrects the effect of inhomogeneous illumination by actinic light of the diffusing disk of the miniature spectrometer.

Typically, the Spec. Cal. Factor is adjusted by comparing the Calibration Factor determined by the PAR sensor of the 2035-B clip with the Calibration Factor determined by the miniature spectrometer in the PAR block (Fig. 9, page 22). The Spec. Cal. Factor is correct if both setups (2035-B clip and miniature spectrometer) result in similar Calibration Factors. Ask Walz for the Spec. Cal. Factor when a 2035-B clip is not at hand.

To calibrate the internal PAR sensor when the MINI-PAM/F1 Miniature Fiberoptics is used, see Section 3.4.14.1 (page 40).

“Internal PAR” opens the menu “Int. PAR Sensor Settings” (Fig. 69, page 96). In this menu, the last two lines display the PAR measured by the internal and by the external PAR sensor ($\mu\text{mol m}^{-2} \text{s}^{-1}$), when internal actinic light is switched on.

Calibration of the internal PAR sensor is done by comparison with a calibrated external sensor. The external sensor is positioned at a defined distance and angle relative to the end of the fiberoptics. Therefore, the readout of the internal PAR sensor applies only for the position of the external PAR sensor. With the external PAR

sensor in calibration position, proper calibration of the internal sensor is apparent from similar readouts of internal and external PAR.

Internal PAR Sensor Settings		Side Panel
Calibration Factor	1136	
Calibrate	→	
Spec. Cal-Factor	1.35	
Internal PAR	415	
External PAR	417	
		▲ , ▼ . Up/down keys move cursor.
		SET. Selects item or open submenu Calibrate.
		EXIT. Return to Main Menu.

Fig. 69: Internal PAR Sensor Settings

If calibration of the internal PAR sensor has previously been carried out but the two PAR values differ clearly, check if:

- (1) External PAR sensor is in the center of the actinic light beam.
- (2) Setup of previous calibration was different from the current one. For instance, if in previous calibration the external PAR sensor was 3 mm away from fiber tip and now the 2035-B leaf clip geometry is used (7 mm instead of 3 mm distance) the internal PAR readout would be much higher than the external one.
- (3) External PAR sensor not connected or incorrect PAR channel selected in window “External PAR Sensor” (Fig. 56, page 85).

PAR sensor compatible with MINI-PAM-II: To calibrate the internal PAR sensor using the PAR sensor of the 2035-B clip connected to the MINI-PAM-II, go to “Calibrate” and establish calibration factor automatically.

MINI-SPEC/MP Miniature Spectrometer: The spectrometer only measures the actinic light intensity properly in the calibration mode. Actinic light in normal operation mode gives a highly fluctuating signal. See Section 3.4.1, page 21, for instructions.

Other PAR sensor: To calibrate the internal PAR sensor using an external PAR sensor which cannot be read by the MINI-PAM-II, adjust calibration factor manually by selecting “Calibration Factor”, pressing **SET** and changing the factor by arrow keys until internal PAR matches that of the external sensor.

MINI-PAM-II Sett. Ft Chart Res. (s) # Auto Off (min:ss) # Backlight (%) Signal LED On/Off Beeper on/off Time/Date → Load Defaults → Time/Date Year # Month # Day # Hour # Minute #	Memory Datasets → New Record → Mark A Record No # Measurement No. # Info MINI-PAM-II → Sensors → Firmware →	MINI-PAM-II Info Hardware Version Serial Number Battery (V) Battery (SAT) (V) Battery (%) Current (A) Battery Type: Temp LED (C) Temp. (Detector) (C) Operation Time (h) Sensor Info Head Keys Mini-Spec Firmware Info Version Number Date Build Date Programmed	Reset Reset Settings → Reset Sys. Settings → Reset Porometer →
---	--	---	--

Fig. 70: Overview: MINI-PAM-II, Memory, Info, and Reset Menus

4.3.5 MINI-PAM-II Settings

The menu “MINI-PAM-II Settings” permits adjusting of device settings, choosing between two time intervals for the Ft chart settings, and retrieving the default configuration for measurements.

Ft Chart Resolution (s)

Ft chart resolution can be either 0.2 or 1.0 s/dot corresponding to 25 or 125 s/total time axis.

MINI-PAM-II Settings		<div>Side Panel</div> <div>▲ , ▼ . Up/down keys move cursor.</div> <div>SET. Change parameter by up/down keys, or switch function on/off, or open submenu (lines with →).</div> <div>EXIT. Return to Main Menu.</div>
Ft Chart Resolution (s)	0.2	
Auto Off (min)	5	
Backlight (%)	30	
Signal LED	on	
Beeper	on	
Time/Date	→	SET
		EXIT

Fig. 71: MINI-PAM Settings

Auto Off (min)

Time interval without saturation pulse analysis after which the MINI-PAM-II powers off.

Backlight (%)

Percentage of maximum intensity of the display’s backlight LED array.

Signal LED

On/off switch for LED on top of the MINI-PAM-II (green flash every 2 s, normal operation; green double flash every 2 s, clock-controlled operation; continuous light, saturation pulse analysis; green flash every 10 s, sleep mode).

Beeper

On/off switch for beeper which acoustically confirms keystrokes and saturation pulse analysis.

Time/Date

Simple menu for setting time and date.

4.3.6 Memory

The “Memory” window provides an overview on stored saturation pulse analyses (Datasets). Importantly, here fluorescence kinetics induced by saturation pulses can be viewed. To see stored induction and light curves, use the **MEM** key of window “Induction Curve” (Fig. 36, page 56) and window “Light Curve” (Fig. 37, page 59), respectively. Also in this window, new records can be started and the mark of saturation pulse data can be changed.

Memory				Side Panel
Datasets	→			▲, ▼ . Up/down keys move cursor. SET . Open submenu (lines with →) or change mark by up/down keys. EXIT . Return to Main Menu.
New Record	→		▲	
Mark	A		▼	
Record No.	68			
Measurement No.	185			
			SET	
			EXIT	

Fig. 72: Memory

Deletion of data from the MINI-PAM-II internal memory is not allowed in the stand-alone mode. However, the memory can be cleared using the software WinControl-3.

4.3.7 Info

“Info” (Fig. 73, page 100) consists of 3 links to submenus of which three list hardware and software information:

- MINI-PAM-II provides hardware information of the fluorometer.
- “Sensors” lists serial numbers of sensors connected to the MINI-PAM-II.
- Firmware shows serial number and date of the firmware of the MINI-PAM-II.

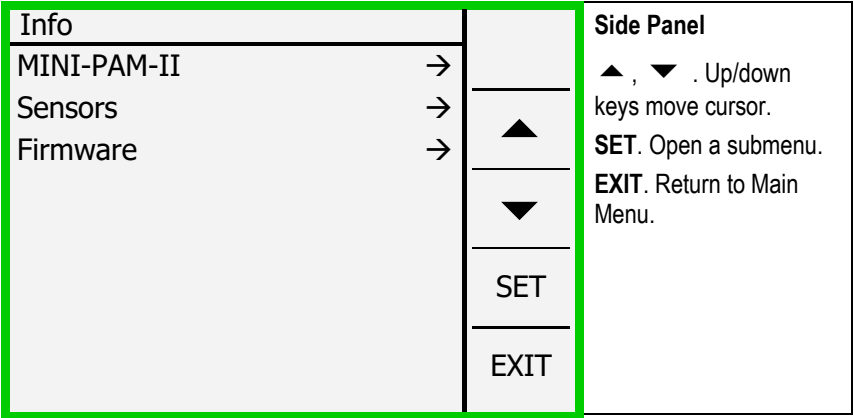


Fig. 73: Info

4.3.7.1 MINI-PAM-II Info

Information available on this window is (1) hardware version of fluorometer, (2) serial number of fluorometer, (3) battery voltage at normal operation, (4) battery voltage during a saturation pulse, (5) charge status in percent, (6) present current consumption (7) temperature (in °C) of actinic LED, (9) temperature (in °C) of detector, and (9) operation time of device.

MINI-PAM-II Info		Side Panel ▲, ▼ . Up/down keys move cursor. EXIT. Return to Info.
Hardware Version	19	
S/N PYAD0191A		
Battery Voltage (V)	7.4	
Battery (SAT) (V)	7.2	
External DC (V)	12.1	
Current (A)	0.04	
Temp. LED (°C)	23.6	
Temp. Detector (°C)	-	
Operation Time (h)	86	EXIT

Fig. 74: MINI-PAM-II Info

4.3.7.2 Sensor Info

Serial number and hardware information of MINI-PAM-II measuring head and accessories.

4.3.7.3 Firmware Info

Version and times of completion of firmware.

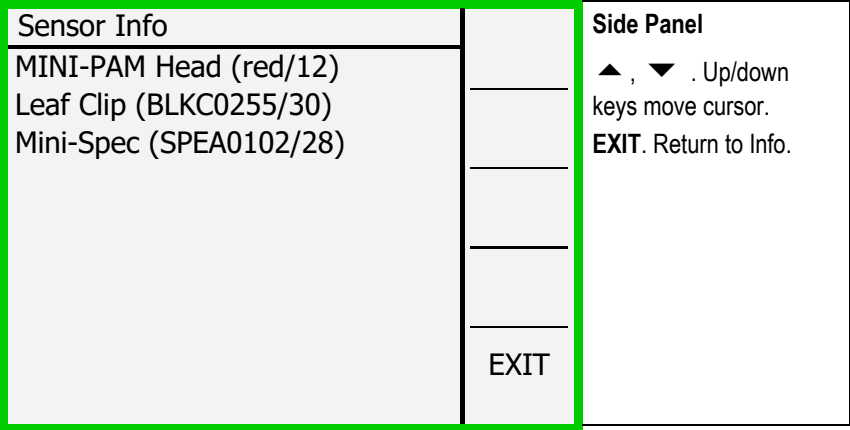


Fig. 75: Sensor Info

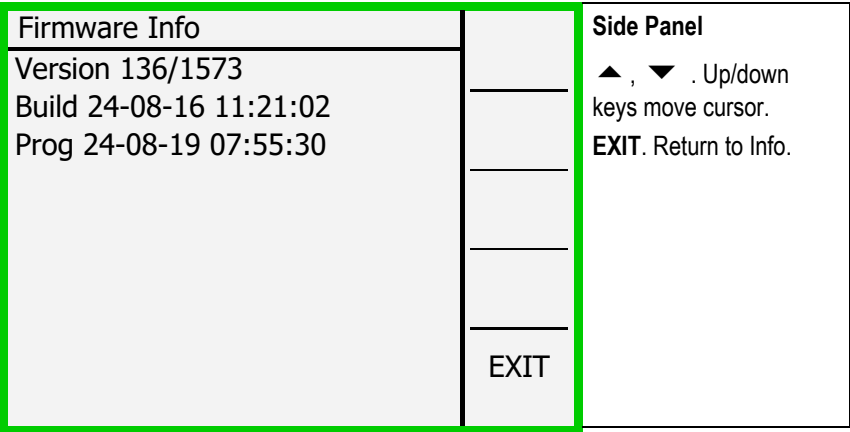


Fig. 76: Firmware Info

4.3.8 Reset

The „Reset“ menu consists of tree items and is used to install default settings:

Reset		Side Panel
Reset Settings →		▲ , ▼ . Up/down
Reset Sys. Set. →		keys move cursor.
Reset Porometer →		EXIT. Return to Info.
	EXIT	

Fig. 77: Reset

4.3.8.1 Reset Settings

Installs the default MINI-PAM-II setting (Table 9, page 103) left and center column. Current settings of some variables are saved when the MINI-PAM-II is shut off. These variables are marked by “Yes” in the rightmost column of Table 9.

Table 9: Default Settings

	Default Setting	Current Setting (saved and restored)
Measuring Light		
Status	On	No
Intensity	6, relative unit	Yes
Frequency	3, see Table 8	Yes

Table 9: Default Settings

	Default Setting	Current Setting (saved and restored)
Frequency high status	Off	No
Actinic Light		
Status	Off	No
Intensity	6, relative unit	Yes
Factor	1.00	Yes
PAM Signal		
Gain	1, relative unit	Yes
Damping	2, relative unit	Yes
Far-red Light		
Width 5	5, s	Yes
Intensity	8, relative unit	Yes
Saturation Pulse		
Intensity	10, relative unit (\triangleq 5000 $\mu\text{mol m}^{-2} \text{s}^{-1}$)	Yes
Width	0.6, s	Yes
Program Actinic Light and Yield		
Actinic light width	30, s	Yes
Initial Pulse	Yes	Yes
Program Induction Curve		
Delay	40, s	Yes
Width	20, s	Yes
Length	12	Yes
Program Light Curve		
Width	20, s	Yes
Intensity	3, relative unit	Yes
Length	8, light steps	Yes
Clock		
Item	Saturation pulse	Yes
Interval	60, s	Yes
Hardware		
Signal LED status	On	Yes
Beeper status	On	Yes

Table 9: Default Settings

	Default Setting	Current Setting (saved and restored)
Automatic power down	15, min	Yes
Background light	60%	Yes
Graphics		
Ft chart time resolution	0.2, s/dot	Yes
External Light		
Status	Off	No
Total intensity	1%	Yes
Red LED	10%	Yes
Green LED	10%	Yes
Blue LED	10%	Yes
White LED	10%	Yes
Stirrer		
Status	Off	No
Speed	10%	Yes
Pre-SAT off	10, s	Yes
Reverse	0, s	Yes
Interval mode	Off	Yes
Interval	2, min	Yes
Stirring interval	5, s	Yes
Stir in program	Off	Yes
F₀' Mode		
Status	Off	Yes
PAR Sensor		
Status	Internal	No
Mark		
Character	A	Yes

4.3.8.2 Reset Sys. Settings

Restores device settings like calibration factor of internal PAR sensor (calibrated for geometry of 2035-B leaf clip), measuring

light current and calibration factors of external devices like those stored on the 2035-B leaf clip

4.3.8.3 Reset Porometer

Restores the factory settings of the porometer. See separate manual.

5 WinControl-3 Installation

The WinControl-3 software is provided on a Walz USB flash drive. The WinControl-3 software is regularly optimized. The latest software version is available on the Walz website:

https://www.walz.com/products/chl_p700/mini-pam-ii/downloads.html

WinControl-3 can be installed from the Walz USB flash drive or using the setup software downloaded from the Walz website.

5.1 Installation process

The installation of WinControl-3 is mostly automatic. Dialog boxes appearing during setup provide advice or allow configuration of WinControl-3. To install WinControl-3, proceed as follows:

- Close other programs as advised by the setup wizard.
- Execute setup file: double-click on file or right-click on file and choose “run” from context menu.
- A pop-up windows must appear which identifies the Heinz Walz GmbH as verified publisher.
- Accept default folder for program installation or choose different folder after clicking **Browse...**.
- Select “Standard” Installation. (The “JUNIOR-PAM Teaching Edition” runs only with JUNIOR-PAM fluorometers.)
- Install USB driver and select optional WinControl-3 links (icon or shortcut).







- Connect MINI-PAM-II to computer and run PAM Firmware Update. If the current firmware* is outdated, PAM Firmware Update will automatically replace it by the recent version. Running PAM Firmware Update after installation of WinControl-3 is important because new software properties may function only in the presence of the latest firmware.

*The term firmware denotes a piece of software residing on a flash memory of the MINI-PAM-II. The firmware is integrated in the WinControl-3 software.

5.2 WinControl-3 Program Group

Setup of WinControl-3 creates the WinControl-3 program group (Table 10, p. 108) in the Windows Start menu. The WinControl-3 program group is comprised of 5 items. The items are introduced in sections 5.2.1 through 5.2.5.

Table 10: WinControl-3 in Windows Start Menu

	WinControl-3
	PAM Firmware Update
	Uninstall WinControl-3
	WinControl-3
	WinControl-3 – Network Mode
	WinControl-3 – Offline

5.2.1 PAM Firmware Update

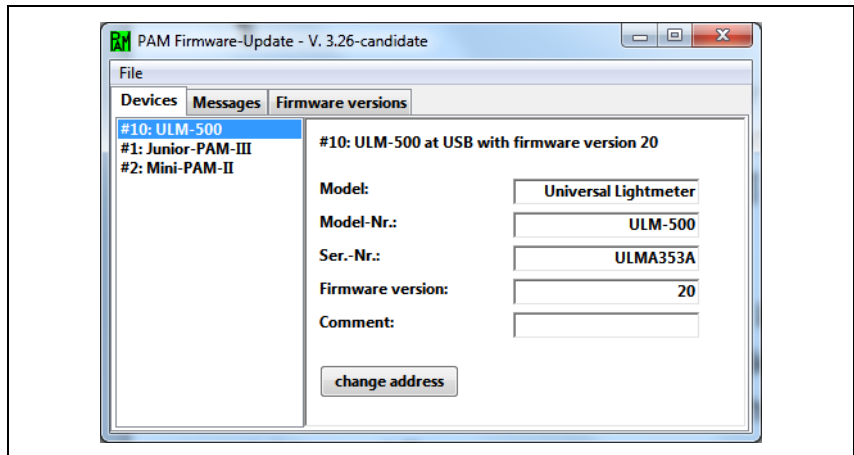


Fig. 78: PAM Firmware Update

Several devices connected to the same computer must have different addresses (channel numbers).

Initialization of PAM Firmware Update triggers a search for PAM devices connected to the computer. The result is displayed in the right panel of window "Devices" (Fig. 78, page 109). Each device name is preceded by its address number (between hash and colon).


"PAM Firmware Update" compares the firmware in the device with the firmware included in the WinControl-3 software. If WinControl-3 includes newer firmware, the device is automatically updated.

"PAM Firmware Update" cannot update firmware of first-generation devices (DIVING-PAM, MICROFIBER-PAM, MICROSCOPY-PAM, MINI-PAM, WATER-PAM). In these devices, firmware resides on an EPROM chip and firmware update requires exchange of this chip.

A device can be selected by mouse click. The currently selected device is highlighted (white letters on blue background). The main panel of the window shows information on the device selected. The first four information lines define the hardware and software state of the device. The last line displays a comment associated with the device and typed in using the WinControl-3 software.

Devices with identical addresses cannot be operated simultaneously. If WinControl-3 detects identical addresses, the window “Device Channel Configuration” (Fig. 79, page 110) pops up offering a working address configuration and the option to change addresses manually. Note that address number is synonymous to channel number in the software WinControl-3.

Fig. 79: Device Channel Configuration

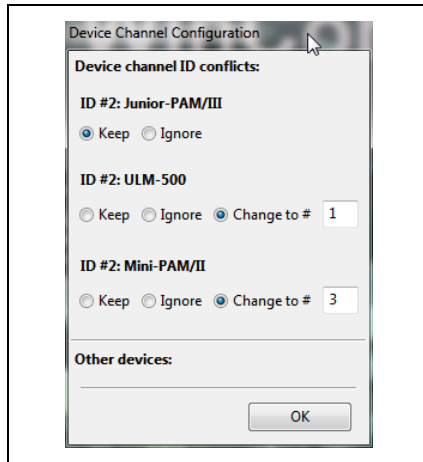

Address numbers can be changed manually via the button **change address**. Then, determine new address by picking a number from the drop-down list “New address:”. If several devices are connected, the drop-down list offers only unused address numbers. The window “Messages” displays the protocol of activities including firmware update of devices. The window “Firmware

Versions” compiles all software version provided by PAM Firmware Update.

5.2.2 Uninstall WinControl-3

This program removes WinControl-3 and all its links. It does not remove the USB driver software.

5.2.3 WinControl-3

This command starts WinControl-3 in the default mode. When devices which are compatible with the WinControl-3-type software are detected, WinControl-3 enters the measure mode. Clicking **Offline-Mode** interrupts the search process and WinControl-3 starts in the offline mode. When the search process fails to find compatible devices, a pop-up window provides three options: **Yes** triggers another search for devices, **No** launches the offline mode of WinControl-3, and **Cancel** ends the whole process. Measuring mode and offline mode, and several instances of WinControl-3 in the offline mode, can run in parallel.

After detection of the MINI-PAM-II, measuring of fluorescence is automatically started. With a green leaf placed in the 2035-B leaf clip, fluorescence values of 400 to 600 are observed (Ft value, bottom of window). Check “Rec. Online” or click **Start Online** to display continuously the Ft on the Chart window. If the Ft is much lower than 400, make sure that the optical fiber is properly inserted. Click **Autoscale** if data are not visible. Trigger saturation pulse analyses by pressing **Fo, Fm** or **SAT**. A healthy leaf, which was kept dark before, should show a value for F_v/F_m of 0.8 or higher.

5.2.4 WinControl-3 Network Mode

The Network mode requires MONITORING-PAM measuring heads and connection via a special interface. It is not available for the MINI-PAM-II.

5.2.5 WinControl-3 Offline Mode

This command launches Wincontrol-3 without the initial search for available PAM devices.

6 WinControl-3 Operation

WinControl-3 offers the same functions as touchscreen operation of the MINI-PAM-II, except that continuous data acquisition and light curve analysis is confined to WinControl-3.


WinControl-3 functions are arranged in 11 Windows (Table 11). The Chart window appears at software start. When more than one device is connected, the additional window “Moni-Bus” appears.

Table 11: Windows of WinControl-3

	Window	Availability	Main panel	Content
1	Chart	Online and Offline	Graphics	Data versus time of all experiments
2	Induct. Curve	Online and Offline	Graphics	Data versus time of fluorescence induction curves
3	Light Curve	Online and Offline	Graphics	Data versus time of light response curves
4	SAT-Chart	Online and Offline	Graphics	Saturation pulse kinetics
5	Spectrum	Online and Offline	Graphics	Spectra
6	Report	Online and Offline	Alphanumerics	Saturation pulse data and data collected at the same time
7	Memory	Online only	Alphanumerics	Information on files stored on the MINI-PAM-II flash memory
8	Batch	Online and Offline	Commands	Site for automatic execution of experiments
9	Control	Online only	Settings	Configuration of external light panel PAR sensor and stirrer control
10	Sensors	Online only	Settings	Control of PAR sensor, spectrometer, and oxygen sensor
11	Settings	Online only	Settings	Device settings

6.1 Chart Window

Fig. 80 divides the Chart window into six areas. Most of the side bars are also present when other windows are viewed.

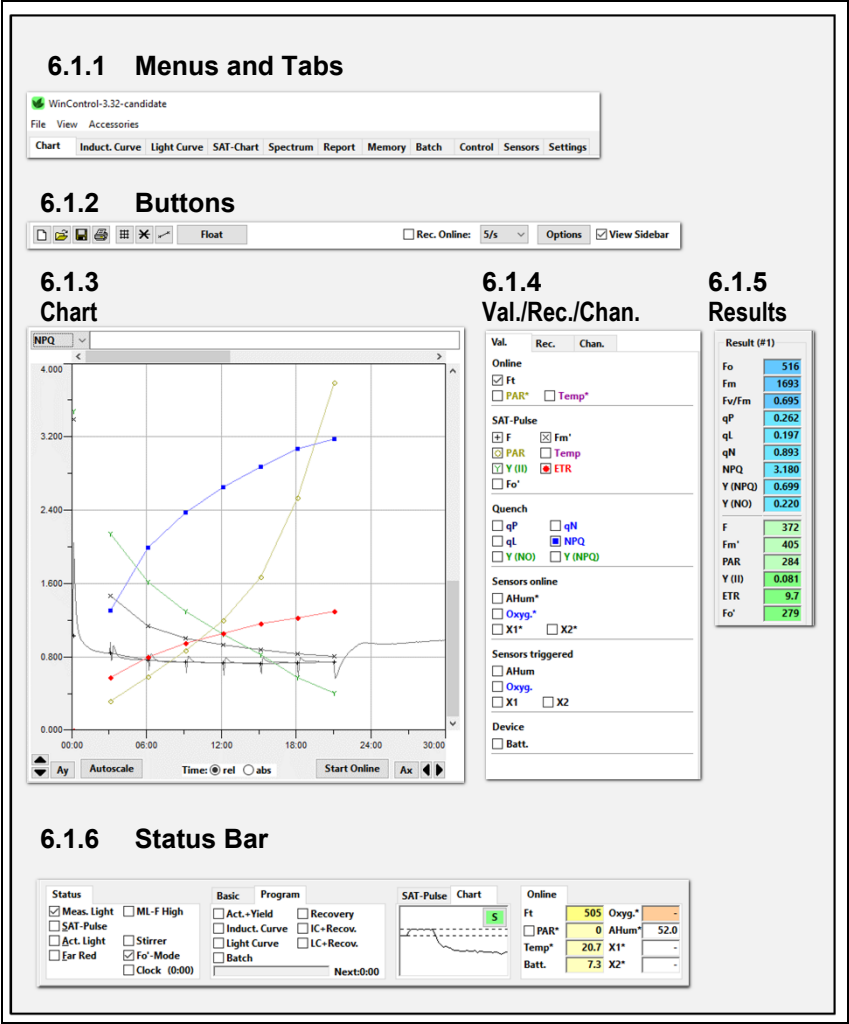

6.1.1 Menus and Tabs

The top bar includes three menus (File, View, and Accessories) and all window tabs (compare Table 11, page 113). The three menus are explained in Table 12.

Table 12: Menu Overview		
Menu	Item	Comment
File	Load Data	Opens WinControl-3 files. Save current data first and switch off online recording of data.
	Save Data	Saves present data in WinControl-3 format.
	Save Settings	Saves all current instrument settings in a batch file. The settings can be restored by executing this batch file.
	Export Report	Exports the data displayed on Report window. For information on configuration of export data see Table 13 (page 116).
	Export Chart Record	Exports the data displayed on Chart.
	Quit	Exit WinControl-3.
View	Results Panel	Switches Results panel (Section 6.1.3) on or off.
	Status Panel	Switches Status panel (Section 6.1.6) on or off.
	Warnings	Prompts the display of the 'program starting time' and non-critical errors.
	Batch Window	Switches Batch File window on or off.
Accessories	Temperature Units	Toggles between °Celsius and °Fahrenheit. The temperature unit affects only the numerical display on the Status bar (6.1.6, p.124).
	Plugins	Does not apply for the MINI-PAM-II.
	Record File	Saves data continuously to reduce data loss in case of program failure. The command prompts for a folder in which the file should be saved. The file name is created automatically using data and time of record start: WinControl-Record-YYYY-MM-DD-hh_mm_ss.pam.

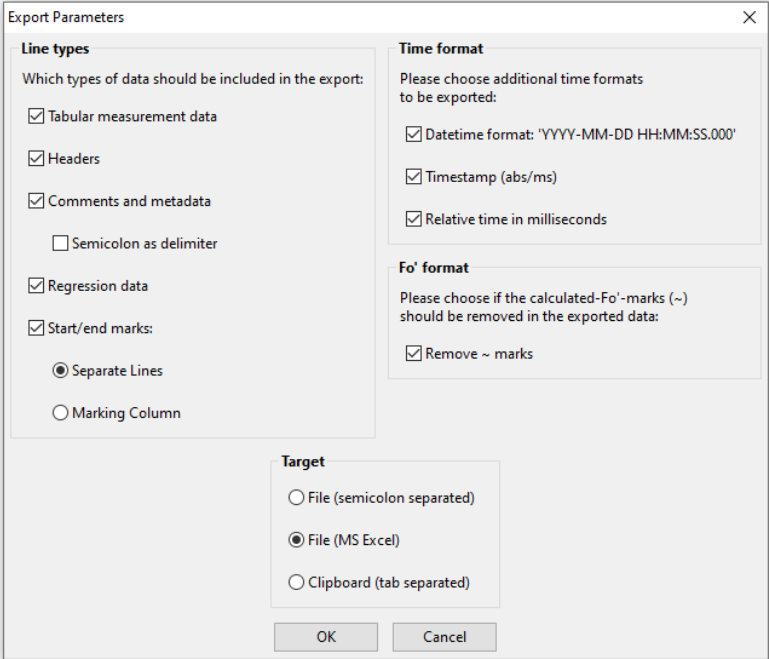


Fig. 81: Export Format for Reports

Table 13: Export Format for Reports

Parameter	Action when checked
<input checked="" type="checkbox"/> Tabular measurement data	Exports data of saturation pulse analysis and all other data recorded at the same time.
<input checked="" type="checkbox"/> Headers	Exports column headers.
<input checked="" type="checkbox"/> Comments and metadata	Exports user and software comments (=“metadata”, e.g., device information).
<input checked="" type="checkbox"/> Semicolon as delimiter	Applies to comments and metadata. Information after a semicolon is written in a new cell.
<input checked="" type="checkbox"/> Regression data	Exports the cardinal numbers of light curves.

Table 13: Export Format for Reports

Parameter	Action when checked
<input checked="" type="checkbox"/> Start/end marks	
<input checked="" type="radio"/> Separate Lines	Writes start and end of an experiment in separate lines.
<input type="radio"/> Marking Column	Writes start and end of an experiment in a separate column.
<input checked="" type="checkbox"/> Datetime format	Exports date and time. Format: day/month/year hours:minutes:seconds.ms. In Excel, custom-formatting may be needed: dd/mm/yyyy hh:mm:ss.000.
<input checked="" type="checkbox"/> Timestamp	Exports UTC (Coordinated Universal Time) time in ms. To convert into date/time use: $= \frac{SourceCell + HoursTimeShift * 3.6 \cdot 10^6}{86.4 \cdot 10^6} + 25569$ Format target cell, e.g.: dd/mm/yyyy hh:mm:ss.000
<input checked="" type="checkbox"/> Relative time in ms	Export time of experiment in ms.
<input checked="" type="checkbox"/> Remove “~” marks	Removes tilde (swung dash) signs. WinControl-3 marks calculated F_0 values by a tilde.
<input checked="" type="radio"/> File (semicolon separated)	Creates a file in which individual data are separated by semicolons.
<input checked="" type="radio"/> File (MS Excel)	Creates an Excel file. Note that columns “Date” and “Time” are differently formatted but contain the same information: dd/mm/yyyy hh:mm:ss.000
<input type="radio"/> Clipboard (tab separated)	Copies the data to the clipboard. Data are separated by tab characters. Use the paste command to copy the data into a spread sheet program.

6.1.2 Buttons

The elements of the buttons bar are introduced in Table 14 and Table 15.

Table 14: Buttons









Icon	Meaning	Comment
	Delete	Deletes all current data.
	Load	Same function as “Load Data” in the File menu (Section 6.1.1.).
	Save	Same function as “Save” in the File menu (Section 6.1.1.).
	Print	Prints current chart view.
	Grid on/off	Controls display of chart grid.
	Lines	Connects data points with lines.
	Symbols	Controls the display of symbols.
Float	Add Chart	Creates an additional chart window with separate view settings.
<input checked="" type="checkbox"/> Rec. Online	Continuous recording	Controls continuous recording of fluorescence, PAR, and temperature.
5/s 	Sampling frequency	Sets sampling frequency for online data. Open drop-down menu by the downward arrow. Available sampling frequencies are 5/s, 1/s, and 1/10s. 5/s is available only for fluorescence.
Options		See Table 15.
Or right-click on chart.		
<input checked="" type="checkbox"/> View Sidebar		Controls display of side bar (see 6.1.4, Val./Rec./Chan.)

Table 15: Options

Menu Item	Present	Comment
Export Record	Always	Exports the data displayed.
Select current record	Always	Highlights data of a Record on the chart, and also on windows SAT-chart and Report.
Split Selection	After selection of interval.*	Put selection in separate record.
Zoom to Selection	After selection of interval.*	Displays the interval enlarged.
Export Selection	After selection of interval.*	Exports the data selected.

* How to select an interval on the chart: keep the left-hand mouse button pressed, move cursor across x-interval and release.

6.1.3 Chart

Fig. 82 outlines the tools to adjust the chart. Section 6.1.4 explains how to select data for display.

The scale of y- and x-axis can be manually changed: a vertical pair of arrows affects the y axis and a horizontal pair of arrows adjusts the x-axis (see lower border of Fig. 82). Right next the two pairs of arrows are buttons for automatic axis scaling.

Manual and automatic y-axis scaling affects only the currently active y-axis unit. In contrast, the command Autoscale adjusts y- and x-axis so that all data fit in the chart area.

The y-axis unit can be picked from a drop-down menu located in the upper left corner of the chart. The active y-axis unit also determines the y-data of the cursor position on the chart.

A way to zoom in to a particular x-axis interval is to mark the x-interval of interest (see Fig. 82), open a context menu by right-click on the chart area, and select the command “Zoom to Selection”.

The selection of data on the chart results in highlighting the corresponding saturation pulse kinetics in the window SAT-Chart and the corresponding data lines in the window Report.

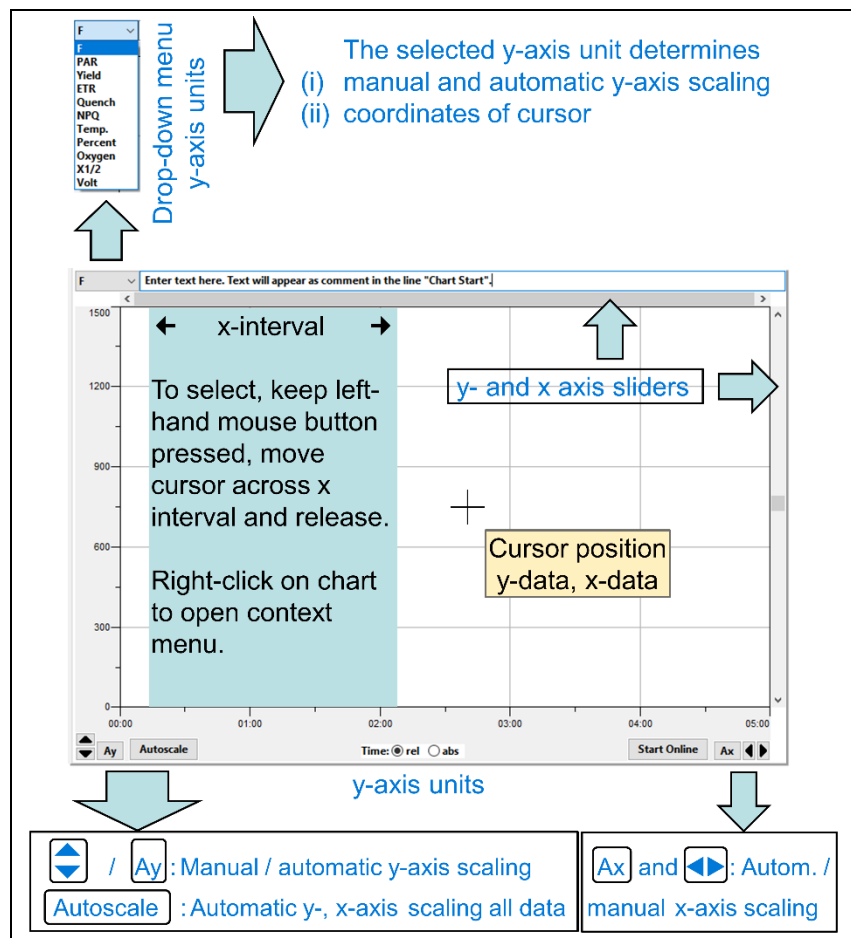


Fig. 82: Chart Tools

6.1.4 Val./Rec./Chan.

The field “Val. Rec. Chan.” includes three different sidebars (Fig. 83). Data which should be displayed on the chart are selected on the Val. (= values) sidebar. The Val. sidebar distinguishes continuously recorded data (data groups: Online and Sensors Online) from saturation pulse data and data taken at the time of a saturation pulse (data groups SAT Pulse, Quench, and Sensors Triggered).

The Rec. (= Record) sidebar lists the number of individual charts of the present data file (= Report). Additional Records can be started by the button **New Record** and existing Records can be deleted by **Delete Record**. The latest Record, or the Record selected by mouse click, will be displayed on chart. The windows “Induction Curve” and “Light Curve” also possess the Rec. sidebar but in these cases the Induction Curve and Light Curve experiments, respectively, are listed.

The Chan. sidebar displays all devices detected by the WinControl-3 software. Only a MINI-PAM-II was connected in Fig. 83.

On the Val. sidebar, the Ft represents a continuously recorded PAM fluorescence signal (online signal). The fluorescence level “F” is the corresponding saturation pulse signal (the F is the Ft measured right before the saturation pulse).

PAR* (photosynthetic active radiation, $\mu\text{mol photons m}^{-2}\cdot\text{s}^{-1}$) and Temp* (Temperature, °C) are two further online signal. They are distinguished by a superscript asterisk from their pendants which are measured together with saturation pulses. In the same way, data of additional sensors are marked (Fig. 83). The additional

sensors available measure humidity (AHum) or oxygen concentration (Oxyg.). The sensor channels X1 and X2 are prepared for future use.

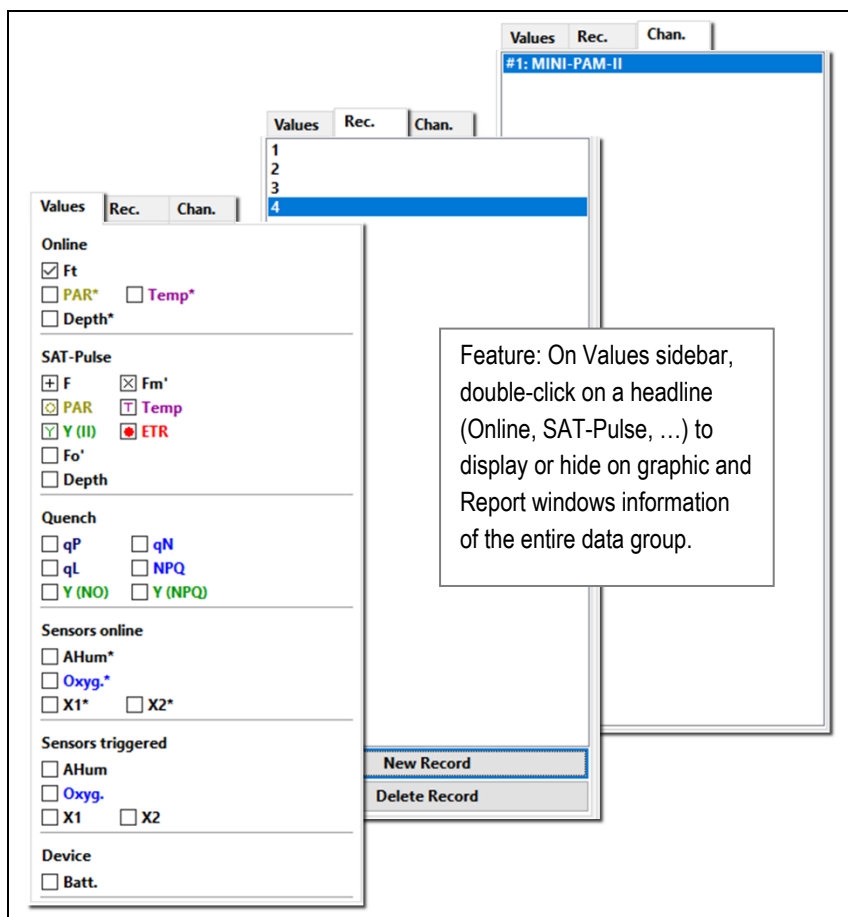


Fig. 83: Val. Sidebar

Besides F level fluorescence, the SAT-Pulse fluorescence levels are maximum fluorescence F_M' (including F_M) and minimum fluorescence F_0' (F_0). Derived from these fluorescence levels is the photochemical quantum yield of photosystem II, $Y(II)$ (including

F_v/F_m). The electron transport rate (ETR, $\mu\text{mol electrons}/(\text{m}^2 \cdot \text{s})$) is calculated with $Y(\text{II})$ and PAR. The section “Quench” of the Val. sidebar includes six more saturation pulse parameters. All fluorescence parameters are explained in Chapter 8 (page 165).

Table 16: Sidebar

Parameter	Sample Data	Comment
Fo	441	Fluorescence properties of the dark-acclimated sample.
Fm	1818	
Fv/Fm	0.757	
qP	0.237	Fluorescence quotients describing the state of the light acclimated sample.
qL	0.147	
qN	0.816	
NPQ	2.212	
Y(NPQ)	0.616	
Y(NO)	0.278	
F	506	Fluorescence properties of the light-acclimated sample.
Fm'	566	
PAR	285	Actinic light intensity, $\mu\text{mol m}^{-2} \text{s}^{-1}$.
Y(II)	0.106	Fluorescence property of the light-acclimated sample.
ETR	12.7	Relative electron transport rate, $\mu\text{mol m}^{-2} \text{s}^{-1}$.
Fo'	313	Fluorescence property of the light-acclimated sample.

6.1.5 Results

The Results sidebar (Table 16) shows numerical values of the current experiment. The upper part of the sidebar displays data of the F_o, F_m determination. These data are maintained throughout the current experiment. All other data are updated with each saturation pulse analysis. All fluorescence levels and ratios are explained in Chapter 8 (page 165).

6.1.6 Status Bar

The “Status” field indicates the state of light sources and special functions (Table 17). The checkboxes function both as indicator and as on/off switches.

The “Basic” and the “Program” fields are stacked and can be selected by tabs. Actinic light and the clock frequency is adjusted on the Basic field, all automatic routines of the WinControl-3 software can be started in the Program field (Table 17).

The stacked “SAT-Pulse” and “Chart” field provide buttons to trigger F_v/F_m or $Y(II)$ analyses, and graphic areas displaying fluorescence kinetics induced by saturation pulses. The “Online” field display live data numerically.

Table 17: Status Bar

Status	
<input checked="" type="checkbox"/> Meas. Light	<input type="checkbox"/> ML-F High
<input type="checkbox"/> SAT-Pulse	
<input type="checkbox"/> Act. Light	<input type="checkbox"/> Stirrer
<input type="checkbox"/> Far Red	<input checked="" type="checkbox"/> Fo'-Mode
	<input type="checkbox"/> Clock (0:00)

Meas. Light: Low frequency PAM measuring light.

ML-F high: High frequency measuring light. Measuring light changes automatically to high frequency when actinic light is switched on.

SAT-Pulse: Saturation pulse analysis to determine $Y(II)$; equivalent to **SAT** button (see below).

Act. Light: Actinic light to drive photosynthesis.

Far-red: Light at wavelengths > 700 nm.

Stirrer: Does not apply for the MINI-PAM-II.

Fo'-Mode: Automatically takes as F_0' fluorescence the minimum fluorescence in a period of far-red-light illumination following a saturation pulse.

Clock: Repetitive trigger of the event specified on

Settings Window (0, page 150). See below on how to adjust the interval between trigger events.

Basic	
Act. Int.	7 μ 285 μ
Clk. Time:	5:00 Memory:
#1: MINI-PAM-II	5 k



Program	
<input type="checkbox"/> Act.+Yield	<input type="checkbox"/> Recovery
<input type="checkbox"/> Induct. Curve	<input type="checkbox"/> IC+Recov.
<input checked="" type="checkbox"/> Light Curve	<input type="checkbox"/> LC+Recov.
<input type="checkbox"/> Batch	
Next:0:11	

Act. Int.: Setting and intensity of MINI-PAM-II light source in $\mu\text{mol m}^{-2} \text{s}^{-1}$.

Clk. Time: Time interval between automatically triggered events in minutes : seconds.

#1: MINI-PAM-II: Device connected (channel).

Memory: Current file size in kilobytes (k).

Experimental Routines

Experimental routines are defined on the

Settings Window (0, page 150). Most routines can be triggered by the clock.

Act.+Yield: Period of actinic illumination terminated by saturation pulse analysis.

Recovery: Dark phase with saturation pulse analyses performed at increasing intervals.

Table 17: Status Bar


Experimental routines continued

IC+Recov.: Induction curve plus dark phase with saturation pulse analyses performed at increasing intervals.

Light Curve: F_0 , F_M determination followed by illumination with stepwise increasing light intensities where each step is terminated by a saturatin pulse analysis.

LC+Recov.: Light Curve followed by a dark phase with saturation pulse analyses.

Batch: Execution of batch file program.

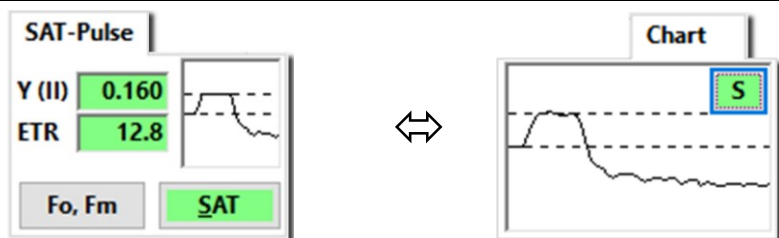

F₀, F_m: Saturation pulse analysis to determine F_0 , F_M , and F_v/F_M .


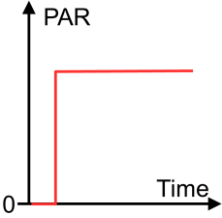
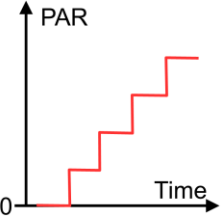
SAT = **S**: Saturation pulse analysis to determine F , F'_0 , F'_M , and $Y(II)$.

Online			
Ft	313	Oxyg.*	-
<input type="checkbox"/> PAR*	0	AHum*	49.0
Temp*	21.3	X1*	-
Batt.	7.4	X2*	-

Online are continuously recorded with the frequency defined (see Table 14, page 118).

Ft, PAR* and Temp*: see 6.1.4, page 121. When PAR* is unchecked, the intensity is taken from the internal PAR list (Section 6.10.3, page 152); when checked, the actually measured value is displayed. Batt., battery voltage.

Table 18: Three Graphics Windows

Chart Window	Induct. Curve Window	Light Curve Window																																													
<u>Rec. Sidebar</u> <table> <tr> <th>Val.</th><th>Rec.</th><th>Chan.</th></tr> <tr><td></td><td>1</td><td></td></tr> <tr><td></td><td>2</td><td></td></tr> <tr><td></td><td>3</td><td></td></tr> <tr><td></td><td>...</td><td></td></tr> </table> <div>Start Online</div>	Val.	Rec.	Chan.		1			2			3			...		<u>Rec. Sidebar</u> <table> <tr> <th>Val.</th><th>Rec.</th><th>Chan.</th></tr> <tr><td></td><td>IC 1</td><td></td></tr> <tr><td></td><td>IC 2</td><td></td></tr> <tr><td></td><td>IC 3</td><td></td></tr> <tr><td></td><td>...</td><td></td></tr> </table> <div>Start IC</div>	Val.	Rec.	Chan.		IC 1			IC 2			IC 3			...		<u>Rec. Sidebar</u> <table> <tr> <th>Val.</th><th>Rec.</th><th>Chan.</th></tr> <tr><td></td><td>LC 1</td><td></td></tr> <tr><td></td><td>LC 2</td><td></td></tr> <tr><td></td><td>LC 3</td><td></td></tr> <tr><td></td><td>...</td><td></td></tr> </table> <div>Start LC</div>	Val.	Rec.	Chan.		LC 1			LC 2			LC 3			...	
Val.	Rec.	Chan.																																													
	1																																														
	2																																														
	3																																														
	...																																														
Val.	Rec.	Chan.																																													
	IC 1																																														
	IC 2																																														
	IC 3																																														
	...																																														
Val.	Rec.	Chan.																																													
	LC 1																																														
	LC 2																																														
	LC 3																																														
	...																																														
<u>Lighting</u> 	<u>Lighting</u> 	<u>Lighting</u> 																																													
<u>Right click on chart</u> <div> Export Record Select current record </div>	<u>Right click on chart</u> <div> Export Record Select current induction curve </div>	<u>Right click on chart</u> <div> Export Record Select current light curve Export Regression Data </div>																																													
<u>Val. Sidebar</u> Quench <input type="checkbox"/> qP <input type="checkbox"/> qN <input type="checkbox"/> qL <input type="checkbox"/> NPQ <input type="checkbox"/> Y (NO) <input type="checkbox"/> Y (NPQ) Sensors online	<u>Val. Sidebar</u> Quench <input type="checkbox"/> qP <input type="checkbox"/> qN <input type="checkbox"/> qL <input type="checkbox"/> NPQ <input type="checkbox"/> Y (NO) <input type="checkbox"/> Y (NPQ) Sensors online	<u>Val. Sidebar</u> Quench <input type="checkbox"/> qP <input type="checkbox"/> qN <input type="checkbox"/> qL <input type="checkbox"/> NPQ <input type="checkbox"/> Y (NO) <input type="checkbox"/> Y (NPQ) Regression <input type="checkbox"/> REG1 <input type="checkbox"/> REG2 Sensors online																																													
<u>x-axis units</u> Time: <input checked="" type="radio"/> rel <input type="radio"/> abs	<u>x-axis units</u> Time: <input checked="" type="radio"/> rel <input type="radio"/> abs	<u>x-axis units</u> <input type="radio"/> Time <input checked="" type="radio"/> PAR																																													

6.2 Induct. Curve/Light Curve Windows

The icons and bars of the Chart window introduced above are also present in the Induction Curve and Light Curve windows. In contrast to the Chart window, which displays all data of a Record, the Induction Curve window displays individual induction curves, and the Light Curve window displays individual Light Curves. The same Record can include several Induction and Light Curves. Use up and down arrow keys to scroll through the list of curves.

Table 18 summarizes further differences between the three windows. The windows have different start buttons and only the Light Curve windows possesses PAR as x-axis unit. The button **Start IC** starts an experiment in which a sample is exposed to a single defined light intensity, the button **Start LC** initiates a routine in which a sample is exposed to incrementally increasing light intensities. The x-axis radio button “Time” of the Light Curve window corresponds to the radio button “rel” (relative time) in the two other windows.

The Light Curve window allows fitting model functions to experimental data. The experimental data are the electron transport rates (ETR) plotted against the photon flux density, PAR. WinControl-3 provides two model functions called REG1 and REG2 which are introduced in Fig. 84 and Fig. 85, respectively. The function REG1 can decrease at high PAR values. Therefore, REG1 can consider photoinhibition of photosynthesis, where β is a photoinhibition parameter. In contrast, REG2 is a rectangular hyperbola which cannot describe photoinhibition.

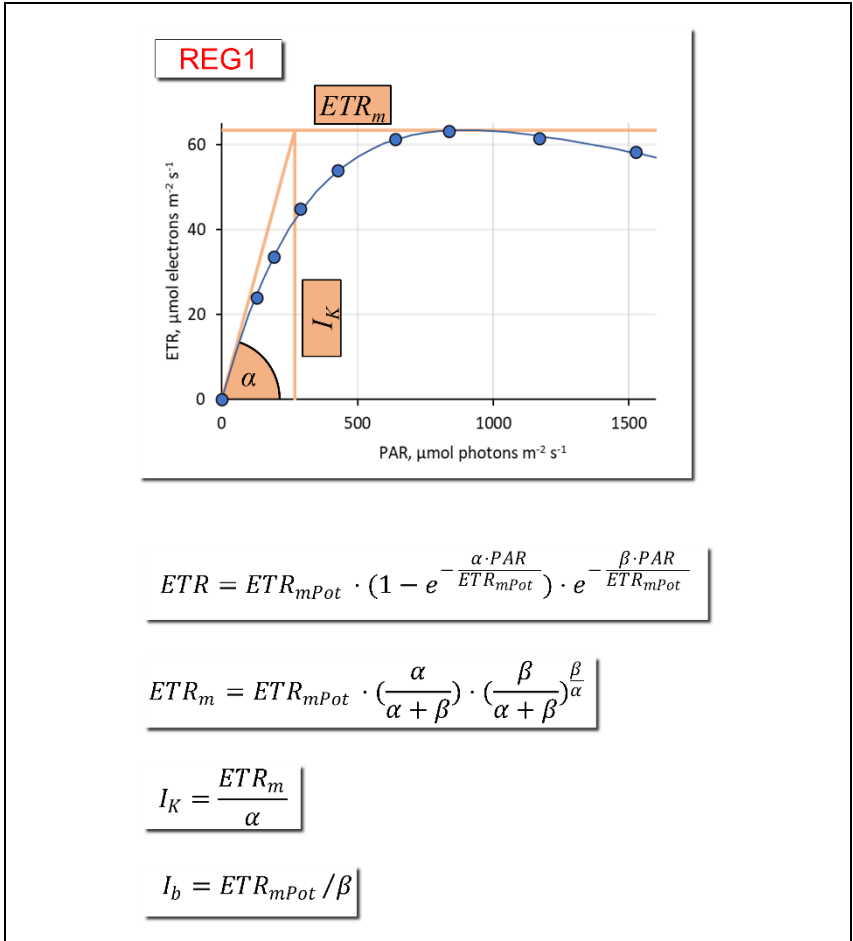


Fig. 84: Model Function REG1

The three cardinal points of the light curve are indicated (α , ETR_m , and I_K). The decrease at $PAR > 1000 \mu\text{mol m}^{-2} \text{s}^{-1}$ is frequently attributed to photoinhibition of photosystem II by strong light. I_b is the theoretical PAR at which the light curve reaches $1/e$ of ETR_{mPot} . ETR_{mPot} is the ETR_m in the absence of photoinhibition. According to: Platt T, Gallegos CL, Harrison WG (1980) Photoinhibition of photosynthesis in natural assemblages of marine phytoplankton. J Mar Res 38: 687-701

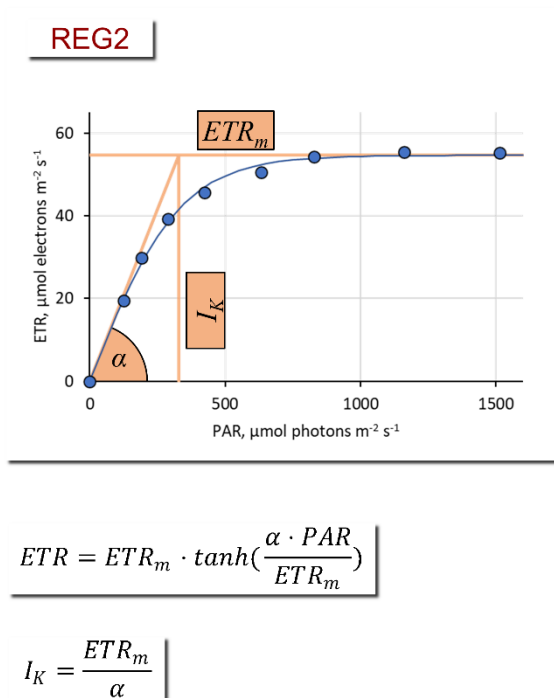


Fig. 85: Model Function REG2

The three cardinal points of the light curve are indicated (α , ETR_m , and I_K). According to Jassby AD, Platt T (1976) Mathematical formulation of the relationship between photosynthesis and light for phytoplankton. *Limnol Oceanogr* 21: 540-547.

Both models calculate the three cardinal parameters of a light curve:

- (i) α , electrons/photons: Initial slope of RLC which is related to the quantum efficiency of photosynthesis.
- (ii) ETR_m , $\mu\text{mol electrons m}^{-2} \cdot \text{s}^{-1}$: Maximum electron transport rate.

(iii) I_K , $\mu\text{mol photons m}^{-2}\cdot\text{s}^{-1}$: Idealized PAR value at which light-limited photosynthesis becomes light-limited.

The cardinal parameters are written into the Report. To export cardinal parameters separately, right click on the chart of the Light Curve window and select from the menu “Export Regression Data” (see Table 18). The item “Select current light curve” in the same menu highlights the data of the currently displayed light curve in the Chart and Report windows. A similar function is available for the window Induction Curve.

6.3 SAT-Chart Window

The SAT-Chart window displays saturation pulse kinetics of fluorescence (Fig. 86). The automatically determined values of F and F_M' are shown as dashed horizontal lines. Fluorescence traces are automatically scaled to fit into the coordinate system. The x-axis covers about 2 s, the saturation pulse starts at 200 ms, and the distance between data points is 50 ms. The mouse wheel scrolls through graphs.

Each graph is accompanied by a protocol panel. The protocol panel lists the pulse number (Nr) and the line number of the Report (Rep. Nr). Select a fluorescence trace by double-click in the protocol window. The selected graph automatically moves to the top of the window.

The **Options** menu (upper right corner of the SAT-Chart window) links the actual view of saturation pulse kinetics to the saturation pulse events selected in one of the three previous windows. The option “Follow Selection” automatically displays on top of the window the saturation pulse kinetics of marked events. When Follow

Selection” is inactive, the option “Jump to Selection” brings the selected kinetics into view.

The Val. sidebar is similarly configured as described in Section 6.1.4 (page 121), except “Online Data” are omitted and the electron transport factor (ETR-F.) is added. The ETR-F. is the fraction of absorbed photons by the sample relative the PAR to which the sample is exposed. The ETR-F. is a factor of the equation for electron transport (Chapter 8, page 165). Each item selected on the Val. sidebar is numerically displayed below the corresponding saturation pulse kinetics.

Single saturation pulse kinetics can be exported by right-click in its chart field. To export a series of saturation pulse kinetics, open the PAM file with a spread sheet program. In the text file, saturation pulse kinetics are preceded by an SG (Table 19). Sort the lines of the file by SG. Create a time scale by considering the neighboring points are spaced apart by 50 ms.

6.4 Spectrum Window

The window is active in the presence of spectral data. With the cursor placed inside the drawing area, or below the x-axis, the mouse wheel has a zoom function for the x-axis. When the cursor is placed left of the y-axis, the mouse wheel has a zoom function for the y-axis.

Right click on the drawing area to open a menu containing zoom and export commands (Fig. 87). The “Zoom to selection” is available after selection of an x-axis interval described before (Fig. 82).

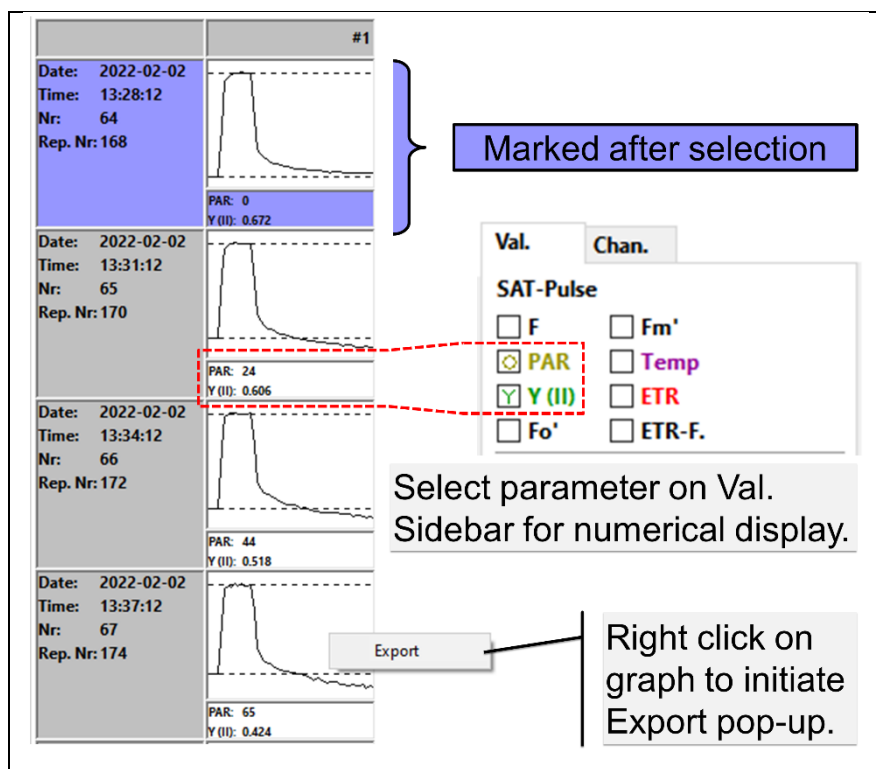


Fig. 86: SAT-Chart Window

Table 19: SAT Kinetics in Raw File

To extract a series of saturation pulse kinetics, open PAM file with spread sheet program, sort by SG and copy data right of SG.

	A	B	C	D	...	AD	AE	AF	AG	AH
1	Date	Datetime	Type	No.						
2	08/03/2022	08/03/2022 09:17:42.997	F	2	...	SG	50	419	419	...
3	08/03/2022	08/03/2022 09:17:44.389	F	3	...	SG	50	383	383	...
						SAT Graph Label	Time interval in ms	Data of Saturation pulse kinetics		

An application menu is available in the top right corner of the window (Fig. 87). To change between the three principal applications (PAR, fluorescence, reflectance), the spectrometer MINI-SPEC/MP has to be reconfigured (see Fig. 10, page 23).

The dark current of the spectrometer at room temperature is measured in the factory and stored on the flash memory of the device. To newly establish the dark current, fully darken the entrance optics (cv. Fig. 10, page 23) and press **Calibr. Dark**.

The standard PAR configuration is “PAR cosine”. This configuration employs a diffusing disk as light entrance (Fig. 10). The configuration shows an approximate cosine response toward incoming radiation. The spectrometer calibration differs between measurements in air and in water. Therefore, choose PAR cosine air or PAR cosine water for proper measurements under the two conditions.

“PAR sphere” employs a spherical light entrance. The sphere configuration measures light from all directions with similar weight. At the time of writing of this manuscript, the spherical sensor is under development.

The setting “PAR Open sensor” is used by service engineers.

“Fluorescence (blue)” and “Fluorescence (red)” measures fluorescence emission spectra excited by blue or red light, respectively. Measuring “Reflectance” requires that the 100% reflectance signal has been established with the highly reflective reference material provided with the spectrometer (Fig. 11, page 25).

The sidebar also display the **Integration time (ms)** which is the integrated measuring time for a single spectrum. ☒ Auto Range optimizes the integration time for a spectrum depending on incoming light. Auto Range ON is the default setting. Auto Range OFF is for special applications. When spectra look unrealistic, switch

on Auto Range. **Average** determines the number of spectra averaged to yield the final spectrum. ☒ **SAT+Spec** measures a spectrum for saturation pulse analysis.

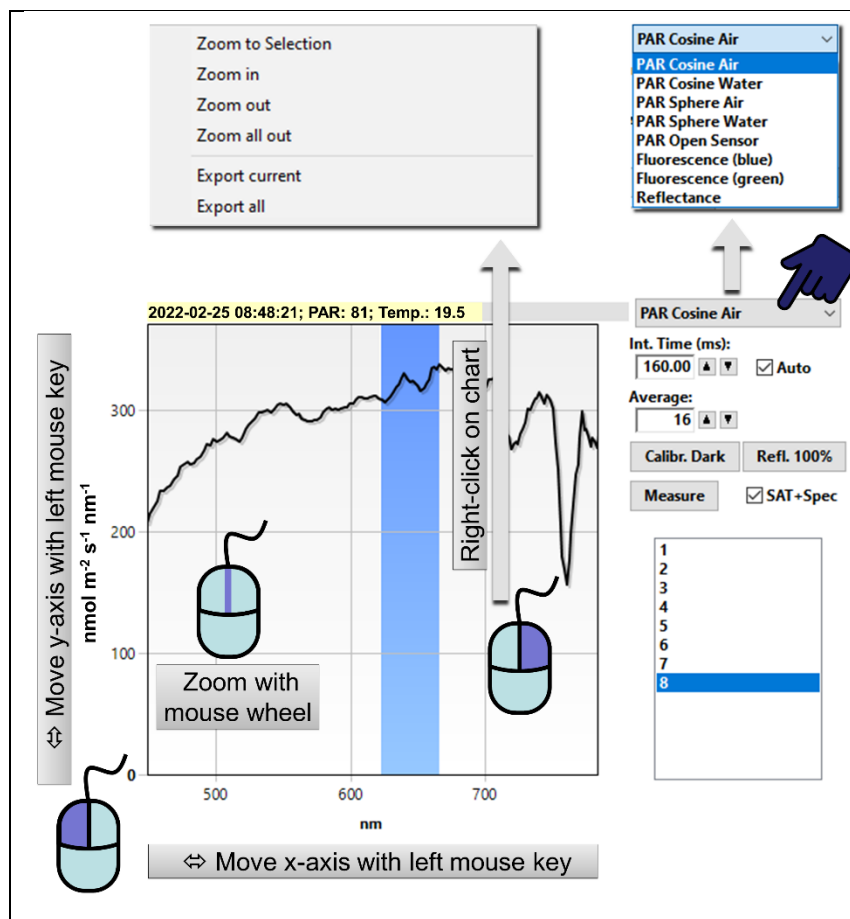


Fig. 87: Spectrum Window

6.5 Report Window

The Report window lists all data associated with saturation pulse analyses. Continuously recorded signals, saturation pulse kinetics and spectra are not listed. The fact that a spectrum was recorded is indicated by “SPEC” in the column “Type” (Table 20). The Report window also displays start and end of induction and light curves, as well as the cardinal points of light curves (Table 20).

The types of data displayed is controlled by the Val. side bar, which is configured as described for the SAT-CHART window (Section 6.3, page 131). Only data displayed on the Report table are exported.

Table 20: Report Table

Type: D, device. FO, F_0 and F_M determination. F, determination of F, F_0' , and F_M' . SPEC, spectrum. SICWS and SICSE, induction curve start and end, respectively. SLCS and SLCE, light curve start and end, respectively. REG1 and REGS, cardinal parameters of regression analysis 1 and 2, respectively (see 6.2). **Mark:** Letters assigned to data (see 4.2.1, page 46). Hide Mark column via option menu. Here, the MINI-PAM-II has the channel number #2. “#2” is preceding the cardinal values of light curve analysis, REG1 and REG2.

	Date	Time	Type	No.	2:Mark	2:F	2:Fm'	2:PAR	2:Y (II)	2:ETR
1			D		Device Nr: #2, MINI-PAM-II (PYAD0191A)					
2	2022-03-08	09:07:11	SCHS							
3	2022-03-08	09:16:20	FO	1	A	461	1932	6	0.761	1.9
4	2022-03-08	09:16:50	F	2	A	680	774	6	0.121	0.3
5	2022-03-08	09:17:24	SPEC		PAR: 12; Temp.: 21.2					
6	2022-03-08	09:17:42	SICS		Induction Curve start					
7	2022-03-08	09:17:48	SICE		Induction Curve end					
8	2022-03-08	09:17:59	SLCS		Light Curve start					
9	2022-03-08	09:18:07	REG1		#2: alpha: 0.362, ETRm: 13.99, Ik: 38.634 (beta: 0.001, ETRmPot: 14.179)					
10	2022-03-08	09:18:07	REG2		#2: alpha: 0.272, ETRm: 13.665, Ik: 50.280					
11	2022-03-08	09:18:05	SLCE		Light Curve end					

The options menu of the Report window can be opened by clicking the **Options** button or by right-click in the Report field. The options menu contains various tools for handling the Report table. All menu items and their function are summarized in Table 21.

Command “Insert Settings” of the options menu writes the current device settings in the report. The abbreviations used for the various settings are compiled in Table 22.

Table 21: Options Menu	
Options Menu Item	Comment
	General Report Management
Follow Selection	Automatically displays the data associated with events marked in other windows
Show Mark	Display marker letters (see “Mark” in Table 20)
Insert Settings	Writes settings of MINI-PAM-II in Report table (see Table 22)
	Manipulation of All Data
Export All	For details see Section 6.1.1, page 115.
Delete All Measure Data	Self-explaining
Page Setup for Printing	Basic configuration of print layout. Checks if output fits on page
Preview Printing	Self-explaining
Print Report	Selects and configures printer, prints current Report
	Manipulation of Selected Data
	Requires that saturation pulse events on a chart or lines of the Report have been selected. To select, move the cursor with left mouse key pressed over events
Export Selected Lines	Confines export to lines of interest

Table 21: Options Menu	
Options Menu Item	Comment
Jump to Selection	Brings data into view, when "Follow Selection" is off
Delete Selected Data	As above. Affects only selected lines.
Preview Print Selection	As above. Affects only selected lines.
Print Selected Data	As above. Affects only selected lines.
Mark as Light Curve	Combines a series of saturation pulse analysis into a light curve and performs regression analysis REG1 and REG2. The series must not contain start or end marks of previous light or induction curves.

Table 22: Abbreviations for Settings

MEA	Measuring Light Intensity
MI	Measuring Light Frequency
DI	Damping
GA	Gain
EF	ETR-Factor
FZ	F Offset
CW	Clock Interval
SI	Saturation Pulse Intensity
SW	Saturation Pulse Width
AI	Actinic Light Intensity
AF	Actinic light factor
FRI	Far-red Intensity
FRW	Far-red Width
AW	Length of exposure of "Act. + Yield" routine
ICD	Delay of Induction Curve
ICW	Interval Between Saturation Pulses of Induction Curve
ICL	Length of Induction Curve
LCW	Step Length of Light Curve
LCI	Initial Intensity Setting of Light Curve
LCL	Length of Light Curve
LO	Offset of External PAR Sensor
LG	Calibration Factor of External PAR Sensor
LO2	Offset of External PAR Sensor 2

Table 22: Abbreviations for Settings

LG2	Calibration Factor of External PAR Sensor 2
TO	Offset of External Temperature Sensor
TG	Gain of External Temperature Sensor
ILO	Offset of Internal PAR Sensor
ILG	Gain of Internal PAR Sensor
PARGAIN_RED	Calibration Factor for Red LED of 2054-L
PARGAIN_GREEN	Calibration Factor for Green LED of 2054-L
PARGAIN_BLUE	Calibration Factor for Blue LED of 2054-L
PARGAIN_WHITE	Calibration Factor for White LED of 2054-L
PARGAIN_CUSTOM	Calibration Factor for Other light Sources
ITO	Offset of Internal Temperature Sensor
ITG	Gain of Internal Temperature Sensor
MLC	Fm Correction Factor 1
MLC2	Fm Correction Factor 2
TRM	Trim Value for Measuring Light
TRA	Trim Value for Actinic Light
TRSA	Tim Value for Saturation Pulse
TRFR	Trim Values for Far-red Light

6.6 Memory Window

The Memory window accesses the data of the internal memory of the MINI-PAM-II. The MINI-PAM-II memory is a circular buffer which can store 512 files or 30 000 data entries maximally. The memory is full when one of the two limits is reached. For instance, if each file contains only two data entries, the full memory contains 512 files but only 1024 data entries, or, if each file contains 1000 data entries, then only 30 files fit into the memory.

When the memory is full, the oldest data are overwritten with the latest ones. Usually, data acquired in the offline mode are downloaded (6.6.2) and the memory is cleared (0) which avoids that all memory is used.

6.6.1 Clear Memory

Device Time:
 Date: 2022-03-03
 Time: 11:14:21
 Sync with PC
 Files:
 Number of files: 30
 Current filename: 30
 Clear Memory

Clear Memory

Are you sure that you want to clear the complete device memory?

OK Cancel

Clear Memory

Deleting the device's internal memory can not be undone and thus all data will be lost. Do you want to abort the operation now?

Yes No

6.6.2 Download Memory

File details:

Create new file Download selected files

Showing page 1 out of 3 pages

001:	2021-01-18 14:04:23 - 2021-11-11 12:23:32	data entries: 258
002:	2021-11-12 10:06:07 - 2021-11-12 10:06:07	data entries: 2
003:	2021-11-12 10:07:04 - 2021-11-26 15:46:37	data entries: 381
004:	2021-11-30 13:49:12 - 2021-11-30 13:49:12	data entries: 2
005:	2021-11-30 13:49:12 - 2021-11-30 13:49:12	data entries: 2
006:	2021-11-30 13:49:12 - 2022-03-03 09:53:50	data entries: 5532
007:	2022-03-03 09:55:50 - 2022-03-03 09:56:33	data entries: 18
008:	2022-03-03 09:56:34 - 2022-03-03 10:07:40	data entries: 14
009:	2022-03-03 10:07:46 - 2022-03-03 10:08:13	data entries: 16
010:	2022-03-03 10:08:16 - 2022-03-03 10:08:48	data entries: 12

Download

Please choose:

☒ Clear existing data in WinControl before Download

☐ Merge with current data in WinControl

Caution: Merging data will destroy any Light Curve or Induction Curve information in the actual data.

OK Cancel

Max. time distance

Please choose the maximum timedistance in which measures should have the same measure number:

☐ 1 second

☐ 10 seconds

☐ 30 seconds

☒ 1 minute

☐ 5 minutes

OK Cancel

Fig. 88: Memory Window

6.6.1 Clear Memory

Clearing the memory is irreversible. To avoid inadvertent deletion of data, two warnings must be passed before the command **Clear Memory** is executed.

Above the Clear Memory button, memory information is given: “Number of Files” is the file count, “Current file number” is the serial number of the file to which data are currently written. The Current file number can be greater than the number of files when the memory is full and old files are overwritten by new ones.

The Clear Memory field display date and time of the internal computer of the MINI-PAM-II. Click **Sync with PC** to apply the computer time settings to the MINI-PAM-II.

6.6.2 Download Memory

A list of files is displayed in the Download field of the Memory window (Fig. 88). You can download individual or groups of files. To pick several files, hold down **Ctrl** key and click on files of interest. To pick consecutive files, hold down **Shift** key and click on first and last files of the row. Then click **Download selected files**.

Usually, data are written in an empty Report window. Selecting “Clear existing data ...” in the download dialogue box, empties the Report window.

When parallel measurements of two different fluorometers should be combined, choose “Merge with current data ...”. The newly imported data will be sorted so that time points of import data matches time points of existing data. Two time points are consid-

ered as matching when their time difference is smaller as the interval selected from the box “Max. time distance” (cf. Fig. 88). The merge operation deletes the report entries for begin and of an induction or a light curve.

6.7 Batch Window

Batch files automatically execute experimental procedures. To activate the Batch window, click icon new batch or open existing batch file (Fig. 89, “Batch Start Buttons”). The click **Edit** and choose between “Add command” and “Record Macro”. The command “Update indentation” is an automatic editing tool to improve readability.

6.7.1 Add Command

Add command leads to five groups of commands (Fig. 89). A command can be inserted into a batch file by left click.

Measure commands

Saturation pulse	Saturation pulse analysis of light-exposed sample
F_0 , F_M determination	Saturation pulse analysis of dark acclimated sample

Light Sources commands

Controls for actinic and measuring light. The light ON command asks for the light intensity: simple enter the light intensity setting in the grey-shaded field. New intensity settings can be chosen using the command “Change intensity”.

Settings commands

“Reset settings to default” installs the settings described in Table 9 (page 103).

Repetition commands

The group contains two commands. “Repeat specific number of times” and “Repeat indefinitely”. Both commands write two lines on the batch file sheet. In case of the first command, these lines are:

Line 1:

```
for $loopvar = 1 to count // Start repetition block with specified number of repetitions
```

Line 2:

```
next // End of repetition block
```

Write commands to be repeated between Line 1 and 2. Define how often the commands should be repeated by the number entered in field `count`.

The second command (Repeat indefinitely) repeats the commands placed between the two lines until the batch program is stopped manually.

Line 1:

```
while 1 // Start block repeating indefinitely
```

Line 2:

```
wend // End of repetition block
```

Timing

The command “Delay” inserts a time interval after the previous command has been terminated. The next command is executed when the time interval ends.

The command “Continue Delay” takes the end of the previous delay phase as starting point. Actions within this time interval are

performed without affecting the interval defined for Continue Delay. A continuous time scale can be built by a series of Continue Delay commands.

“Wait until time of day” delays start of the batch program until the time specified.

6.7.2 Record Macro

The function “Record Macro” converts your manually entered commands into batch file lines. Simply click “Record Macro”, perform experiment, and click “Record Macro” again.

6.7.3 Options

All items of the options menu of the Batch window are related to the MONI-DA data acquisition system.

Check MONI-DA Compatibility

Some batch file commands cannot be executed by the MONI-DA. This command searches for such incompatibilities.

Upload to Device

Transfers a batch file to the MONI-DA memory.

Export

Export the current batch file as “WinControl-3 Compiled Batch File” (*.wccb). This file format is for future use. Upcoming MONI-DA versions will be enabled to directly download wccb files, that is, without being connected to WinControl-3. The final goal is to install batch files by remote control.

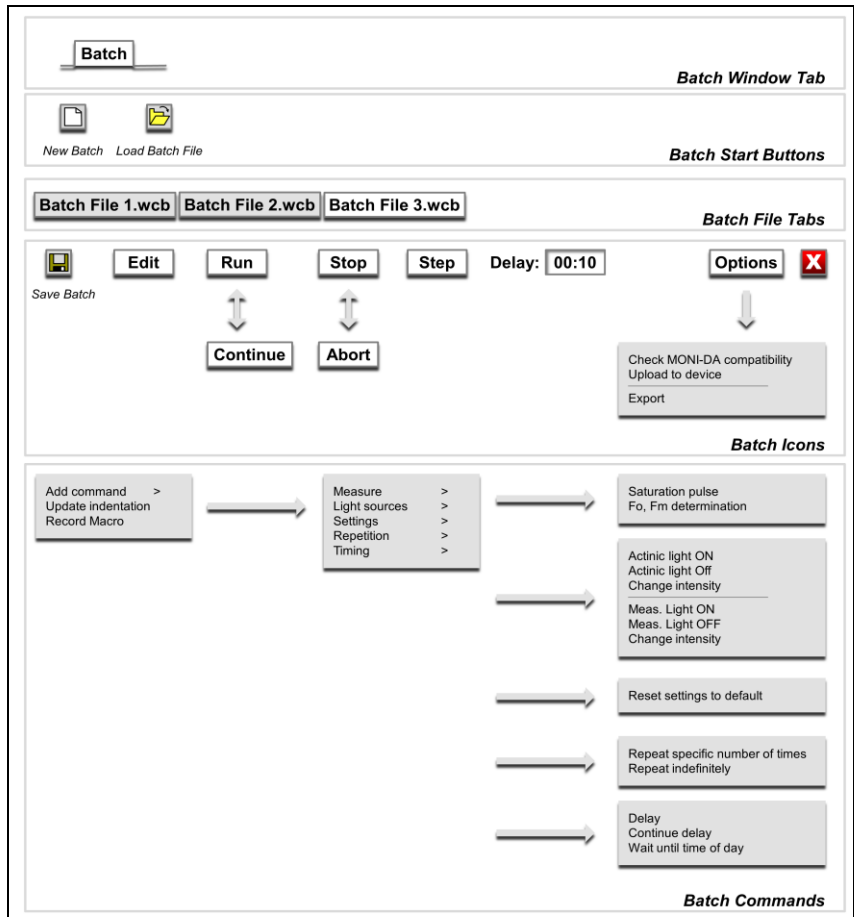


Fig. 89: Batch Window Overview

6.8 Control Window

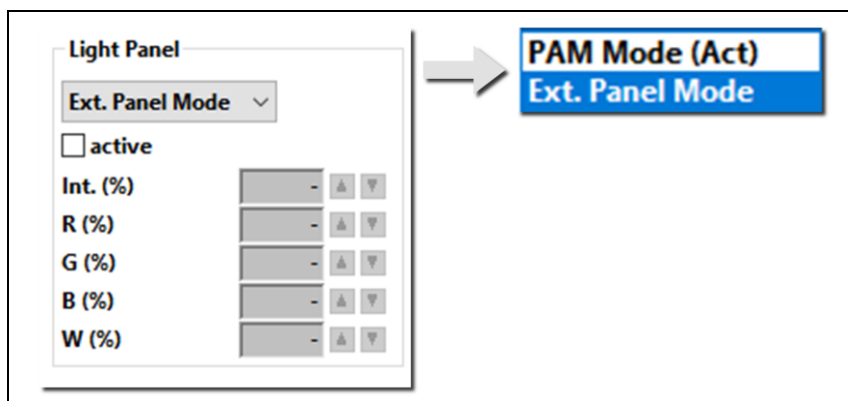


Fig. 90: Control Window

Fig. 90 shows the section of the Control Window which is important for the MINI-PAM-II fluorometer. The drop-down menu on top of the Light Panel area consists of two points: “PAM Mode (Act.)” and “Ext. Panel Mode”. Ext. Panel Mode is the standard mode in which the fluorometer’s internal light source is used as actinic light to which an external light source may be added (see Table 23). Activating PAM Mode (Act.) selects the external light source as the sole actinic light. The meaning of numerical inputs (Int. (%) to W (%) varies depending on the mode chosen (Table 23).

Typically, the maximum intensity of all four LED groups is 5000 $\mu\text{mol m}^{-2} \text{s}^{-1}$ or higher. Operation of the 2054-L External LED Source requires line power.

Table 23: External Light Source

	Ext. Panel Mode	PAM Mode (Act.)
Internal light source of MINI-PAM-II		
Function	Main actinic source	Not available
Intensity setting	Actinic light intensity (Settings window)	Not available
External LED Light Source 2054-L		
Requirement	Not mandatory	Required
Function	Background illumination	Main actinic source
Intensity setting	Int %	Actinic light intensity (Settings window)
Control		
Int. (%)	Relative intensity (0-100%). The number multiplies the intensity of all light sources of the External LED Light Source 2054-L	Not active
R (%), G (%), B(%), W (%)	Absolute intensity in percent (is multiplied by Int. (%))	Factor defining the intensity with which a light source contributes to total intensity.

6.9 Sensors Window

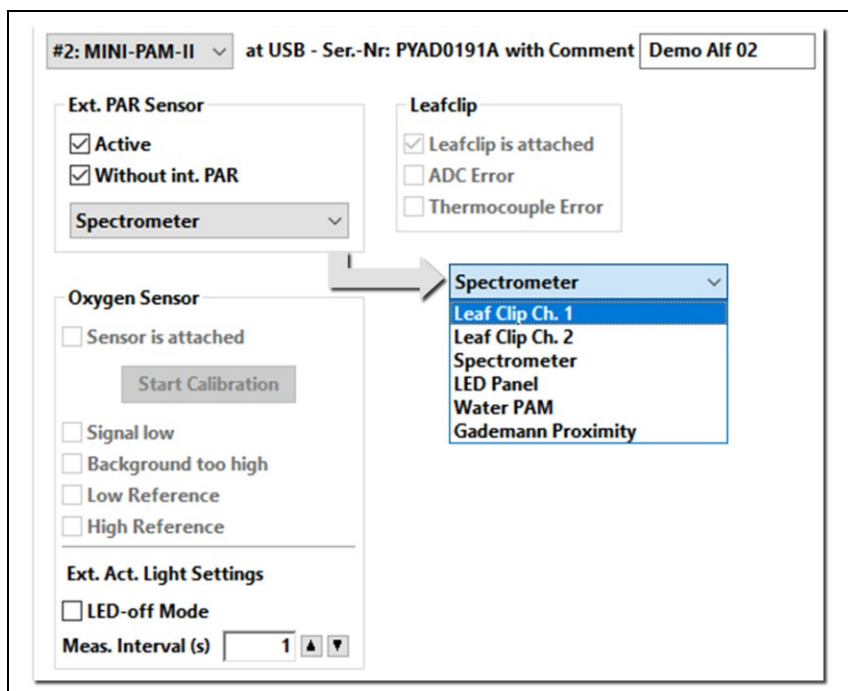


Fig. 91: Sensors Window

External PAR Sensor

The internal PAR sensor measures the light intensity of internal light at sample level. Internal actinic light is used for short-term illumination (Act. + Yield), or when induction and light curve experiments are carried out. External PAR sensors are designed to measure the intensity of external light at sample level. External PAR sensors are employed to estimate electron transport rates driven by external artificial or natural light.

The checkbox “Active” selects sensing of external PAR. Checking “Without int. PAR” ignores measurements by the internal PAR sensor. The drop-down menu below these checkboxes lists external PAR sensors.

Drop-Down Menu

“Leaf Clip Channel 1” corresponds to the micro quantum sensor of the 2035-B clip (Fig. 13, page 26). “Leaf Clip Channel 2” is the signal of a PAR sensor connected to the side port of the 2035-B clip. “Spectrometer” uses PAR spectra integrated over the visible range as PAR value. “LED panel” is the PAR sensor connected to the LED-Panel RGBW-L084. (Does not apply for External LED Light Source 2054-L). “WATER-PAM” is available when working with this PAM fluorometer. “Gademann Proximity” is a sensor under development.

Leaf Clip

The checkbox of “Leafclip is attached” automatically indicates that a 2035-B clip is connected to the MINI-PAM-II. Erroneous conversion of the analogue measuring signal to digital information is signaled by “ADC Error”. A defective thermocouple activates the checkbox of “Thermocouple error”.

Oxygen sensor

This section, including “External Actinic Light Settings” is dedicated to the accessory Fiber-Optic Oxygen Meter FireStingO2 (see separate manual).

6.10 Settings Window

6.10.1 Title Bar

#2: MINI-PAM-II at USB - Ser.-Nr: PYAD0191A with Comment Demo Alf 02

6.10.2 Light Intensity

Measuring Light		SAT-Pulse	
Int.	6	Int.	10
Freq.	3	PAR	
		Width	0.6
			μ
		Actinic Light	
		Int.	3
		PAR	
		Factor	1.00
			65μ
		Far Red Light	
		Int.	8
		Width	5

6.10.3 Actinic Light List

Act. Light List	
0:	0
1:	25
2:	45
3:	65
4:	90
5:	125
6:	190
7:	285
8:	420
9:	625
10:	820
11:	1150
12:	1500
<input type="button" value="Calibrate"/>	

6.10.4 F-Offset

F-Offset
 F-Offs. 19

6.10.5 System Parameter

System Parameter
 Damp. 2
 Gain* 1
 ETR-F. 0.84

6.10.6 Programs

Act.+Yield
 Width 0:30
☒ With Initial Pulse

Induct. Curve
 Delay 0:40
 Width 0:20
 Length 12
☒ With FoFm Pulse

Light Curve
 Width 0:20
 Int. 3
 Length 8
☒ With FoFm Pulse

6.10.7 Clock

Clock
 Time 5:00
 1: SAT-Pulse
 1: SAT-Pulse
 2: Act.+Yield
 3: Light Curve
 4: LC+Rec.
 5: Induct. Curve
 6: IC+Rec.

6.10.8 Indicators

Set
☒ Sign. LED active
☐ Beeper active

Fig. 92: Settings Window

6.10.1 Title Bar

The Reset button restores factory values for all settings of the current window. Located right of the Reset button is a drop-down menu listing all devices connected. The settings displayed on the current window belong to the device selected, and also the serial number which is displayed right of the drop-down menu. In the text box right of the serial number, up to 20 characters can be entered. Serial number and text box content are written in the first line of each record file.

6.10.2 Light Intensity

When actinic light is switched off, μ s measuring pulses are delivered at 5 to 25 Hz depending on frequency setting "Freq." (Table 8, page 72). Switching on actinic light automatically increases the measuring light frequency to 100 Hz. The measuring light frequency can also be manually set to 100 Hz by checking "MF-F High" on the Status Bar (Table 17, page 125).

The intensity setting "Int." adjusts measuring light intensity. Twelve intensity levels are available. The numerical value of the setting is proportional to the measuring light intensity. How to estimate the integrated measuring light intensity for the frequency and intensity settings used is described in Section 4.3.1.2 (page 70).

Note that an increase in flash intensity increases the F_0 fluorescence level (the fluorescence intensity, in general is, proportional to the flash intensity). An increase in the number of flashes/pulse increases the F_0 level only if some PS II reaction centers become closed by the increased integrated measuring light intensity (the measuring light becomes actinic). If this is the case, the integrated

measuring light intensity must be reduced either by reducing intensity or by reducing frequency.

Like measuring light, the intensity of saturation pulses, actinic light and far-red light can be adjusted in 12 levels. The PAR of saturation pulses is indicated in newer versions of WinControl-3. The PAR of actinic light is taken from the light list (see below). Most of far-red emission is not absorbed by chlorophylls. Therefore, intensity data for far-red radiation (PAR or photon flux density) is not given as it would strongly overestimate its actual actinic effect.

6.10.3 Actinic Light List

Twelve actinic light levels are listed. These values are target PAR values. As delivered from the factory, the light emission of the LED is adjusted so that these target values are present at the sample level of a 2035-B Leaf Clip Holder. The target values can be varied by changing the “Factor” in the field “Actinic Light” (Section 6.10.2).

Calibrate newly adjusts the emission of the LED. Calibration must be performed when the sample level is closer or further away than the sample level of a 2035-B Leaf Clip Holder (see Section 4.2.11, page 61). Calibration is also necessary when the MINI-PAM/F1 Miniature Fiberoptics is used. (See Section 3.4.14.1, page 40.)

6.10.4 F-Offset

See Section 4.3.1.7 (page 73).

6.10.5 System Parameter

Damping (Damp.) Function for data smoothing. The smoothing effect increases with increasing numerical value, at the same time, but instrument response decreases. See 4.3.1.4 (page 72).

Gain Factor (Gain). See 4.3.1.3 (page 71). The Gain factor amplifies signal and noise. At low signal levels, increasing the gain improves digital resolution.

ETR-Factor (ETR-F.). The ETR-Factor is the percentage of light absorbed by the sample and is a parameter for calculating the electron transport rate (Section 8.3, page 172).

6.10.6 Programs

Act. + Yield: Light exposure followed by saturation pulse analysis

“Width” is the duration of light exposure. The actinic light intensity adjusted in the field “Actinic Light” applies (Section 6.10.2, page 151). A saturation pulse analysis is always carried out at end of illumination. Checking “With Initial Pulse” places an additional saturation pulse analyses before illumination. The fluorescence levels of the initial saturation pulse analysis are defined as F_0 and F_M . All fluorescence ratio parameters requiring these two fluorescence levels (Table 29, page 174) are only available with initial saturation pulse.

Induct. Curve: Fluorescence curve following the onset of light

Three parameters determine the sequence of events. “Delay” indicates the dark time interval between F_0 and F_M determination and onset of light. The delay time can be adjusted between 5 seconds and 10 minutes. Default value is 40 seconds. “Width” is the

time interval between neighboring saturation pulses in the light period. Length is the number of saturation pulse analyses carried out. Time interval of actinic illumination is approximately “Width” times “Length”. A dark period can be appended to an induction experiment (Table 17, page 125 and Table 6, page 61).

Light Curve: Saturation pulse analyses after exposure to different light intensities

The “Light Curve” program exposes a sample to consecutively increasing actinic light intensities. Three input values define the procedure. “Width” is the duration of exposure to a particular light intensity. “Int.” is the first light level of the light curve. “Length” is the number of light steps. The duration of a light curve is determined by “Width” times “Length”. A dark period can be appended to illumination (see above).

☒ **With FoFm pulse:** Both Induction and Light Curves can be performed without initial saturation pulse (F_0 and F_M measurement).

6.10.7 Clock

The clock utility repetitively triggers one out of 6 different measuring routines. functions: Saturation pulse analysis, Act.+Yield, Induct. Curve, IC+Recov., Light Curve and LC+Recov (sse Above). The interval time between two events (Clk. Time) can be adjusted between 10 s and 50 min. The time interval must be longer than the time required by the triggered event.

6.10.8 Indicators

The signal LED located above the touchscreen indicates the operational state of the MINI-PAM-II as listed in Table 24. The

beeper acoustically marks begin and end of a saturation pulse analysis.

Table 24: LED Signal Code

LED Color	Frequency	Process
M e a s u r i n g	m o d e	
Green	1 Hz	Normal operation
Green double flash	1 Hz	Normal operation and clock running
Green	continuous	Normal operation and saturation flash
Additional codes during firmware update		
Green/red alternating	high	Waiting for software
Red	continuous	Update running

System information is displayed on the bottom of the Settings and the System Settings windows (see Table 25 below)

Table 25: System Information

System information displayed on bottom of Systems and Systems Settings window.

Headline	Model	Model Number	Serial Number
Display	MINI-PAM-II (blue) or	MINI-PAM-II/B or	PYAC####(A) or
	MINI-PAM-II (red)	MINI-PAM-II/R	PYAD####(A)
Information	Color version	Order code	S/N convention

6.11 System Settings Window

Different from all other windows of WinControl-3, the window System Settings is not represented in the row of tabs. To access system settings, open Settings window and click “System Settings” (Fig. 93). The functional elements of System Settings are shown in Fig. 94 together with their section numbers.

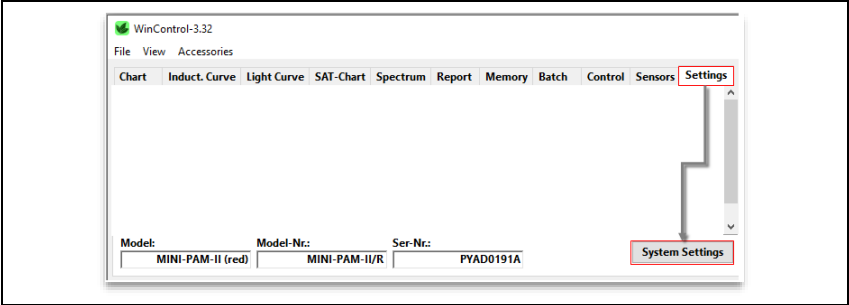


Fig. 93: System Settings Button
Click to open System Settings window.

Reset

#2: MINI-PAM-II

at USB - Comment

Demo A1f 02

Reset complete System

Do you really want to load the default system settings?
This is only recommended for advanced users!

OK Abbrechen

Int. PAR-Sensor

☒ Active

Offset -

Calib. 1044

Calibrate

PAR sensor calibration

Do you really want to calibrate the internal PAR sensor of the PAM?

OK Abbrechen

6.11.2 Custom PAR Gain

Cust. PAR Gain

Sunlight

Red 142

Green 166

Blue 190

White 155

Custom 150

Sunlight

Red

Green

Blue

White

Custom

Mixed

6.11.4 Ext. PAR/T sensors

Ext. PAR-Sensor

☐ Active

Offset 0

Calib. 152

Channel 2:

Offset 0

Calib. 150

Ext. Temp. Sensor

Offset -0.7

Gain 1.02

6.11.5 PAM Trim Val.

PAM Trim Values

Fm-Corr. 0

Fm-Corr. 2 0

Meas. Light 186

Act. Light 200

SAT Light 200

FR Light 153

6.11.6 Internal Temperature Sensor

Int. Temp. Sensor

Offset -

Gain -

Fig. 94: System Settings

156

6.11.1 Title Bar

The **Reset** button of the System Settings page recalls all calibration data established at the factory. This information is taken from the flash memories of the MINI-PAM-II and of the external sensor (e.g., the 2035-B Leaf-Clip Holder). The original calibration data cannot be changed by the user. The command Reset on the System Settings window is equivalent to the command “Reset System Settings” of the MINI-PAM-II menu (Section 4.3.8.2, page 105). The command Reset on the Settings window (Section 6.10.1, page 151) is equivalent to the command “Reset Settings” of the MINI-PAM-II menu (Section 4.3.8.1, page 103).

The other elements of the Title Bar are described in Section 6.10.1, page 151.

6.11.2 Custom PAR Gain

Different calibration factors for different light sources are provided. The calibration factor varies because of slight spectral variations in sensitivity of the PAR sensor. Table 26 summarizes the factors available.

6.11.3 Int. PAR sensor

The internal PAR sensor received a small fraction of the actinic light emitted by the internal LED of the MINI-PAM-II. The calibration factor of the internal PAR sensor is displayed in the “Calib.” numerical field (Fig. 94).

The calibration of the internal PAR sensor is done by comparison with a calibrated external PAR sensor. The reading of the external PAR sensor depends on its position relative to the end of the optical fiber of the MINI-PAM-II. Factory calibration was established with a 2035-B leaf clip in which the fiber was fully inserted. For other configurations, the internal sensor must be newly calibrated.

PAR sensor compatible with MINI-PAM-II: To calibrate the internal PAR sensor using the PAR sensor of the 2035-B clip, connect clip to MINI-PAM-II, position light sensor relative to fiber identical to the experimental situation (with the light sensor in the sample plane) and click Calibrate.

Other PAR sensor: To calibrate the internal PAR sensor using an external PAR sensor which cannot be read by the MINI-PAM-II, proceed as described above but adjust calibration factor manually until the readout of the internal sensor matches that of the calibrated external sensor.

Table 26: Custom PAR Gain

The table compiles the types of gain factors with their application, which are available for the PAR sensor of the 2035-B leaf clip or the PAR sensor of the 2065-M device.

Gain factor	Application
Sunlight	Outdoor natural light. Factor displayed in "Ext. PAR Sensor/Calib.".
Red, green, blue, white	Red, green, blue, white LED light from MINI-PAM-II internal light source or from 2054-L external light source.
Mixed	2054-L external light source. The active factor is dynamically adjusted depending on the color mixture selected.
Custom	User-defined factor.

6.11.4 Ext. PAR/T sensors

The boxes “Ext. PAR Sensor” and “Ext. Temp. Sensor” (Fig. 94) display the factory-established calibration factors of the 2030-B clip or the 2065-M Mini Quantum/Temp.-Sensor, depending on which device is connected. In box “Ext. PAR Sensor”, the “Channel 2” calibration applies to a second PAR sensor connected to the 2035-B leaf clip. Calibration factors must be manually entered when a PAR sensor does not support automatic installation of its calibration data.

6.11.5 PAM Trim Val.

Trim values have been set at the factory so that the MINI-PAM-II meets its specifications.

Table 27: PAM Trim Values

The table compiles the types of gain factors with their application, which are available for the PAR sensor of the 2035-B leaf clip or the PAR sensor of the 2065-M device.

Type	Application
Fm-Corr.	Factor to compensate a signal decrease during a saturation pulse. The factor is determined with a Walz fluorescence standard foil. Signal compensation is obsolete (value=0) for the latest generation of LEDs.
Fm-Corr.2	Second factor to compensate a signal decrease during a saturation pulse. Two factors are required when a bi-phasic signal decrease occurs. Signal compensation is obsolete (value=0) for the latest generation of LEDs.
Meas. Light Act. Light SAT Light FR Light	Factors adjusting the intensities of measuring light/actinic light/saturation pulse light/far-red light to meet the respective specification.

6.11.6 Internal Temperature Sensor

Does not apply to the MINI-PAM-II device.

7 Hints & Troubleshooting

7.1 Instrument Settings

Instrument settings were adjusted at the factory for optimum performance of the MINI-PAM-II. For example, LED currents have been adjusted to meet target PAR values for the measuring setup of the 60° distance clip or the 2035-B leaf clip (Fig. 6 and Fig. 13). For a different geometry, the internal PAR sensor requires recalibration to correctly measure internal actinic light (see Fig. 69, page 96). Estimation of PAR according to Fig. 7 (page 20) is rather inaccurate.

Also, the fluorescence offset (F-Offset) of your system has been measured and was saved on the MINI-PAM-II memory. That means that the MINI-PAM-II should show a fluorescence value close to zero with the fiber fully inserted and in the absence of a sample. If this signal deviates clearly from zero, newly adjust F-Offset (see Fig. 45, page 69).

7.2 Default settings

For fluorescence measurements with most healthy green leaves, default settings for measuring light and saturation pulses are well suited. Some samples, however, require special settings. The following sections will provide some hints to adjust settings properly.

7.3 F₀ Fluorescence

Usually, measuring light intensity is adjusted to reach F₀ fluorescence levels around 500 mV (for a definition of F₀ see Chapter 8, page 165). Theoretically, the F₀ should stay below 640 mV. The latter upper value is derived from the assumption that the maximum F_V/F_M of any sample is 0.84 und from the fact that signal saturation occurs at 4000 mV (see equation below, Table 28). If required, measuring light can be adjusted (Fig. 45, page 69).

Table 28: Maximum F₀ of a Dark-acclimated Leaf

$\left(\frac{F_V}{F_M}\right)_{Max} = \frac{(F_M)_{Max} - (F_0)_{Max}}{(F_M)_{Max}}$	(F _M) _{MAX} , maximum possible F _M value = 4000.
with $(F_V/F_M)_{Max} = 0.84$	(F ₀) _{MAX} , unknown maximum F ₀ value (the F _M associated with this F ₀ , or with smaller F ₀ , is not saturating).
$(F_0)_{Max} = 640$	(F _V /F _M) _{MAX} , assumed maximum possible PS II photochemical yield.

At low signal levels, signal height can be increased by increasing measuring light intensity. At too high intensities, the measuring light might drive photosynthesis to some degree. Therefore, test if switching on measuring light results in a stable signal or if any signal increase occurs. In the latter case, the effective measuring light intensity must be decreased either by reducing the amplitude of measuring light (measuring light intensity) or by reducing measuring light frequency or both (Section 4.3.1, page 69, and Section 6.10.2, page 151).

7.4 F_M Fluorescence

The F_M and F_M' levels are determined as the maximum of the fluorescence signal induced by a saturation pulse. Factory settings of saturation pulse width and intensity are adjusted to arrive at a plateau with normal green leaves (Fig. 95A). Some samples (for instance, high light grown leaves) do not reach a plateau with standards settings (Fig. 95B). In this case, saturation pulse intensity or/and length should be increased. Also, fluorescence kinetics can reach its maximum clearly before end of the saturation pulse (Fig. 95C). The latter does not result in erroneous F_M or F_M' values because these values correspond to the maximum of fluorescence kinetics. In case of Fig. 95C, saturation pulse intensity or/and length might be decreased.

Some samples, particularly low light grown or senescing plants, exhibit with standard settings somewhat decreased F_V/F_M values but show normal fluorescence kinetics. These samples increase the F_V/F_M with decreasing saturation pulse intensity. Therefore, testing the F_V/F_M at saturation pulse intensities also below and above standard settings is important to optimize your saturation pulse settings.

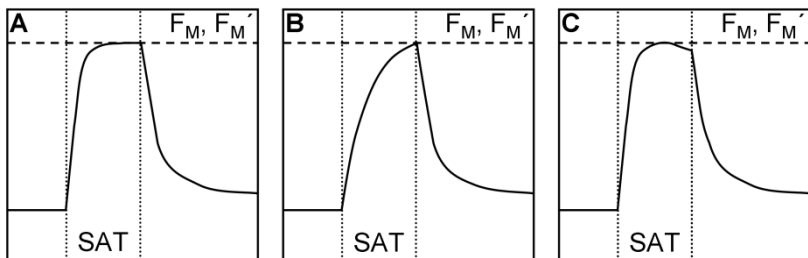


Fig. 95: Fluorescence Kinetics Induced by Saturation Pulse

7.5 Signal Noise

Check if fluctuating light sources (fluorescent tubes, computer screens) affect the signal.

Exclude that automatic scaling of Y axis at low Y(II) has amplified signal noise.

7.6 System Hangs

Disconnect line power and remove batteries. Power on again.

7.7 External PAR Sensor is not Responding

Check the following:

External sensor activated (Fig. 54, page 83. Section 6.11.4, page 159)?

Correct external PAR sensor selected (see Section 4.3.4.1, page 85)?

Sunlight profile selected when measuring external light (Fig. 66, page 93. Section 6.9, page 148)?

When the miniature spectrometer is used, make sure that the spectrometer is in the light measuring mode (not fluorescence or reflectance).

7.8 Intensity of external lamp cannot be adjusted

Repeat lamp calibration.

8 Saturation Pulse Analysis

8.1 Pulse-amplitude Modulated (PAM) Fluorescence

The PAM principle is illustrated by Fig. 96. The top part shows the total fluorescence of a sample. μ s-measuring flashes are given throughout the experiment starting with “Pulse on”. These flashes cause the spikes in the fluorescence trace. From left to right, an external effect induces a “False Signal” of continuous fluorescence in the darkened sample. Then the sample is exposed to a period of actinic illumination (“Actinic on” and “Actinic off”), and, finally, the sample is kept in the dark again.

During actinic illumination, an effect of stray light on the fluorescence signal is additionally assumed. The fluorescence level at onset of stray light plus actinic light is denoted “Actinic F_0 ”. The further increase of continuous fluorescence during illumination is denoted “Actinic F_V ”, where the V stands for variable fluorescence. The “Actinic F_V ” reflects changes of the fluorescence yield in the sample because stray light and actinic light are constant during the illumination period.

In Fig. 96, not only continuous fluorescence varies but also the amplitude of fluorescence spikes. PAM fluorometers ignore the changes of continuous fluorescence and measure only the amplitude of fluorescence spikes. This is achieved by subtracting the fluorescence level just before the μ s-measuring flash from the fluorescence level at the μ s-measuring flash. In Fig. 96, the PAM fluorescence amplitude during the initial dark phase is denoted “Pulsed F_0 ”, and the maximum variable fluorescence at the end of actinic illumination is denoted “Pulsed F_V ”.

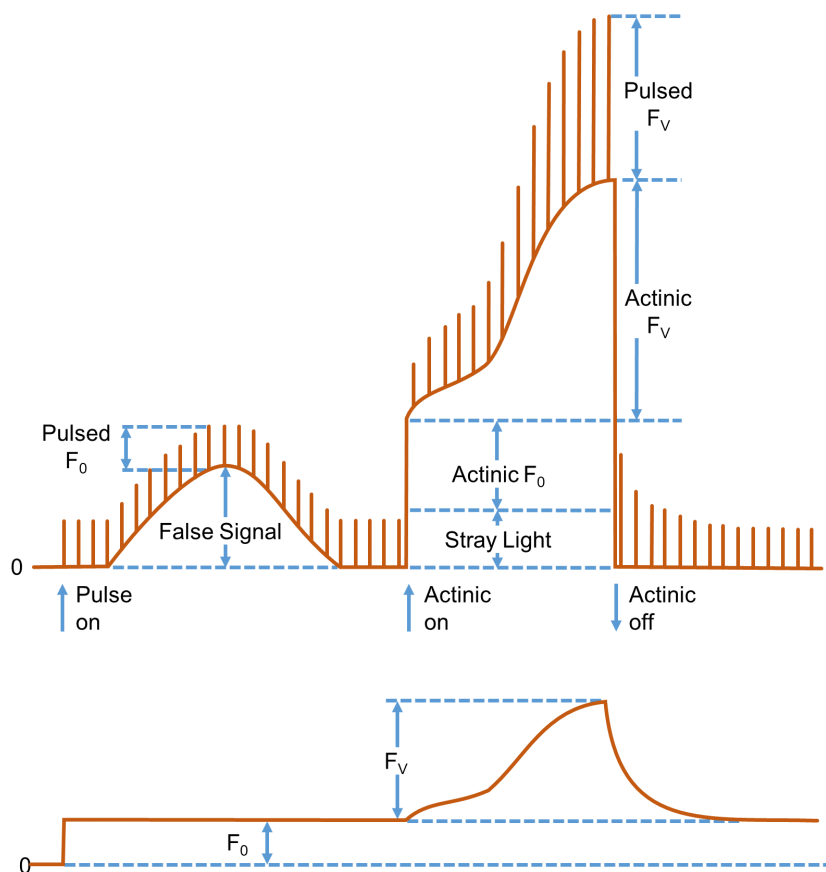


Fig. 96: Illustration of the PAM measurement principle

Figure redrawn from Dr. Ulrich Schreiber, Manual for PAM-101.

Because the μ s-measuring flashes have constant amplitude, the varying amplitudes of fluorescence spikes is a measure of how efficient excitation light is converted into fluorescence. In other words, PAM fluorescence is proportional to the fluorescence yield.

The lower trace in Fig. 96 outlines the PAM fluorescence trace. Obviously, PAM fluorescence irons out the “False Signal” of total

fluorescence at the beginning of the experiment, and also the fluorescence jumps when actinic light is switched on and off. The course of continuous fluorescence within the range “Actinic F_v ” resembles the corresponding trace of PAM fluorescence, because both measuring light and actinic illumination are constant.

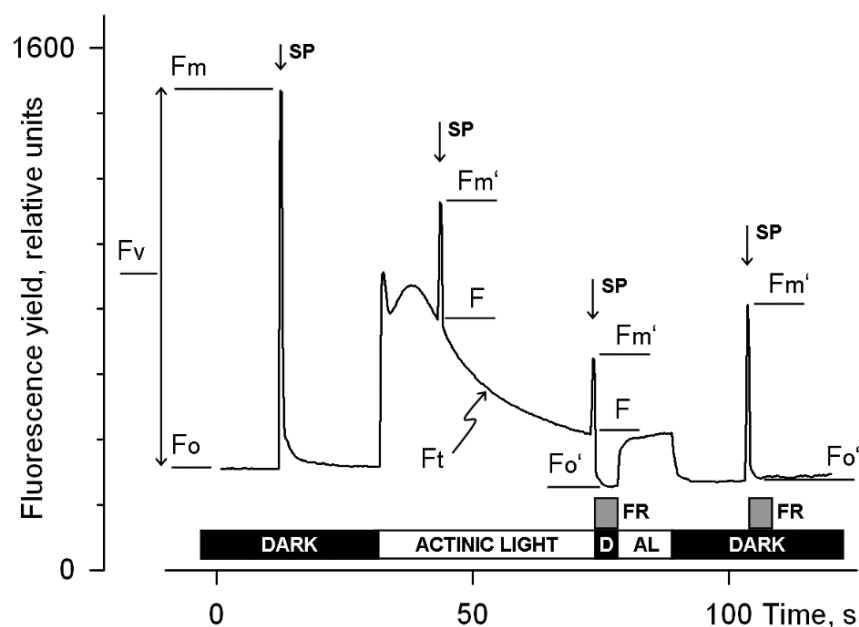


Fig. 97: Fluorescence Levels of Saturation Pulse Analysis

Y-axis (Fluorescence yield) corresponds to PAM fluorescence, see 8.1. AL, Actinic Light; D, dark; SP, Saturation Pulse; F_t , continuously recorded PAM fluorescence; FR, far-red illumination.

8.2 Saturation Pulse Analysis

The five principal levels of PAM fluorescence which are used for saturation pulse analysis are shown in Fig. 97. Two of these levels (F_0 and F_M) must be measured with the dark-acclimated sample. The three other levels (F_0' , F , and F_M') are measured with the actinic light-exposed sample or in a dark period following this light treatment. Some parameters of saturation pulse analysis require fluorescence measurement of the same sample in both the dark-acclimated and light-exposed state (Table 29, page 174).

Because PAM fluorescence is excited by μs pulses of constant amplitude, variations between fluorescence levels are usually interpreted as variation in chlorophyll fluorescence yield. This applies for variations between different types of fluorescence levels (e.g. between F_0 and F_M) and for variations of the same type of fluorescence level (e.g. the change of F_M' during a fluorescence induction curve).

Measurements with Dark-Acclimated Samples

- F_0** Minimum fluorescence level excited by very low intensity of measuring light to keep photosystem II reaction centers open.
- F_M** Maximum fluorescence level elicited by a pulse of saturating light (Saturation Pulse) which closes all photosystem II reaction centers.

Measurements with Illuminated Samples

- F_0'** Minimum fluorescence level of illuminated sample. The F_0' is lowered relative to F_0 by non-photochemical quenching. The measuring routine for F_0' (see 4.3.1.6, page 72) determines the F_0' level during a dark interval following a Saturation Pulse. In this dark interval, far-red

light is applied which selectively drives photosystem I. As a consequence, electrons are removed from the intersystem electron transport chain and opening of photosystem II reaction centers is efficiently accelerated

If the F_0' Mode is switched off, the F_0' will be calculated according to Oxborough and Baker:

$$F_0' = \frac{1}{\frac{1}{F_0} - \frac{1}{F_M} + \frac{1}{F_M'}}$$

Oxborough K, Baker NR (1997) Resolving chlorophyll a fluorescence images of photosynthetic efficiency into photochemical and non-photochemical components - calculation of qP and F_v'/F_m' without measuring F_0' . *Photosynth Res* 54 135-142. <https://doi.org/10.1023/A:1005936823310>

In the Record table, the calculated value F_0' is preceded by a tilde sign (~).

- F_M'** Maximum fluorescence level of the illuminated sample. The F_M' is induced by a Saturation Pulse which temporarily closes all photosystem II reactions centers. F_M' is decreased relative to F_M by non-photochemical quenching.
- F** The F corresponds to the momentary fluorescence level (F_t) of an illuminated sample shortly before application of a Saturation Pulse.

To quantify photochemical use and non-photochemical losses of absorbed light energy, fluorescence ratio expressions have been derived which use two or more of the five relative fluorescence yields introduced above. Table 29 (page 174) compiles the fluorescence ratio parameters available in the DualPAM software. Below, these parameters will be explained briefly.

F_V/F_M and $Y(II)$ Maximum and effective photochemical quantum yields of photosystem II

The F_V/F_M and $Y(II)$ estimate the fraction of absorbed quanta used for photosystem II photochemistry. F_V/F_M corresponds to the maximum photochemical yield of photosystem II, $Y(II)$ is the effective photochemical yield of photosystem II. Measurements of F_V/F_M require that samples are acclimated to darkness or dim light so that all reactions centers are in the open state and non-photochemical dissipation of excitation energy is minimal.

In algae and cyanobacteria, however, the dark-acclimated state often is not showing maximal photosystem II quantum yield, as the photosystem II acceptor pool may be reduced in the dark by stromal reductants and, consequently, the so-called state 2 is formed exhibiting low photosystem II quantum yield. In this case, preillumination with moderate far-red light should precede determinations of F_0 and F_M .

The $Y(II)$ value estimates the photochemical use of excitation energy in the light. It is lowered with respect to F_V/F_M by partial closure of photosystem II centers and various types of non-photochemical energy losses induced by illumination.

 q_P and q_L Coefficients of photochemical fluorescence quenching

Both parameters estimate the fraction of open photosystem II reaction centers. The q_P is based on the concept of separated photosystem II antenna units (puddle model), whereas the q_L assumes interconnected photosystem II antenna units (lake model) which was assumed to be present in leaves (*cf. Kramer et al.*, 2004). Determinations of q_P and q_L do not require fluorescence measurements with the dark-acclimated sample, except the F_0'

mode is switched off and F_0' is calculated according to Oxborough and Baker (1997).

q_N and NPQ Parameters of non-photochemical quenching

Both parameters are associated with non-photochemical quenching of excitation energy, mainly involving a low thylakoid lumen pH- and a zeaxanthin-dependent quenching mechanism. The q_N and the NPQ parameters require fluorescence measurements with the sample in the dark-acclimated and in the light-exposed states (cf. Table 29, page 174).

Calculation of NPQ (or SV_N ; Gilmore and Yamamoto, 1991) corresponds to the Stern-Volmer equation for fluorescence quenching which predicts proportionality between fluorescence quenching (NPQ) and the concentration of fluorescence-quenching centers in the photosynthetic antennae (e.g. zeaxanthin).

$Y(NO)$, $Y(NPQ)$ and $Y(II)$ Complementary photosystem II yields

Genty *et al.* (1996) and Kramer *et al.* 2004 have presented expressions describing the partitioning of absorbed excitation energy in photosystem II between three fundamental pathways the sum of which adds up to one:

$Y(NO)$ non-regulated losses of excitation energy including heat dissipation and fluorescence emission,

$Y(NPQ)$ regulated energy losses of excitation energy by heat dissipation involving ΔpH - and zeaxanthin-dependent mechanisms, and

$Y(II)$ use of excitation energy for charge separation.

This concept of "complementary photosystem II quantum yields" is useful to analyze the partitioning of absorbed light energy in photosynthetic organisms. For instance, in the presence of strong

light, a much higher $Y(NPQ)$ than $Y(NO)$ indicates that excess excitation energy is safely dissipated at the antenna level and that photosynthetic energy fluxes are well-regulated.

In variance, high values of $Y(NO)$ would signify that excess excitation energy is reaching the reaction centers, resulting in strong reduction of photosystem II acceptors and photodamage, e.g. via formation of reactive oxygen species.

8.3 Relative Electron Transfer Rate (ETR)

Relative electron transfer rates for photosystem II are calculated according to:

$$ETR(II) = PAR \cdot ETR\text{-Factor} \cdot P_{PS2}/P_{PS1+2} \cdot Y(II).$$

The basic idea of the ETR equation is to multiply $Y(II)$, the effective photochemical quantum yield of photosystem II, by an estimate for the photon flux density absorbed by all photosystem II in the sample. The latter estimate is derived from three numbers:

- (1) **PAR** Quantum flux density of photosynthetically active radiation (PAR) impinging on the sample.
- (2) **ETR-Factor** Sample absorbance (= 1 – transmittance)

The ETR-Factor describes the fraction of incident photons absorbed by the sample. The most frequently used default value for green leaves is 0.84 meaning that 84% of incoming light is absorbed. The ETR-Factor can be lower in bleached leaves or leaves containing considerable amounts of non-photosynthetic pigments like anthocyanins.

- (3) P_{PS2}/P_{PS1+2} Relative distribution of absorbed PAR to photosystem II

The default P_{PS2}/P_{PS1+2} is 0.5 which assumes the photosystem II contributes 50% to total sample absorbance. The P_{PS2}/P_{PS1+2} may deviate from the idealized factor of 0.5 depending on wavelength of light and acclimation status of the sample.

8.3.1 Reviews on Saturation Pulse Analysis of Photosystem II

Baker NR (2008) Chlorophyll fluorescence: a probe of photosynthesis in vivo. *Annu Rev Plant Biol* 59: 89–113. <https://doi.org/10.1146/annurev.arplant.59.032607.092759>

Kalaji HM, Schansker G, Ladle RJ, Goltsev V, Bosa K, Allakhverdiev SI, Brestic M, Bussotti F, Calatayud A, Dąbrowski P, Elsheery NI, Ferroni L, Guidi L, Hogewoning SW, Jajoo A, Misra AN, Nebauer SG, Pancaldi S, Penella C, Poli DB, Pollastrini M, Romanowska-Duda ZB, Rutkowska B, Seródio J, Suresh K, Szulc W, Tambussi E, Yannericari M, Zivcak M (2014) Frequently asked questions about *in vivo* chlorophyll fluorescence: practical issues. *Photosynth Res* 122: 121–158. <https://doi.org/10.1007/s11120-014-0024-6>

Krause GH, Weis E (1991) Chlorophyll fluorescence and photosynthesis: the basics. *Annu Rev Plant Physiol Plant Mol Biol* 42: 313–349. <https://doi.org/10.1146/annurev.pp.42.060191.001525>

Maxwell K, Johnson GN (2000) Chlorophyll fluorescence – a practical guide. *J Exp Bot* 51: 659–668 <https://doi.org/10.1093/jexbot/51.345.659>

Murchie EH, Lawson T (2013) Chlorophyll fluorescence analysis: a guide to good practice and understanding some new applications. *J Exp Bot* 64: 3983–3998. <https://doi.org/10.1093/jxb/ert208>

Schreiber U (2004) Pulse-amplitude-modulation (PAM) fluorometry and saturation pulse method: an overview. In: Papageorgiou GC, Govindjee (eds) *Chlorophyll a Fluorescence: A Signature of Photosynthesis*. Springer, Dordrecht, pp 279–319. https://doi.org/10.1007/978-1-4020-3218-9_11

Table 29: Fluorescence Ratio Parameters

Source	Equation	Sample State	Range [Theory] [Experiment]
Maximum photochemical quantum yield of PS II (Kitajima and Butler, 1975)	$\frac{F_V}{F_M} = \frac{F_M - F_0}{F_M}$	Dark	[0, 1] [0, ~0.84]
Effective photochemical quantum yield of PS II (Genty <i>et al.</i> , 1989)	$Y(II) = \frac{F'_M - F}{F'_M}$	Light	[0, 1] [0, ~0.84]
Quantum yield of light-induced (Δ pH- and zeaxanthin-dependent) non-photochemical fluorescence quenching (Genty <i>et al.</i> 1996, Kramer <i>et al.</i> 2004)*	$Y(NPQ) = \frac{F}{F'_M} - \frac{F}{F_M}$	Dark and Light	[0, 1] [0, ~0.9]
Quantum yield of non-regulated heat dissipation and fluorescence emission: this type of energy loss does not involve the action of a trans-thylakoid Δ pH and zeaxanthin (Genty <i>et al.</i> 1996, Kramer <i>et al.</i> 2004)*	$Y(NO) = \frac{F}{F_M}$	Dark and Light	[0, 1] [0, ~0.9]
Stern-Volmer type non-photochemical fluorescence quenching (Bilger and Björkman, 1990; Gilmore and Yamamoto, 1991))	$NPQ = \frac{F_M}{F'_M} - 1$	Dark and Light	[0, ∞] [0, ~4]
Coefficient of photochemical fluorescence quenching (Schreiber <i>et al.</i> 1986 as formulated by van Kooten and Snel, 1990)	$q_P = \frac{F'_M - F}{F'_M - F'_0}$	Light. If F'_0 calculated, Dark and Light	[0, 1] [0, 1]
Coefficient of photochemical fluorescence quenching assuming interconnected PS II antennae (Kramer <i>et al.</i> 2004)	$q_L = q_P \cdot \frac{F'_0}{F}$	As q_P .	[0, 1] [0, 1]
Coefficient of non-photochemical fluorescence quenching (Schreiber <i>et al.</i> 1986 as formulated by van Kooten and Snel, 1990)	$q_N = 1 - \frac{F'_M - F'_0}{F_M - F_0}$	Dark and Light	[0, 1] [0, ~0.95]

* Kramer *et al.* (2004) have derived more complex equations for $Y(NO)$ and $Y(NPQ)$. Klughammer and Schreiber (2008) have transformed the equations by Kramer *et al.* (2004) into the simple equations of Genty *et al.* (1996).

Table 30: References Cited in Table 29

- Bilger W, Björkman O (1990) Role of the xanthophyll cycle in photoprotection elucidated by measurements of light-induced absorbance changes, fluorescence and photosynthesis in leaves of *Hedera canariensis*. *Photosynth Res* 25:173-185. <https://doi.org/10.1007/BF00033159>
- Genty B, Briantais J-M, Baker NR (1989) The relationship between the quantum yield of photosynthetic electron transport and quenching of chlorophyll fluorescence. *Biochim Biophys Acta* 990: 87–92. [https://doi.org/10.1016/S0304-4165\(89\)80016-9](https://doi.org/10.1016/S0304-4165(89)80016-9)
- Genty B, Harbinson J, Cailly AL and Rizza F (1996) Fate of excitation at PS II in leaves: the non-photochemical side. Presented at: The Third BBSRC Robert Hill Symposium on Photosynthesis, March 31 to April 3, 1996, University of Sheffield, Department of Molecular Biology and Biotechnology, Western Bank, Sheffield, UK, Abstract P28
- Gilmore AM, Yamamoto HY (1991) Zeaxanthin formation and energy-dependent fluorescence quenching in pea chloroplasts under artificially mediated linear and cyclic electron transport. *Plant Physiol* 96:635-643. <https://doi.org/10.1104/pp.96.2.635>
- Kitajima M, Butler WL (1975) Quenching of chlorophyll fluorescence and primary photochemistry in chloroplasts by dibromothymoquinone. *Biochim Biophys Acta* 376:105-115. [https://doi.org/10.1016/0005-2728\(75\)90209-1](https://doi.org/10.1016/0005-2728(75)90209-1)
- Klughammer C and Schreiber U (2008) Complementary PS II quantum yields calculated from simple fluorescence parameters measured by PAM fluorometry and the Saturation Pulse method. *PAM Application Notes* 1: 27-35 (<https://www.walz.com/files/downloads/pan/PAN078007.pdf>)
- Kramer DM, Johnson G., Kiirats O, Edwards GE (2004) New flux parameters for the determination of Q_A redox state and excitation fluxes. *Photosynth Res* 79: 209-218. <https://doi.org/10.1023/B:PRES.0000015391.99477.0d>
- Schreiber U, Schliwa U, Bilger W (1986) Continuous recording of photochemical and non-photochemical chlorophyll fluorescence quenching with a new type of modulation fluorometer. *Photosynth Res* 10: 51–62. <https://doi.org/10.1007/BF00024185>
- van Kooten O, Snel J (1990) The use of chlorophyll fluorescence nomenclature in plant stress physiology. *Photosynth Res* 25: 147–150. <https://doi.org/10.1007/BF00033156>
-

9 Specifications MINI-PAM-II

9.1 Basic System

9.1.1 Optoelectronic Unit

MINI-PAM-II/B (Blue Version)

Measuring light: Blue (470 nm) LED, standard modulation frequencies 5 to 25 Hz, adjustable in increments of 5 Hz, and 100 Hz, measuring light PAR at standard settings = $0.05 \mu\text{mol m}^{-2} \text{s}^{-1}$. Fluorescence at wavelengths greater than 630 nm is measured

Actinic light: Same blue LED as for measuring light, maximum actinic PAR = $3000 \mu\text{mol m}^{-2} \text{s}^{-1}$, maximum PAR of saturation pulses = $6000 \mu\text{mol m}^{-2} \text{s}^{-1}$ adjustable at increments of $500 \mu\text{mol m}^{-2} \text{s}^{-1}$.

MINI-PAM-II/R (Red Version)

Measuring light: Red (655 nm) LED, modulation frequencies and PAR as described for MINI-PAM-II/B. Fluorescence at wavelengths greater than 700 nm is measured

Actinic light: Same red LED as for measuring light, maximum PAR of actinic light and saturation pulses as described for MINI-PAM-II/B

MINI-PAM-II/B and MINI-PAM-II/R

Far-red light: Peak emission at 735 nm

Signal detection: PIN photodiode protected by long-pass and a short-pass filters

Data memory: Flash memory, 8 MB, providing memory for more than 27,000 saturation pulse analyses

Display: Backlit 160x104 dots (78 x 61 mm) transfective B/W LCD display with resistive touchscreen

Ports: Ports for fiberoptics, USB cable, external light source, 2035-B leaf clip, auxiliaries and 12 V DC power supply


Power supply: 6 AA (Mignon) rechargeable batteries (Eneloop 1.2 V/2 Ah), providing power for up to 1000 yield measurements; 6 spare batteries, automatic power/off, battery charger (100 to 240 V AC, 50-60 Hz, 0.35 A) for 1 to 8 AA/AAA NI-MH/NI-CD batteries, 12 V 5,5 A power supply MINI PAM-II/N

Operating temperature: -5 to +45 °C, non-condensing

Dimensions: 17.2 cm x 11.2 cm x 7.6 cm (L x W x H)

Weight: 1.5 kg (incl. battery)

Table 31: Pin Assignments. View from Outside.

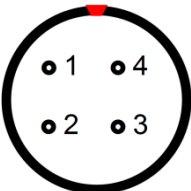	LEAF CLIP/AUX 1/AUX 2
1	GND
2	Serial Clock
3	Serial Data
4	+ 5 V Output

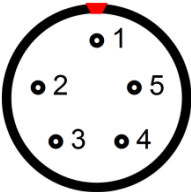
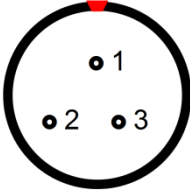
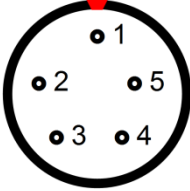
	
1	SYNC
2	GND
3	SYNC
4	RS 485 A
5	RS 485 B
6	Voltage output of current power supply (line or battery)

Table 31: Pin Assignments. View from Outside.

	<p>EXT. DC</p> <table> <tr> <td>1</td><td>Not used</td></tr> <tr> <td>2</td><td>GND</td></tr> <tr> <td>3</td><td>+ 12 V Input</td></tr> </table>	1	Not used	2	GND	3	+ 12 V Input				
1	Not used										
2	GND										
3	+ 12 V Input										
	<p>COMP 2/ COMP 1</p> <table> <tr> <td>1</td><td>+ 5 V Output (= 4 of LEAF CLIP)</td></tr> <tr> <td>2</td><td>Voltage output of current power supply (= 5 of SYNC)</td></tr> <tr> <td>3</td><td>RS 232 TxD</td></tr> <tr> <td>4</td><td>RS 232 RxD</td></tr> <tr> <td>5</td><td>GND</td></tr> </table>	1	+ 5 V Output (= 4 of LEAF CLIP)	2	Voltage output of current power supply (= 5 of SYNC)	3	RS 232 TxD	4	RS 232 RxD	5	GND
1	+ 5 V Output (= 4 of LEAF CLIP)										
2	Voltage output of current power supply (= 5 of SYNC)										
3	RS 232 TxD										
4	RS 232 RxD										
5	GND										

9.1.2 Fiberoptics MINI-PAM/F

Design: Randomized 70 μm glass fibers forming single plastic shielded bundle with stainless steel adapter ends

Dimensions: Active diameter 5.5 mm, outer diameter 8 mm, length 100 cm

Weight: 180 g

9.1.3 Power Supply MINI-PAM-II/N

Input: 100 to 240 V AC, 50 to 60 Hz

Output: 12 V DC, 5.5 A

Operating temperature: -5 to +45 °C, (non-condensing)

Dimensions: 13 cm x 5.5 cm x 3 cm (L x W x H)

Weight: 350 g including cables

9.1.4 Battery Charger 000190101101

Input: 100 to 240 V AC, 50 to 60 Hz

Output: 12 V DC, 1.0 A

Operating temperature: -5 to +45 °C, (non-condensing)

Dimensions: 17.5 cm x 10.5 cm x 3 cm (L x W x H)

Weight: 300 g including cable

9.1.5 Distance Clip 60° 2010-A

Design: Metal clip with fiber holder and 11 mm diameter sample hole: 5.5 cm x 1.4 cm (L x W)

9.1.6 Complementary Items

Sloped Plexiglas rack for convenient desktop operation. Stylus for touchscreen. Carrying strap for optoelectronic unit

9.1.7 Software WinControl-3

Program: WinControl-3 System Control and Data Acquisition Program (Windows XP/Vista, Windows 7+8 32-bit and 64-bit) for operation of measuring system via PC, data acquisition and analysis

Saturation Pulse Analysis: Measured: F_t , F_0 , F_M , F , F_0' (also calculated), F_M' . PAR, leaf temperature and relative humidity using 2035-B Leaf-Clip Holder. Calculated: F_0' (also measured), F_V/F_M

and $Y(II)$ (maximum and effective photochemical yield of PS II, respectively), q_L , q_P , q_N , NPQ, $Y(NPQ)$, $Y(NO)$ and ETR (electron transport rate)

Fitting Routines: Two routines for determination of the cardinal points α , I_k and ETR_{max} of light curves

Programmed Features: Automatic determination of signal offset for all light intensities and all gain levels. Automatic calibration of internal PAR sensor against an external PAR sensor connected to the MINI-PAM-II

Communication Protocol: USB

Computer Requirements: Processor, 0.8 GHz. RAM, 512 MB. Screen resolution, 1024 x 600 pixels. Interface, USB 2.0/3.0.

9.1.8 Transport Case MINI-PAM/T

Design: Aluminum case with custom foam packing

Dimensions: 50 cm x 34 cm x 20 cm (L x W x H)

Weight: 3.8 kg

9.2 Accessories

9.2.1 2035-B Leaf-Clip Holder

Design: Consisting of a port to position the MINI-PAM/F Fiberoptics and a clip to hold the sample. A circular 1 cm diameter hole of the upper clip part defines the measuring area. Standard distance between fiberoptics tip and measuring area is 8 mm. The port aligns the fiberoptics at an angle of 60° relative to the measuring

plane. A mini quantum sensor is positioned on the sample level by a movable Perspex arm, a leaf temperature sensor is mounted below the sample within an up and down movable Plexiglas tube, and a humidity sensor is mounted at 3 cm distance from the sample. The electronics of the 2035-B unit stores calibration factors of sensors. Saturation pulses can be released by remote trigger button. An input socket for an additional light sensor is provided

Micro quantum sensor: LS-C sensor for selective PAR measurement, range 0 to 7000 $\mu\text{mol m}^{-2} \text{s}^{-1}$, cosine corrected for light incident at an angle between -30° to $+30^\circ$ from surface normal, internal preamplifier

Thermocouple: Ni-CrNi, wire diameter 0.1 mm, -20 to $+60^\circ\text{C}$

Humidity sensor: Humidity and temperature sensing integrated circuit, 0 – 100% relative humidity

Power supply: MINI-PAM-II leaf clip socket (5 V/10 mA)

Cable length: 180 cm

Dimensions: 17 cm x 5.7 cm (max.) x 8 cm (max.) (L x W x H)

Weight: 250 g (excluding cable)

9.2.2 **Fiberoptics Adapter 90° 2030-B90**

Accessory to Leaf Clip Holder 2035-B. Positions MINI-PAM-II fiberoptics at a 90° angle relative to the leaf surface

9.2.3 **External LED Source 2054-L**

Accessory to Leaf Clip Holder 2035-B to illuminate the sample plane at an angle of 45° with red, green, blue or white light, or

mixtures of these four light qualities. The intensity of each light quality is adjusted individually and the maximum PAR per light quality is $1500 \mu\text{mol m}^{-2} \text{s}^{-1}$.

Design: LED holder for 4 RGBW LEDs made of anodized aluminum with cooling ventilator on the back side. The device can be fastened by two 7 cm long aluminum rods whose ends fit in drilled holes of the 2035-B clip

Dimensions: 9 cm (max.) x 4.5 cm x 12 cm (max.) (L x W x H)

Weight: 210 g without cable

9.2.4 Dark Leaf Clip DLC-8

Design: Clip made of aluminum with felt contact areas and sliding shutter

Dimensions: 6.5 cm x 2 cm (max.) x 1.5 cm (max.) (L x W x H)

Weight: 3.6 g

9.2.5 Arabidopsis Leaf Clip 2060-B

Design: Aluminum clip with 3.2 mm diameter viewing area designed to position small leaves below the fiberoptics of the MINI-PAM-II, prepared to accommodate PAR and temperature sensors of the Mini Quantum/Temp.-Sensor 2065-M

Dimensions: 7.6 cm x 3.0 cm (max.) x 5.2 cm (max.) (L x W x H)

Weight: 55 g

9.2.6 Conversion Kit for 2030-B/2035-B Leaf Clips 2035-B/RLC

Design: Duct and mounting bar, made of Plexiglas, replacing the standard 60° fiberoptics duct of the 2030-B clip (PAM-2500) or the 2035-B clip (MINI-PAM-II). The duct positions the fiberoptics 2010-F (2030-B clip) or fiberoptics MINI-PAM/F (2035-B clip) perpendicularly to the sample surface. With lateral set screw. Including two circular light-shielding elements made from cellular rubber and a 3 cm x 2 cm (L x W) black plastic base plate with cellular rubber rim

Dimensions, max: 7 cm x 2.6 cm x 2.7 cm (L x W x H)

Weight: 12 g

9.2.7 Fiberoptics Holder for Surfaces 2060-A

Aluminum plate (6.0 x 3.3 cm max.) with 11 mm diameter circular hole (measuring area) and aluminum port to position fiber at an angle of 60° relative to the aluminum plate. With port for temperature sensor of 2065-M unit to measure surface temperature and thread to mount the PAR sensor of the 2065-M unit. Connected to a 10 x 0.8 cm (L x Ø) steal rod with two lateral aluminum supports (12 cm x 1 cm x 1 cm, L x W x H) which are lockable by knurled screws

Dimensions (without aluminum supports): 15 cm x 3.3 cm x 2.5 cm (L x W x H)

Weight: 125 g

9.2.8 Mini Quantum/Temp.-Sensor 2065-M

Mini quantum and temperature sensors connected by 30 cm cables to an electronic unit for signal amplification, digitization and storage of calibration factors. A 10 cm steel rod can be laterally screwed-on to the electronic unit. A 110 cm cable connects the 2065-M unit with the fluorometer

Mini quantum sensor: LS-C sensor for selective PAR measurement, range 0 to 7000 $\mu\text{mol m}^{-2} \text{s}^{-1}$, cosine corrected for light incident at angles between -30 to +30 from surface normal

Thermocouple: Ni-CrNi, wire diameter 0.1 mm, -20 to +60 C

Dimensions of electronic unit: 15 cm x 3.3 cm x 2.5 cm (L x W x H)

Weight: 125 g (excluding cable)

9.2.9 Miniature Spectrometer MINI-SPEC/MP

Design: POM tube, at one side, port for light detection, port for fluorescence excitation by blue (452 nm max) or green (525 nm max) LEDs, and port for white light from a tungsten lamp for reflectance measurements; at the opposite side, 4-pole underwater socket.

Spectrometer: Hamamatsu micro-spectrometer, spectral range: 400 to 800 nm, spectral resolution: between 8 and 10 nm.

Maximum PAR: 4000 $\mu\text{mol m}^{-2} \text{s}^{-1}$ for illumination having spectral characteristics similar to sunlight

Dimension: 3.25 cm diameter, 175 cm length max

Weight: 135 g

9.2.10 Flat Entrance Optics SPEC/P

Design: Hard-anodized aluminum rod of 10 mm diameter and 50 mm length, at one end with lateral light entrance through a 5 mm diameter diffusor and the opposite end inserted in a mounting plate (diameter 33 mm, height 5 mm). Aluminum rod with internal light guide

9.2.10.1 Fluorescence and Reflectance Optics SPEC/R

Design: Spectrometer cap consisting of POM: maximum diameter 35 mm, height 13 mm, weight 16 g. With central 5 mm x 16 mm groove which accommodates at one end a Perspex light guide for fluorescence excitation by blue or green light, and at the other end a Perspex light guide for white light for reflectance measurements. With 3 mm diameter central drilled hole as light channel to the detector window of the spectrometer. The sample is fixed between the cap part and another disk (maximum diameter 40 mm, height 10 mm, weight 8 g). The sample side of cap and disk is padded with foam rubber, both parts have magnets build in to attract each other and, thus, hold the sample. Including a 10 mm thick Zenith Polymer reflectance standard

9.2.11 PAR Calibration Block 000160101439

Design: POM block with drill hole for entrance optics of the Miniature Spectrometer. Oriented at an angle of 60° and 90° relative to the spectrometer port are drill holes for the Fiberoptics MINI-PAM-F

Dimensions: 4.15 cm x 2 cm x 5 cm (L x W x H)

Weight: 40 g

9.2.12 Suspension Cuvette KS-2500

Cuvette: Round stainless-steel cuvette (7.5 mm wide, 9.0 mm deep) with top window adapter for connecting the fiberoptics of the MINI-PAM-II; embedded in POM body with injection port for microliter syringes and hose nozzles for connecting an external flow-through water bath (not included). Including three 6.0 x 1.5 mm magnetic stir bars

Dimensions: 11.5 cm x 9 cm x 8.5 cm (L x W x H, including ports)

Weight: 252 g

9.2.13 Magnetic Stirrer with Fiberoptics Holder MKS-2500

Magnetic stirrer: To drive the magnetic stir bar in the Suspension Cuvette KS-2500; with Plexiglas ring for centering the cuvette and miniature stand to fix the fiberoptics on top of the cuvette

9.2.14 Compact Tripod ST-2101A

Stand for 2035-B Leaf-Clip Holder.

Height: Adjustable between 24 cm and 87 cm

Weight: 400 g

9.2.15 Barcode Scanner BCS-9590

Design: Single-line handheld laser scanner with trigger button and 1 m, partially coiled cord; to be connected to the Photosynthesis Yield Analyzer MINI-PAM-II. Bar codes are stored together with fluorescence data

Dimensions: 9 cm (max.) x 6 cm (max.) x 16 cm (max.) (L x W x H)

Weight: 335 g

9.2.16 MINI-PAM/F1 Miniature Fiberoptics

Active diameter 2 mm, length 1.5 m. Including adapter for attachment to Leaf-Clip Holder 2035-B.

Subject to change without prior notice

10 Guarantee

10.1 Manufacturer's Guarantee

Under this Manufacturer's Guarantee ("Guarantee"), subject to the Conditions and Instructions below, Heinz Walz GmbH, Germany ("Manufacturer"), guarantees (§443 BGB) to the end customer and user ("Customer") that all products supplied by it shall substantially conform in material respects to the Specifications for 24 months from the delivery date (date on invoice). In this Guarantee, "Specifications" means the product's features (as may be amended by Manufacturer from time to time), which are set out under the headings "specifications" and/or "technical specifications" within the product's respective brochure, data sheet, or respective tab on the Manufacturer's website for such product, and which may be included with the documents for the product when delivered. In case of an eligible guarantee claim, this Guarantee entitles the Customer to repair or replacement, at the Manufacturer's option, and this Guarantee does not include any other rights or remedies.

10.2 Conditions

This Guarantee shall not apply to:

- Any defects or damage directly or indirectly caused by or resulting from the use of unauthorized replacement parts and/or service performed by unauthorized personnel.
- Any product supplied by the Heinz Walz GmbH, Germany which has been subjected to misuse, abuse, abnormal use, negligence, alteration or accident.
- Damage caused from improper packaging during shipment or any acts of God.

- Batteries, cables, calibrations, fiberoptics, fuses, gas filters, lamps (halogen, LED), thermocouples, and underwater cables.
- Defects that could reasonably have been detected upon inspection of the product when received by the Customer and not promptly noticed within ten (10) days to Heinz Walz GmbH.
- Submersible parts of the DIVING-PAM or the underwater version of the MONITORING-PAM have been tested to be watertight down to the maximum operating depth indicated in the respective manual. Guarantee shall not apply for diving depths exceeding the maximum operating depth. Further, guarantee shall not apply for damage resulting from improper operation of devices, in particular, the failure to properly seal ports or sockets.

10.3 Instructions

- To obtain guarantee service, please follow the instructions below:
- The Walz Service Information Form available at https://www.walz.com/support/repair_service.html must be completed and returned to Heinz Walz GmbH, Germany.
- The product must be returned to Heinz Walz GmbH, Germany, within 30 days after Heinz Walz GmbH, Germany has received written notice of the defect. Postage, insurance, and/or shipping costs incurred in returning equipment for guarantee service are at customer expense. Duty and taxes are covered by Walz.
- All products being returned for guarantee service must be carefully packed and sent freight prepaid.

-
- Heinz Walz GmbH, Germany is not responsible or liable for missing components or damage to the unit caused by handling during shipping. All claims or damage should be directed to the shipping carrier.

10.4 Applicable law

- This Guarantee is governed by German law. Place of jurisdiction is Bamberg, Germany.

11 Index

2

2030-B90 90° Fiberoptics Adapter .27, 33

2035-B Leaf-Clip External Light Source29

2035-B Leaf-Clip Holder26, 181

2035-B Leaf-Clip Signal LED.....29

2035-B/RLC Conversion Kit.....31, 184

2054-L External LED Source 8, 29, 30, 76

2060-A & 2065-M, Surface Holder with Quantum/Temp.-Sensor34

2060-A Fiberoptics Holder34

2060-B & 2065-M Arabidopsis Clip with Quantum/Temp.-Sensor.....33

2065-M Mini Quantum/Temp.-Sensor35

9

90° Fiberoptics Adapter 2030-B90 .27, 33

A

Abbreviations for Settings.....138

Accessories115, 181

Accessories Menu115

Act. Int.125

Actinic + Yield 54, 55, 80, 125, 153

Actinic + Yield Settings 80

Actinic Factor 74

Actinic Intensity 74

Actinic Light..... 74, 177

Actinic Light List 61, 152

ADC Error..... 149

Address 109

Adjust F-Offset 73

Alpha 130

Arabidopsis Leaf Clip 2060-B..... 183

Asterisk 122

Auto Off 98

Auto Range 135

AUX 1 15

Average..... 135

B

Back Light 98

Barcode Scanner 38, 188

Basic Data..... 46, 47

Basic System 177

Batch Window Overview 145

Battery..... 10, 19, 126, 178

Battery Charger..... 180

BCS-9590 Barcode Scanner..... 39

Beeper..... 99, 155

Bilger	175	Connection of MINI-SPEC/MP	21
Blue Version	177	Control Window	146, 148
Butler	175	Conversion Kit	31, 184
Buttons	118	Cosine Response	28
C		Cursor Mode	46, 54
CAL Command	64	Custom PAR Gain	158
Calibr. Dark	134	Custom PAR Gain	157
Calibrate	159	Cuvette	37
Calibrate Light List	63	D	
Calibration	152	Damping	72, 153
Calibration Factor	66	Dark Current	134
Cardinal Parameters	130	Dark Leaf Clip DLC-8	183
Carrying Belt	14	Date	99
Carrying Belt Fastening	16	Default Settings	103
Change Address	109	Delete Data	118
Change Marker	47	Delete Record	121
Channel	109	Device Channel Configuration	109, 110
Chart Area	119	Distance Clip 60° 2010-A	19, 180
Chart Tools	120	DLC-8 Dark Leaf Clip	36
Chart Window	114	Download Memory	141
Chart Window	114	E	
Circular Buffer	139	Electron Transport Rate	172
Clear Memory	141	Emission Spectra	18, 24
Clk. Time	125	Emission Spectra of Miniature Spectrometer Light Sources	25
Clock	78, 154	Emission Spectra of MINI-PAM II LEDs	18
Clock Item	79	Emission spectra of the 2054-L External LED Source	30
COMP 1	15		
Compact Tripod ST 2101A	27		
Compact Tripod ST-2101A	187		
Components And Setup	13		

Emission Spectrum.....	52	F ₀ ' Measurement.....	168
EPROM.....	109	Far-Red Light	125
Equations.....	174	Far-Red Settings	76
ETR.....	122	Fiberoptics.....	8, 17, 179
ETR Factor	72, 153	Fiberoptics Adapter 90°	27
ETR(II)	172	Fiberoptics Adapter 90° 2030-B90	182
ETRm.....	130	Fiberoptics Adapter 90° 2030-B90 - Mounting	32
Excel.....	115	Fiberoptics Holder for Surfaces 2060- A.....	184
Export	115	File Menu	115
Export Chart Record	115	FireStingO2 Oxymeter	86, 149
Export Format for Reports	116	Firmware	100
Export Record.....	119	Firmware Info	101, 102
Export Regression Data.....	131	First Three Submenus in Detail.....	68
Export Selection.....	119	Flat Entrance Optics SPEC/P	186
Ext. Panel Mode	76, 147	Float	118
Extent of Delivery.....	13	Fluorescence Detection	177
External Light Source	147	Fluorescence Emission	23
External Light Source	29, 62	Fluorescence Kinetics Induced by Saturation Pulse.....	163
External Light Source	182	Fluorescence Levels of Saturation Pulse Analysis.....	167
External PAR	95	Fluorescence Ratio Parameters....	174
External PAR Sensor.....	85, 148, 159	Fluorescence Reviews.....	173
External PAR Sensor Is Not Responding.....	164	Fluorescence Spectra	24, 134
External Temperature Sensor.....	159	F _M Definition	168
External Temperature Sensor Settings	94	F _M Determination.....	163
F		F _M Kinetics	163
F ₀ Definition	168	F _M ' Definition	169
F ₀ Mode	125	F ₀ Maximum Level.....	162
F ₀ ' Calculation.....	169	F ₀ ' Mode	72
F ₀ ' Definition	168		

F-Offset.....	73, 152	Internal PAR Sensor Settings	96
Freezing of MINI-PAM-II.....	164	Internal Temperature Sensor	160
Ft	167	Introduction	9
Ft Chart.....	52	J	
Ft Chart Resolution.....	98		
F _V /F _M Definition.....	170	JUNIOR-PAM Teaching Edition	107
G		K	
Gain.....	71, 153	KS-2500 Suspension Cuvette.....	37
Genty	175	KS-2500 Suspension Cuvette and	
Gilmore + Yamamoto.....	175	MKS-2500 Magnetic Stirrer.....	37
GPS-Status.....	45	L	
Graphics Windows.....	127	Leaf Clip / Ext. PAR Sett.	91
Grid View	118	Leaf Clip 2035-B	26, 91
Guarantee.....	189	Leaf Clip 60°	19
H		Leaf Clip Channel 1 Settings	92
Hardware	100	Leaf Clip Channel 2 Settings	94
Hints	161	Leaf Clip Connector	27
Humidity Sensor	29, 49	Leaf Positioning Setup DUAL-BA ...	34
I		Leaf Temperature	28
Ik	130	Leaf-Clip 2035-B	94
Indicators.....	154	LED	177
Induction Curve	56, 81, 128, 153	LED Signal Code	29, 155
Induction Curve Settings	81	LED Spectra.....	18
Info.....	100	Light Curve.....	57, 59, 82, 128, 154
Int. PAR sensor	157	Light Curve Settings.....	82
Integration Time.....	135	Light Intensity.....	151
Internal PAR	62, 95	Light List Calibration	63
Internal PAR Sensor.....	65	Light Panel	146
		Light Panel Settings	75, 76

Light Sources.....	74, 75, 76	MINI-PAM/F1 Miniature Fiberoptics	40
Limits	91	MINI-PAM-II Info	101
Line View	118	MINI-PAM-II Settings	97
Load Data	115, 118	MINI-SPEC/MP	21, 23, 25, 134
M		MKS-2500	187
Main Menu	66	ML-F High	71
Main Menu with all eight submenus.	67	Model Function REG1	129
Mark.....	47, 48	Model Function REG2.....	130
Maximum F_0 of a Dark-acclimated		Moni-Bus.....	113
Leaf.....	162	MQS-B/A Adapter Set.....	41
Measuring Light	69, 177	N	
Measuring Light and Signal	20	New Record	48, 121
Measuring Light Frequency.....	71, 72, 125	Non-photochemical Quenching.....	171
Measuring Light Intensity.....	71	NPQ Definition	171
Measuring Light Settings ...	70, 71, 152	O	
Measuring Light Trim Value.....	160	O ₂ System Overview	38
Memory.....	99, 125, 141	ON Button	15
Memory Window.....	139, 140	Online Data	122, 126
Menu and Tabs.....	115	Operating Temperature.....	178
Menu Overview.....	115	Options.....	119
Merge Data	142	Options Menu.....	137
Micro Quantum Sensor.....	182	Optode	86, 149
Micro-Quantum-Sensor	27	Optoelectronic Unit.....	177
Mini Quantum/Temp.-Sensor 2065-M		Optoelectronic Unit LED.....	18
.....	35, 185	Optoelectronic Unit Rear Panel.....	14
Miniature Fiberoptics	40, 95, 188	Optoelectronic Unit Sockets.....	15
Miniature Fiberoptics MINI-PAM/F1	39, 40	Output Format.....	87
Miniature Spectrometer	21, 23, 25, 88, 185	Overview Windows.....	44, 67, 68, 84
MINI-PAM Settings	98	Oxygen Sensor	86, 149

Oxygen Sensor Settings..... 86, 87

P

PAM Firmware Update 108, 109

PAM Fluorescence 165

PAM Mode (Act.) 76, 147

PAM Settings..... 69

PAM Trim Values..... 159

PAR Calibration 94

PAR Calibration Block 22, 24, 186

PAR Cosine 134

PAR Display 49

PAR Quantum Sensor 27

PAR Sensor 148

PAR Spectrometer..... 21

PAR Sphere..... 134

Photosynthetic Photon Flux Density 27

Pin Assignment..... 178

Plugins..... 115

Porometer..... 50

Power Supply 178, 179

PPFD 27

Primary Data..... 49

Principal Screen Layout of Top-level

Windows 45

Program Group..... 108

Program Settings..... 78

Program/Clock Settings..... 78

Programs 153

Pulse-amplitude Modulated

Fluorescence 165

PyroScience 86

Q

q_L Definition..... 170

q_N Definition 171

q_P Definition..... 170

Quench Data..... 122

Quenching Analysis 50, 51

Quit 115

R

Rapid Light Curve 31

Rear Panel of MINI-PAM-II 14

Rec. Online 118

Record 48

Record File..... 115

Record Macro 144

Recovery 60, 61

Recovery Curve 58

Red Version 177

References..... 175

Reflectance 23, 52

Reflectance Spectra..... 24

Reflectance Standard 24

REG1 128, 129

REG2 128, 130

Regression Data 116

Remove Marks..... 117

Report Table 136

Report Window 136

Reset..... 103

Reset Porometer..... 106

Reset Settings..... 103, 151

Reset Sys. Settings	105	Sensors Window	148
Results Sidebar	123, 124	Serial Number	100
Review Articles	173	Settings	115, 152
RGBW Light Source	29	Settings Window	150, 155
Ring Buffer	139	Setup	90
S		Sidebar	123
Safety Instructions	7	Signal LED	98
Sampling Frequency	118	Signal Noise	164
SAT Kinetics In Raw File	133	Sockets	15
SAT Pulse Data	122	Software Installation	107
SAT+Spec	135	SPEC Command	52
SAT-Chart Export	132	Specification Memory	178
SAT-Chart Options	132	Specification Touchscreen	178
SAT-Chart Window	131, 133, 135	Specifications	177
SAT-Pulse	125	Spectro. Oper. Mode	89
Saturation Pulse Analysis	166, 174	Spectrometer	52, 53, 134
Saturation Pulse Settings	77	Spectrometer Settings	88
Save Data	115, 118	Spectrum Window	132
Save Settings	115	Start Menu WinControl	108
Screen Calibration	43	Start/End Marks	117
Select Current Induction Curve	131	Status Bar	124, 125
Select Current Light Curve	131	Surface Clip	34
Select Display	83	Suspension Cuvette KS-2500	37, 187
Self-Locking Connector	17	Symbol View	118
Semicolon As Delimiter	116	SYNC	15
Sensor Info	101, 102	Sync with PC	141
Sensor Settings	83	System Hangs	164
Sensor Settings Menu and its		System Information	155
Submenus	84	System Parameter	153
Sensors	83	System Settings	156
		System Settings Button	156

T

Temperature	28, 49
Temperature Calibration	94
Temperature Units	115
Thermocouple	28
Thermocouple Error	149
Time	99
Time Format	117
Time Stamp	117
Title Bar	151, 157
Top Level Windows	44, 45
Touch Screen Calibration	43
Touchscreen Operation	43
Troubleshooting	161

U

Uninstall WinControl-3	111
Unterwater Cable	21
Update Indentation	142
USB Flashdrive	107
USB Socket	15
Used Cal. Profile	93
UTC Time	117

V

Val. Sidebar	122
Val./Rec./Chan.	121
van Kooten + Snel	175
View Menu	115

W

Warnings	115
Watertight Connection	21
WinControl-3	180
WinControl-3 Default Mode	111
WinControl-3 Download	107
WinControl-3 Installation	107
WinControl-3 Network Mode ..	112, 114
WinControl-3 Offline Mode	112
WinControl-3 Program Group	108
WinControl-3 Uninstall	111
WinControl-3 Windows	113

X

X-Axis Scaling	120
----------------------	-----

Y

Y(II) Definition	170
Y(NO) Definition	171
Y(NPQ) Definition	171
Y-Axis Scaling	120
Yield + # Poro	82

Z

Zoom to Selection	119
-------------------------	-----

A

α	130
----------------	-----

12 List of Figures

Fig. 1:	Rear Panel of MINI-PAM-II	14
Fig. 2:	Carrying Belt Fastening	16
Fig. 3:	Self-locking Connector.....	17
Fig. 4:	Fiberoptics	17
Fig. 5:	Emission Spectra of MINI-PAM II LEDs	18
Fig. 6:	Distance Clip 60° 2010-A	19
Fig. 7:	Measuring Light and Signal	20
Fig. 8:	Connection of MINI-SPEC/MP	21
Fig. 9:	PAR Calibration Block	22
Fig. 10:	Miniature Spectrometer MINI-SPEC/MP: Configuration	23
Fig. 11:	MINI-SPEC/MP	25
Fig. 12:	Emission Spectra of Miniature Spectrometer Light Sources	25
Fig. 13:	2035-B Leaf-Clip Holder	26
Fig. 14:	2054-L External LED Source	30
Fig. 15:	Emission spectra of the 2054-L External LED Source.....	30
Fig. 16:	2035-B/RLC Conversion Kit.....	31
Fig. 17:	Fiberoptics Adapter 90° 2030-B90: Mounting	32
Fig. 18:	2060-B & 2065-M Arabidopsis Clip with Quantum/Temp.-Sensor.....	33
Fig. 19:	Leaf Positioning Setup DUAL-BA	34
Fig. 20:	2060-A & 2065-M, Surface Holder with Quantum/Temp.-Sensor.....	34
Fig. 21:	Mini Quantum/Temp.-Sensor 2065-M	35
Fig. 22:	DLC-8 Dark Leaf Clip.....	36
Fig. 23:	KS-2500 Suspension Cuvette and MKS-2500 Magnetic Stirrer	37
Fig. 24:	O ₂ measurement. System Overview.....	38
Fig. 25:	BCS-9590 Barcode Scanner	39
Fig. 26:	Miniature Fiberoptics MINI-PAM/F1	39
Fig. 27:	MQS-B/A Adapter Set.....	41
Fig. 28:	Touch Screen Calibration	43
Fig. 29:	Principal Screen Layout of Top-level Windows	45
Fig. 30:	Basic Data, Change Marker.....	47
Fig. 31:	Primary Data.....	49

Fig. 32:	Quenching Analysis	51
Fig. 33:	Ft Chart.....	52
Fig. 34:	Spectrometer	53
Fig. 35:	Actinic + Yield	55
Fig. 36:	Induction Curve.....	56
Fig. 37:	Light Curve	59
Fig. 38:	Recovery.....	60
Fig. 39:	Actinic Light List.....	61
Fig. 40:	Calibrate Light List.....	63
Fig. 41:	PAR Error	63
Fig. 42:	Main Menu	66
Fig. 43:	Main Menu with all eight submenus.....	67
Fig. 44:	First Three Submenus in Detail	68
Fig. 45:	PAM Settings.....	69
Fig. 46:	Measuring Light Settings	70
Fig. 47:	Light Sources.....	75
Fig. 48:	Light Panel Settings.....	75
Fig. 49:	Program/Clock Settings.....	78
Fig. 50:	Clock Item.....	79
Fig. 51:	Actinic + Yield Settings	80
Fig. 52:	Induction Curve Settings.....	81
Fig. 53:	Light Curve Settings	82
Fig. 54:	Sensor Settings	83
Fig. 55:	Sensor Settings Menu and its Submenus.....	84
Fig. 56:	External PAR Sensor.....	85
Fig. 57:	Oxygen Sensor Settings	86
Fig. 58:	Settings.....	87
Fig. 59:	Output Format.....	87
Fig. 60:	Spectrometer Settings	88
Fig. 61:	Spectro. Oper. Mode	89
Fig. 62:	Setup	90
Fig. 63:	Limits	91
Fig. 64:	Leaf Clip / Ext. PAR Sett.....	91
Fig. 65:	Leaf Clip Channel 1 Settings	92

Fig. 66:	Used Cal. Profile	93
Fig. 67:	Leaf Clip Channel 2 Settings	94
Fig. 68:	External Temperature Sensor Settings.....	94
Fig. 69:	Internal PAR Sensor Settings	96
Fig. 70:	Overview: MINI-PAM-II, Memory, Info, and Reset Menus	97
Fig. 71:	MINI-PAM Settings	98
Fig. 72:	Memory	99
Fig. 73:	Info	100
Fig. 74:	MINI-PAM-II Info	101
Fig. 75:	Sensor Info	102
Fig. 76:	Firmware Info	102
Fig. 77:	Reset	103
Fig. 78:	PAM Firmware Update	109
Fig. 79:	Device Channel Configuration	110
Fig. 80:	Chart Window	114
Fig. 81:	Export Format for Reports	116
Fig. 82:	Chart Tools	120
Fig. 83:	Val. Sidebar	122
Fig. 84:	Model Function REG1	129
Fig. 85:	Model Function REG2	130
Fig. 86:	SAT-Chart Window	133
Fig. 87:	Spectrum Window.....	135
Fig. 88:	Memory Window	140
Fig. 89:	Batch Window Overview.....	145
Fig. 90:	Control Window	146
Fig. 91:	Sensors Window	148
Fig. 92:	Settings Window	150
Fig. 93:	System Settings Button	156
Fig. 94:	System Settings.....	156
Fig. 95:	Fluorescence Kinetics Induced by Saturation Pulse.....	163
Fig. 96:	Illustration of the PAM measurement principle	166
Fig. 97:	Fluorescence Levels of Saturation Pulse Analysis	167

13 List of Tables

Table 1:	Rear Panel of MINI-PAM-II	15
Table 2:	Signal Code of LED on 2035-B Leaf Clip Holder	29
Table 3:	Miniature Fiberoptics MINI-PAM/F1	40
Table 4:	Overview of Top-level Windows	44
Table 5:	GPS-Status	45
Table 6:	Sequence of Saturation Pulse Analyses in a Recovery Curve	61
Table 7:	The CAL Command	64
Table 8:	Measuring Light Frequencies	72
Table 9:	Default Settings	103
Table 10:	WinControl-3 in Windows Start Menu	108
Table 11:	Windows of WinControl-3	113
Table 12:	Menu Overview	115
Table 13:	Export Format for Reports	116
Table 14:	Buttons	118
Table 15:	Options	119
Table 16:	Sidebar	123
Table 17:	Status Bar	125
Table 18:	Three Graphics Windows	127
Table 19:	SAT Kinetics in Raw File	133
Table 20:	Report Table	136
Table 21:	Options Menu	137
Table 22:	Abbreviations for Settings	138
Table 23:	External Light Source	147
Table 24:	LED Signal Code	155
Table 25:	System Information	155
Table 26:	Custom PAR Gain	158
Table 27:	PAM Trim Values	159
Table 28:	Maximum F_0 of a Dark-acclimated Leaf	162
Table 29:	Fluorescence Ratio Parameters	174
Table 30:	References Cited in Table 29	175
Table 31:	Pin Assignments. View from Outside	178
